Gebrauchsanleitung | Operating manual |  
Mode d'emploi | Instrucciones de manejo |  
Istruzione | Instruções | 操作手册

## macro

Pipettierhelfer | Pipette controllers

# Impressum

BRAND GMBH + CO KG

Otto-Schott-Str. 25

97877 Wertheim (Germany)

T +49 9342 808 0

F +49 9342 808 98000

[info@brand.de](mailto:info@brand.de)

[www.brand.de](http://www.brand.de)

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# Table of Contents

<b>1</b>	<b>Terms of use</b> .....	<b>28</b>
1.1	Hazard levels .....	28
1.2	Operating manual symbols .....	28
1.3	Format .....	29
<b>2</b>	<b>Safety</b> .....	<b>30</b>
2.1	Function.....	31
2.2	Limitations of use.....	31
2.3	Operating exclusions .....	31
2.4	Storage Conditions .....	31
<b>3</b>	<b>Functional and operating elements</b> .....	<b>32</b>
<b>4</b>	<b>Pipetting</b> .....	<b>33</b>
<b>5</b>	<b>Cleaning</b> .....	<b>37</b>
5.1	Autoclaving.....	37
5.2	Replace filter and clean adapter .....	37
5.3	Replace valve system .....	39
5.4	Check tightness .....	40
<b>6</b>	<b>Malfunction – what to do?</b> .....	<b>41</b>
<b>7</b>	<b>Product markings</b> .....	<b>42</b>
<b>8</b>	<b>Ordering Information</b> .....	<b>43</b>
<b>9</b>	<b>Spare parts</b> .....	<b>44</b>
<b>10</b>	<b>Repairs</b> .....	<b>46</b>
10.1	Sending for repair .....	46
<b>11</b>	<b>Warranty</b> .....	<b>48</b>
<b>12</b>	<b>Disposal</b> .....	<b>49</b>

# 1 Terms of use

- Please carefully read the operating manual before using the device for the first time.
- The operating manual is part of the device and must be kept in an easily accessible place.
- Be sure to include the operating manual if you transfer possession of this device to a third party.
- You can find up-to-date versions of the operating manual on our website: [www.brand.de](http://www.brand.de).

## 1.1 Hazard levels

The following signal words identify possible hazards:

Signal word	Meaning
DANGER	Will lead to serious injury or death.
WARNING	May lead to serious injury or death.
CAUTION	May lead to minor or moderate injuries.
NOTICE	May lead to property damage.

## 1.2 Operating manual symbols

Symbol	Meaning	Symbol	Meaning
	Danger area		Wear protective gloves
	Wear eye protection		

## 1.3 Format

Format	Meaning	Icon	Meaning
<b>1. Task</b>	Indicates a task.	>	Indicates a condition.
a., b., c.	Indicates the individual steps of a task.	⇒	Indicates a result.

## 2 Safety

### **Please read carefully!**

The macro can be used in combination with hazardous materials, work processes and equipment. However, the operating manual cannot cover all of the safety issues that may occur in doing so. It is the user's responsibility to ensure compliance with the safety and health regulations and to specify the corresponding restrictions before use.

1. Every user must read and observe this operating manual before using the device.
2. Follow the general hazard instructions and safety regulations (e.g., wear protective clothing, eye protection, and protective gloves).
3. Follow the instructions given by the reagent manufacturer.
4. Use the device only for pipetting liquids and only within the defined limits of use. Comply with the operating exclusions; see Operating exclusions, p. 31 . In case of doubt, contact the manufacturer or dealer.
5. Always perform work in a manner that does not endanger the user or other people. Avoid splattering. Use only suitable vessels.
6. Never use force.
7. Use only original accessories and original replacement parts. Do not make any technical modifications. Do not disassemble the device further than described in the user manual!
8. Always check that the device is in proper working condition before use. If the device malfunctions, stop pipetting immediately and refer to the section Malfunction – what to do?, p. 41 . If necessary, contact the manufacturer.

## 2.1 Function

The device serves as an aid for pipetting liquids with graduated and volumetric pipettes made of glass or plastic in the volume range from 0.1 ml to 200 ml and with a suction tube outer diameter < 9.2 mm. When the device is handled correctly, the pipetted liquid comes into contact with the pipette only.

## 2.2 Limitations of use

This device is intended for pipetting liquids within the following physical limitations:

- from 10 to 40°C (from 50 to 104°F) (device and reagent)
- vapor pressure up to 500 mbar. Absorb slowly above 300 mbar in order to prevent the liquid from boiling.
- Density up to 2.4 g/cm<sup>3</sup>

## 2.3 Operating exclusions

The device must not be used for liquids with vapors that can attack the materials silicone or PP.

The device is not suitable for use with Pasteur pipettes.

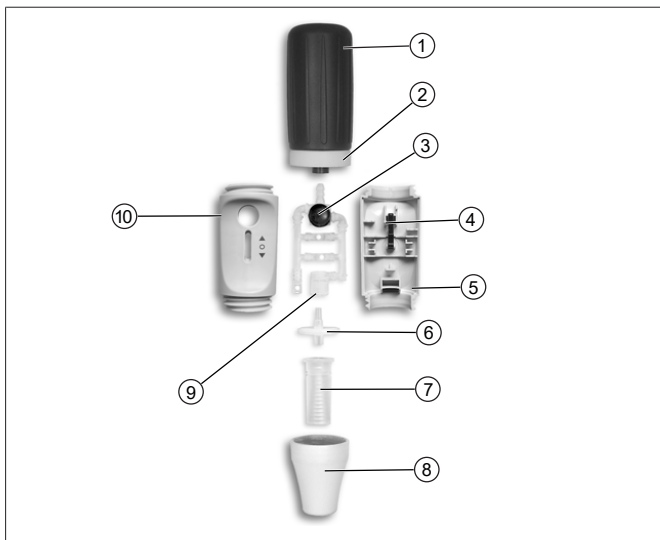
## 2.4 Storage Conditions

Store the unit and accessories in a cool and dry place in cleaned condition only.

Storage temperature from -20 to 50°C (-4 to 122°F).

Relative humidity: 5 % to 95 %.

## 3 Functional and operating elements

1	Suction bellows (SI)	6	Membrane filter (PP/PTFE)
2	Screw ring (PP)	7	Adapter (SI)
3	Blow-out bubble (SI)	8	Adapter housing (PP)
4	Pipetting lever	9	Valve system (PP; SI, PTFE)
5	Housing (PP)	10	Housing (PP)

Abbreviation	Material
PP	Polypropylene
PTFE	Polytetrafluoroethylene
SI	Silicone

## 4 Pipetting

### 1. Vent the suction bellows

1.

- a. Press the suction bellows together before inserting the pipette.

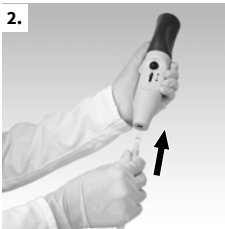
### 2. Attach pipette

#### WARNING

Ensure that the pipette is firmly in place. Never use force! Especially with thin pipettes, there is a risk of glass breakage!  
Risk of injury!

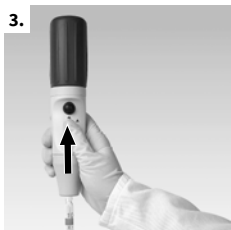
2.

- a. Always hold the pipette as close as possible to the upper end and push it carefully into the adapter until it is firmly in place.
- b. Always hold the device with the pipette inserted vertically, tip down.

### 3. Fill pipette

3.

- a. Immerse pipette tip into the test liquid.
- b. Carefully push the pipetting lever upwards, and fill the pipette slightly above the desired mark.

#### NOTICE

Make sure that no liquid gets into the device! The function of the membrane filter is impaired, and the suction capacity decreases. In this case, replace membrane filter; see Replace filter and clean adapter, p. 37 .

#### NOTICE

The further the lever is moved upwards, the stronger the suction effect (i.e., the faster the meniscus rises)!

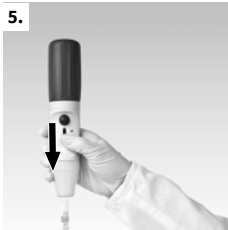
With large pipettes (> 50 ml), the vacuum in the suction bellows is not sufficient to draw up the entire amount of liquid at once. In this case: Press the suction bellows together again and continue to pull it up.

### 4. Set volume

- a. If necessary, wipe the tip of the pipette with a suitable lint-free cloth.
- b. Press the pipetting lever down gently until the meniscus is precisely adjusted.

## 5. Empty pipette

5.

- a. Hold collecting vessel at an angle. Place the pipette tip against the inner wall and press the pipetting lever down.

**For pipettes with waiting time** (imprint e.g., 'Ex + 5 s'):

As soon as the meniscus in the pipette comes to a standstill, observe the waiting time indicated on the pipette (here 5 s).

Pull up the tip of the pipette a few millimeters along the wall of the vessel.

### NOTICE

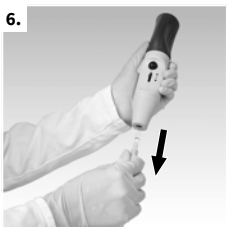
#### When using blow-out pipettes manufactured by others

As soon as the meniscus in the pipette tip has come to a standstill, press the small bellows once. Pull up the tip of the pipette a few millimeters along the wall of the vessel.

## 6. After pipetting

6.

- a. Grip the pipette as high up as possible, and pull it out of the adapter.

## 5 Cleaning

Occasionally clean the housing parts with a damp cloth. When used properly, the device is maintenance-free.

### 5.1 Autoclaving

The device must be carefully cleaned before it is autoclaved. Clean the housing parts with a damp cloth. The device can be autoclaved at 121°C (250°F), 2 bar, and a holding time of at least 15 minutes according to DIN EN 285.

### 5.2 Replace filter and clean adapter

#### ⚠ WARNING

#### Risk of contamination and infection

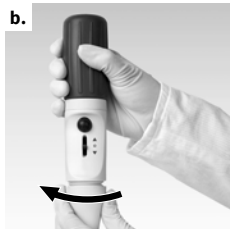
Wear protective gloves and goggles.

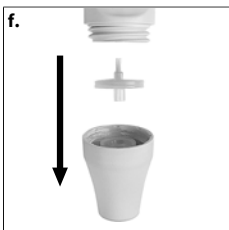
#### NOTICE

Only necessary if a pipette has been overfilled or the suction capacity is impaired.

**b.**

- a.** Pull out pipette.
- b.** Unscrew the adapter housing.
- c.** Remove defective membrane filter.
- d.** Press the adapter upwards out of the adapter housing.

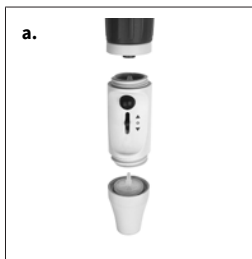
- e.** For cleaning, rinse the adapter carefully with a spray bottle filled with a suitable cleaning liquid (e.g., water). Then blow out and dry.
- f.** First press the new membrane filter into the adapter with the thicker connection facing downwards (the thinner, conical side facing upwards).
- g.** Unscrew the adapter by hand.
- h.** The complete device, including membrane filter, can be autoclaved at 121°C (2 bar) according to DIN EN 285.
- i.** Unscrew the suction bellows for autoclaving.

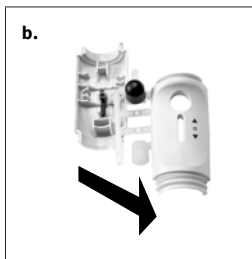
#### NOTICE

Not every membrane filter is suitable! Use only original accessories. The device comes standard with filters of 3 µm pore width. The membrane filter can be autoclaved up to five times at 121°C. The effectiveness of autoclaving must be verified by the user.

## 5.3 Replace valve system

- a.** Unscrew suction bellows and adapter housing.
- b.** Remove the upper part of the housing for replacement. Carefully remove and replace the valve system.
- c.** Reassemble the unit in reverse order and check for leaks.

## 5.4 Check tightness

After adjusting the meniscus, no liquid should drip from the pipette without operating the valves. If the pipette drips; see Malfunction – what to do?, p. 41

## 6 Malfunction – what to do?

Fault	Possible cause	What to do
Suction capacity impaired	Filter dirty or valve system defective	Replace filter, Replace filter and clean adapter, p. 37 .
Pipette drips	Filter not placed correctly or valve system defective	Insert filter correctly, or replace valve system; see Replace filter and clean adapter, p. 37 .
Pipette does not hold	Silicone adapter damaged	Unscrew the lower part of the device, and replace the silicone adapter
Suction bellows ventilates independently	Suction bellows or valve system defective	Replace upper screw connection with suction bellows or valve system

### NOTICE

A visual inspection of the device is often sufficient to detect a defective part. This can then be easily replaced with a spare part. The advantages of self-repair are obvious: Cost and time savings.

## 7 Product markings

Symbol or number	Meaning
	General warning signs
	Follow the operating instructions
XXZXXXXX	Batch number

## 8 Ordering Information

### macro pipette aid

Description	Color*	Order no.
macro pipetting aid, including. Replacement membrane filter (3 µm) and operating instructions.	gray	262 00
	Green	262 01
	Blue	262 02
	magenta	262 03

\* The colors refer to the adapter housing and screw ring.

## 9 Spare parts

### Adapter housing

Color*	Length (mm)	Material	Order no.
gray	49	PP	262 20
Green	49	PP	262 23
Blue	49	PP	262 21
magenta	49	PP	262 22

### Suction bellows with screw ring

Color*	Order no.
gray	262 25
Green	262 28
Blue	262 26
magenta	262 27

\* The colors refer to the adapter housing and screw ring.

### Replacement membrane filter (3 µm)

Packaging unit	Order no.
1 pc in blister packaging	260 52

### Replacement membrane filter (3 µm)

Packaging unit	Order no.
10 pc in blister packaging	260 56

## Silicone adapter

Length (mm)	Order no.
44	261 46

## Valve system

Description	Order no.
Valve system	261 28

# 10 Repairs

## 10.1 Sending for repair

### NOTICE

Transporting of hazardous materials without a permit is a violation of federal law.

---

### Clean the device thoroughly and decontaminate!

- When returning products, please enclose a general description of the type of malfunction and the media used. The device cannot be repaired if information about the media used is not provided.
- The device is returned at the risk and expense of the sender.

### Outside USA and Canada

Fill out the “Declaration of Health Safety” and send it together with the device to the manufacturer or dealer. Pre-printed forms can be requested at the dealer or manufacturer, or are available for download at [www.brand.de](http://www.brand.de).

### Within USA and Canada

Please clarify the requirements for the return delivery with BrandTech Scientific, Inc **before** sending the device in for service.

Send only cleaned and decontaminated devices to the address, which you received together with the return number. Attach the return number in a clearly visible place on the package.

### Contact addresses

**Germany:**  
BRAND GMBH + CO. KG

**USA and Canada:**  
BrandTech® Scientific, Inc.

Otto-Schott-Str. 25  
97877 Wertheim (Germany)  
T +49 9342 808 0  
F +49 9342 808 98000  
info@brand.de  
www.brand.de

**India:**

BRAND Scientific Equipment Pvt. Ltd.  
303, 3rd Floor, 'C' Wing, Delphi  
Hiranandani Business Park,  
Powai  
Mumbai-400 076 (India)  
T +91 22 42957790  
F +91 22 42957791  
info@brand.co.in  
www.brand.co.in

11 Bokum Road  
Essex, CT 06426-1506 (USA)  
T +1-860-767 2562  
F +1-860-767 2563  
info@brandtech.com  
www.brandtech.com

**China:**

BRAND (Shanghai) Trading Co., Ltd.  
Guangqi Culture Plaza  
Room 506, Building B  
No. 2899, Xietu Road  
Shanghai 200030 (P.R. China)  
T +86 21 6422 2318  
F +86 21 6422 2268  
info@brand.com.cn  
www.brand.cn.com

## 11 Warranty

We shall not be liable for the consequences of improper handling, use, servicing, operating or unauthorized repairs of the device or for the consequences of normal wear and tear, especially of wearing parts such as pistons, seals, valves and the breakage of glass. The same applies for failure to follow the instructions of the operating manual. We are not liable for damage resulting from disassembly beyond that described in the operating manual or if non-original spare parts or components have been installed.

### **USA and Canada:**

Find more warranty information on [www.brandtech.com](http://www.brandtech.com).

## 12 Disposal

Before disposal, observe the relevant national disposal regulations and ensure that the product is disposed of properly.

Subject to technical changes, errors, and misprints.

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BRAND Scientific Equipment Pvt.Ltd.