[Gebrauchsanleitung](#) | [Operating manual](#) | [Mode d'emploi](#) | [Instrucciones de manejo](#) | [Istruzione](#) | [Instruções de utilização](#) | [操作手册](#) | [Руководство по эксплуатации](#) | [사용 지침](#) | [Bruksanvisning](#) | [使用説明書](#) | [Használati utasítás](#) | [Návod k použití](#) | [Gebruiksaanwijzing](#) | [Instrukcja użytkowania](#) | [Kullanım Talimatları](#)

accu-jet[®] S

Pipettierhelfer | Pipette controllers

Impressum

BRAND GMBH + CO KG

Otto-Schott-Str. 25
97877 Wertheim (Germany)

T +49 9342 808 0
F +49 9342 808 98000
info@brand.de
www.brand.de

Do you need more operating manuals and translations?

Please refer to <http://www.brand.de/om> or use the following

QR Code:

The original operating manual is written in German. Other languages are translations of the original operating manual. Technical changes, errors or misprints reserved.

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This document declares the accordance with the named harmonized regulations, but does not assure specific properties.

UK Declaration of Conformity

Product name: Pipetting controller accu-jet® S
Product variations: anthracite, berry, petrol, amethyst
Cat.-No. 26350, 26351, 26352, 26353, 26360, 26361, 26362, 26363

The product described above meets the basic requirements that are specified in the harmonization legislation listed below:	Applied harmonized standards:
The Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic Equipment Regulations 2012 (SI 2012 No. 3032)	EN IEC 63000:2018
Electromagnetic Compatibility Regulations 2016 (SI 2016 No. 1091, as amended by SI 2019 No. 696)	EN 61326-1:2013
Other applied standards:	Scope:
EN 61010-1:2010, EN 61010-1:2010/A1:2019/AC:2019-04 EN/IEC 62368-1:2014 (angewendet auf Fremdherstellernetzteil / applied to third-party power supply)	LVD LVD

The manufacturer is solely responsible for issuing this declaration of conformity with regard to meeting the essential requirements and preparing the technical documentation.

Manufacturer	Address
BRAND GMBH + CO KG	Otto-Schott-Str. 25 97877 Wertheim Germany

Wertheim, 18 February 2022

Patrick Ziemeck
Managing Director Technology

i.A. Siegfried Ott
Regulatory Affairs

This document declares the accordance with the named harmonized regulations, but does not assure specific properties.

Declaration of conformity – China RoHS 2

BRAND GMBH + CO KG has made reasonable efforts to ensure that restricted hazardous substances are not used in BRAND products above the limits specified by China RoHS 2.

In order to determine the concentration of hazardous substances in all homogeneous materials of the subassemblies, a “Product Conformity Assessment” (PCA) procedure was performed. As defined in GB 26572-2025 the “Maximum Concentration Value” limits (MCV) apply to these restricted substances:

+ Lead (Pb):	0.1 %	+ Hexavalent chromium (Cr(+VI)):	0.1 %
+ Mercury (Hg):	0.1 %	+ Polybrominated biphenyls (PBB):	0.1 %
+ Cadmium (Cd):	0.01 %	+ Polybrominated diphenyl ether (PBDE):	0.1 %
+ Bis(2-ethylhexyl) phthalate (DEHP):	0.1 %	+ Dibutyl phthalate (DBP):	0.1 %
+ Butyl benzyl phthalate (BBP):	0.1 %	+ Diisobutyl phthalate (DIBP):	0.1 %

Environmental Friendly Use Period (EFUP)

EFUP defines the period in years during which the hazardous substances contained in electrical and electronic products will not leak or mutate under normal operating conditions. During normal use by the user such electrical and electronic products will not result in serious environmental pollution, cause serious bodily injury or damage to the user's assets.

The Environmental Friendly Use Period for BRAND instruments is 40 years.

Material Content Declaration for BRAND Products

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Apart from the disclosures in the above table, the subassemblies are not intentionally manufactured or formulated with lead (Pb), mercury (Hg), cadmium (Cd), hexavalent chromium (Cr+VI), polybrominated biphenyls (PBB), polybrominated diphenyl ethers (PBDE), bis(2-ethylhexyl) phthalate (DEHP), butyl benzyl phthalate (BBP), dibutyl phthalate (DBP) and diisobutyl phthalate (DIBP).

Products manufactured by BRAND may enter into further devices or can be used together with other appliances. With these third party products and appliances in particular, please note the EFUP labeled on these products. BRAND will not take responsibility for the EFUP of those products and appliances.

Place, date: Wertheim, 15|07|2025

Patrick Ziemeck
(Managing Director Technology)

i.A. Siegfried Ott
(Regulatory Affairs)

Table of contents

1 Introduction	45
1.1 Using the operating manual.....	45
2 Safety	46
2.1 Function.....	48
2.2 Limitations of use.....	48
2.3 Operating exclusions	48
2.4 Storage conditions.....	48
3 Scope of delivery	50
4 Functional and operating elements	51
5 Connecting the battery	53
6 Charging the battery	56
7 Turning on the instrument	58
8 Selecting the operating mode	59
9 Aspirating or dispensing liquid	60
10 Setting the motor speed	61
11 Changing the pipetting speed.....	62
12 Resting the instrument, mounting the wall mount.....	63
13 Pipetting	64
14 Cleaning.....	66
14.1 Replace filter and clean adapter	66
14.2 Check tightness.....	68
14.3 UV sterilization	68
15 Troubleshooting	69
16 Technical data	70
16.1 Product markings	70

17 Ordering Information	72
18 Repairs.....	74
18.1 Sending for repair	74
19 Warranty	76
20 Disposal	77

1 Introduction

1.1 Using the operating manual

- Please carefully read the operating manual before using the device for the first time.
- Keep the operating manual in an easily accessible place. It is part of the instrument.
- Be sure to include the operating manual if you transfer possession of this device to a third party.

1.1.1 Signal words and their meaning

⚠ WARNING The signal word WARNING indicates a hazardous situation that, if not avoided, could result in death or serious injury.
or **⚠WARNING! ...**

⚠ CAUTION The signal word CAUTION indicates a hazardous situation that, if not avoided, could result in minor or moderate injury.
or **⚠CAUTION! ...**

NOTICE The signal word NOTICE is used to address practices not related to physical injury. Example: Possible property damage.
or **NOTICE ...**

1.1.2 Operating manual symbols

Symbol	Meaning	Symbol	Meaning
	Hazardous area		Explosion hazard
	Wear eye protection		Wear protective gloves

1.1.3 Presentation of descriptions of actions

Format	Meaning
1. Task	Indicates a task.
a., b., c.	Indicates the individual steps of a task.
>	Indicates a prerequisite for a task.
⇒	Indicates a result of a completed task.

2 Safety

Please read carefully!

The accu-jet® S can be used in combination with hazardous materials, work processes and equipment. However, the operating manual cannot cover all of the safety issues that may occur in doing so. It is the user's responsibility to ensure compliance with the safety and health regulations and to specify the corresponding restrictions before use.

- Every user must read and observe the operating manual enclosed with this laboratory device before and during use. The device may only be used by trained and qualified personnel.
- Follow the general hazard instructions and safety regulations (e.g. wear protective clothing, eye protection and protective gloves).
- When working with infectious or hazardous samples/media (e.g. hazardous substances), the general safety rules in the laboratory must be observed and regulations for handling the samples/media must be followed. The information provided by the media manufacturer (e.g. safety data sheets) must be observed.
- Use the laboratory device only for pipetting or dispensing media within the defined limitations and restrictions of use. Comply with the operating exclusions.
- If working with inflammable media, make sure to take the necessary precautions to avoid the buildup of static electricity (e.g., do not dispense into plastic vessels and do not wipe instruments with a dry cloth). Do not use the device in potentially explosive atmospheres. If in doubt, contact the manufacturer or supplier.
- Always check that the instrument is in proper working condition before use. If device malfunctions are indicated (e.g. sluggish pistons, leaks or power supply faults), stop working with the device immediately and refer to the "Troubleshooting" section in the operating manual. Contact the manufacturer, if necessary.

- Always perform work in a manner that does not endanger users or other people. Avoid splashes. Only use suitable vessels. Never use unnecessary force when operating, cleaning or servicing the laboratory device.
- If the laboratory device is supplied with power through a power adapter, batteries or accumulators, the components and connections on the device must be regularly checked for proper working condition. Do not operate the laboratory device and its accessories in an unprotected, humid or wet environment.
- Do not make any technical modifications. Use only original replacement parts of the manufacturer; even power adapters or accumulators of identical size and specification from other manufacturers are not permitted. Do not disassemble the laboratory device and its accessories (e.g., power adapters, cables, racks, accumulators or batteries) any further than described in the operating manual!
- Only autoclave the laboratory device when this is permissible according to the operating manual.
- Use only the original power adapter to charge the nickel-metal hydride battery.
- Protect the power adapter from moisture and only use it in conjunction with this device.
- Stop using the power adapter if the connection cable is damaged.
- Only authorized service technicians are permitted to carry out repairs and internal maintenance on the instrument.

WARNING

Potential risk of explosion due to damaged battery

In extreme cases, improper handling of the instrument or the battery (short circuit, mechanical damage, overheating, etc.) may cause the battery to explode.

2.1 Function

The device serves as an aid for pipetting liquids with graduated and volumetric pipettes made of glass or plastic in the volume range from 0.1 ml to 200 ml and with a suction tube outer diameter < 9.2 mm. When the device is handled correctly, the pipetted liquid comes into contact with the pipette only.

2.2 Limitations of use

This device is intended for pipetting liquids within the following physical limitations:

- from 10 to 40°C (from 50 to 104°F) (device and reagent)
- vapor pressure up to 500 mbar. Absorb slowly above 300 mbar in order to prevent the liquid from boiling.
- Density up to 9 g/cm³

2.3 Operating exclusions

The instrument must not be used for liquids with vapors that can cause corrosion or attack the materials silicone or EPDM.

The instrument is not suitable for use with Pasteur pipettes.

WARNING

Operating the instrument in a potentially explosive atmosphere

The instrument must never be operated or charged in a potentially explosive atmosphere. Pipetting highly flammable media (flash point under 0 °C (32 °F), e.g. ether, acetone) is prohibited.

2.4 Storage conditions

Store the unit and accessories in a cool and dry place in cleaned condition only.

Storage temperature from -20 to 50°C (-4 to 122°F).

Relative humidity: 5 % to 95 %.

3 Scope of delivery

Pipette controller, nickel-metal hydride battery, power adapter optional (note different item numbers), 2 replacement membrane filters 0.2 µm, sterile, wall mount, operating manual

4 Functional and operating elements

- | | |
|---|---|
| 1 Upper housing shell | 2 Upper housing shell base |
| 3 Motor speed adjustment wheel | 4 Operating mode selection switch |
| 5 Power adapter | 6 Country adapter |
| 7 Upper pipetting button (aspirate liquid) | 8 Bottom pipetting button (dispense liquid) |
| 9 Handle base | 10 Charging socket |
| 11 Adapter housing (PP) | 12 Silicone adapter |
| 13 Membrane filter (PP/PTFE) | 13 NiMH battery pack |
| 15 Pressure compensation opening | 14 For voltage and current consumption specifications, see Technical data, p. 70 . |

4 Functional and operating elements

Abbreviations	Material
PP	Polypropylene
PTFE	Polytetrafluoroethylene
SI	Silicone

5 Connecting the battery


a.

a. Remove the instrument from the packaging. Unscrew the adapter housing (1).

⇒ The adapter and the membrane filter remain in the adapter housing.

b.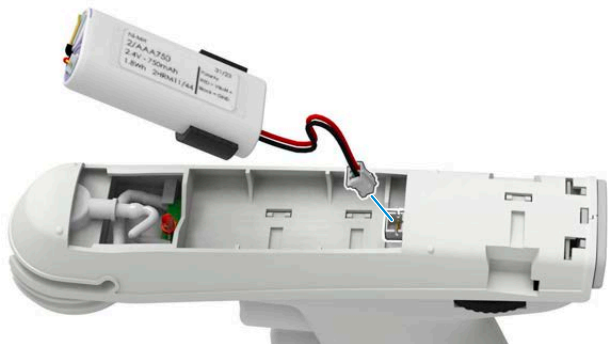

- b. Remove the battery from the packaging. Grip battery plug by the wire, confirm alignment of the plug and insert it into the socket. Push it firmly into the socket until you can hear it snap into place. Insert the battery into the compartment.

c.

- c. Remove the housing upper section from the packaging. Attach it carefully. Attach it carefully. While doing so, make sure that the upper housing section engages in the designated slots above the adjustment wheel and in the screw connection of the adapter housing. Then carefully slide the housing upper section toward the selection switch, as shown in the image.

d.

- d. Screwing on the adapter housing

- e. Charge the battery until the LED charge indicator is flashing green. Charging time is approx. 4 hours, see Charging the battery, p. 56 .

6 Charging the battery

⚠WARNING! Risk of injury from electric shock due to a damaged power adapter, charging cable, or country adapter. Do not use damaged power adapters, charging cables, or country adapters.

NOTICE Using another power adapter can damage the instrument and the battery.

- a. Place the country adapter onto the power adapter.
 - b. Connect the power adapter to the instrument.
 - c. Plug the power adapter into the power outlet.
- ⇒ The instrument is charging. This is indicated by the LED charge indicator on the upper housing section. For operation and charging duration, see Technical data, p. 70 .

LED charge indicator status

Flashing green. Battery is charged. Trickle charging is in progress. This minimizes the lazy battery effect. To utilize maximum battery capacity, only recharge when the LED charge indicator lights up.

Solid green. The instrument is charging. You can continue working while it charges.

Flashing red. Battery is almost empty. Charge the battery. You can also continue to work while charging.

Solid red. Battery is empty. Charge the battery.

7 Turning on the instrument

The instrument is ready for operation when a charged battery is inserted. There is no ON/OFF switch.

8 Selecting the operating mode

Selector switch, see Functional and operating elements, p. 51 .

Select between the following modes to dispense fluid:

Operating mode	
Gravity delivery	slide the selection switch upward. Press the lower pipette button and fluid will leak out of the pipette.
“Blow-out with motor support” operating	slide the selection switch downward. If you press the lower pipette button, fluid is blown out of the pipette using the motor.

9 Aspirating or dispensing liquid

Pipetting button, see Functional and operating elements, p. 51

To fill the pipette, press the upper pipetting button; to dispense, press the lower pipetting button.

10 Setting the motor speed

Adjustment wheel, see Functional and operating elements, p. 51

The motor speed can be infinitely adjusted from maximum (+) to minimum (-) by turning the adjustment wheel. We recommend using the maximum motor speed for large volume pipettes and a lower motor speed for small volume pipettes.

11 Changing the pipetting speed

The speed is limited by the maximum motor speed set. The further the upper and lower pipetting buttons are pressed in, the faster the instrument aspirates or dispenses liquid.

12 Resting the instrument, mounting the wall mount

Resting on a level surface

The accu-jet® S can be set down on an even surface with an inserted pipette.

Mounting the wall mount

- Clean the mounting surface with a suitable cleaner (no moisturizing household cleaners) and a lint-free cloth, and allow it to dry thoroughly.
- Remove the wall mount (1) from the packaging. Peel the protective film off the adhesive strip on the back of the wall mount.
- Firmly** press the wall mount onto the mounting surface.

13 Pipetting

1. Attach the pipette

1.

- a. Always hold the pipette as close as possible to the upper end and push it carefully into the adapter until it is firmly in place.

⚠WARNING! Risk of injury!

Never use force! Especially with thin pipettes, there is a risk of glass breakage! Ensure that the pipette is firmly in place.

- b. Always hold the instrument with the pipette inserted vertically, tip down.

2. Fill the pipette

2.


- a. Immerse pipette tip into the test liquid.
- b. Press the upper pipetting button slowly and fill the pipette so that the meniscus lies slightly above the desired marks.
- c. You can aspirate or dispense liquid with greater precision by pressing both pipetting buttons. This is particularly helpful when adjusting the meniscus or when using small volume pipettes.

NOTICE Do not overfill the pipette.

3. Set volume

3.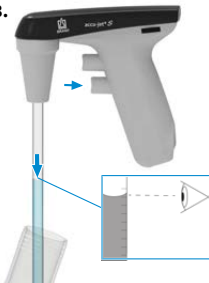

- a. If necessary, wipe the tip of the pipette with a suitable lint-free cloth.
- b. Press the lower pipetting button slowly and dispense liquid until the meniscus is adjusted exactly to the desired volume.

4. Empty the pipette

- a. Press the lower pipetting button slowly. The pipetting speed depends on how far the pipetting buttons have been pressed in. In the “Blow-out with motor support” operating mode, the speed also depends on the set motor speed.

5. After pipetting

- a. Grasp the pipette as high as possible and pull it out of the adapter.

14 Cleaning

NOTICE accu-jet® S is not autoclavable

Occasionally clean the housing parts with a damp cloth. When used properly, the instrument is maintenance-free.

14.1 Replace filter and clean adapter

⚠WARNING! Risk of contamination and infection when replacing the filter and cleaning the adapter

Wear protective gloves and eye protection.

NOTICE Filter replacement or cleaning is only necessary if a pipette has been overfilled or the suction capacity is impaired.

b.

- a. Pull out pipette.
- b. Unscrew the adapter housing.
- c. Pull out the upper housing.

d.

- d. Push out the adapter and membrane filter from the tapered side of the adapter housing.

e.

e. Remove defective membrane filter.


f.

f. For cleaning, rinse the adapter carefully with a wash bottle filled with a suitable cleaning liquid (e.g., water). Then blow out and dry.

g.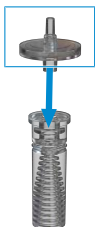

g. First press the new membrane filter into the adapter with the thicker connection facing downwards (the thinner, conical side facing upwards) and insert it into the adapter housing.

h.

h. Press the upper part of the housing into the housing shells and screw on the adapter housing.

NOTICE Autoclavable filters:

The effectiveness of autoclaving must be verified by the user. Not every membrane filter is suitable! Use only original accessories:

- The instrument comes equipped with a 0.2 μm membrane filter (red color code). The membrane filter can be autoclaved up to five times at 121°C (250 °F). Class B volume tolerances can be achieved with 0.2 μm membrane filters (for cell cultures).
- 3 μm membrane filters are required to comply with Class A volume tolerances. Due to the larger pore size, these have a reduced impact on the free delivery time of the pipettes, see .

14.2 Check tightness

After adjusting the meniscus, no liquid should drip from the pipette.

- Adjust the meniscus.
- Hold the instrument vertically.
- If a drop forms after a short waiting time (30 s), follow the remedy under Malfunction - What to do?

14.3 UV sterilization

The device is resistant to normal exposure to a UV disinfection lamp. The effects of the UV exposure may cause some color change.

15 Troubleshooting

Error	Cause	Corrective action
Suction capacity impaired	Motor speed is reduced	Increase motor speed, see Setting the motor speed, p. 61
Pipette dripping	Filter not properly positioned	Insert filter correctly, see Replacing the filter and cleaning the adapter, p. 66 .
	Adapter or pipette damaged	Replace the adapter or pipette, see Replacing the filter and cleaning the adapter, p. 66 .
Pipette not held in adapter	Adapter dirty or damaged	Clean adapter, allow to air dry or replace if necessary, see Replacing the filter and cleaning the adapter, p. 66 .
LED charge indicator does not light up/blink after the power adapter is connected and power is supplied.	Battery plug not firmly inserted	Firmly insert the battery plug, see Connecting the battery, p. 53 .
	Wrong power adapter	Use only original power supply unit, see Functional and operating elements, p. 51
	Battery or power adapter defective	Replace battery/power adapter
Low battery capacity	Instrument defective	Send instrument in for repair, see Repairs, p. 74 .
	Battery defective	Replace the battery, see Connecting the battery, p. 53 .
The instrument does not show any function	Internal error (perform reset)	Remove the battery plug and hold a pipetting button for 10 seconds. Then firmly insert the battery plug again, see Connecting the battery, p. 53 and Pipetting, p. 64

If the troubleshooting table above does not solve the problem, contact your authorized dealer or the manufacturer directly.

16 Technical data

Technical data	Values
Charging socket	5 VDC
Power adapter	Input: 100 – 240 VAC, 50/60 Hz Output: 5 VDC
Battery	Battery pack NiMH 2.4 V, 730 mAh Charging time: 4 h Maximum operating time: approx. 8 h Charging temperature: +10 °C to +35 °C

16.1 Product markings

Symbol or number	Meaning
XXZXXXXX	Serial number
Quick Response Code	Contains serial number
	With this mark, we confirm that the product complies with the requirements set out in the EC Directives and has been subjected to the specified testing procedures.
	UKCA: United Kingdom Conformity Assessed With this mark, we confirm that the product complies with the requirements set out in the UK Designated Standards.
www.brand.de/ip	Patent information
 (here: 40 years)	China RoHS (EFUP) EFUP defines the period in years during which the hazardous substances contained in electrical and electronic equipment do not leak or change under normal operating conditions. Under normal use by the user, such electrical and electronic products do not cause severe environmental pollution, serious personal injury or damage to the user's property.
	KC mark Korea (on battery pack)

Symbol or number	Meaning
	This electrical instrument may not be disposed of in household waste.

17 Ordering Information

Color of the upper housing section	accu-jet® S with universal power adapter and country adapter Item no.	accu-jet® S without universal power adapter and without country adapter Item no.
Anthracite	26350	26360
Berry	26351	26361
Petrol	26352	26362
Amethyst	26353	26363

Scope of delivery, see . Note the item numbers, with or without power adapter.

Housing top cover

Description	Packaged unit	Item no.
Housing top cover, anthracite	1 pc.	26780
Housing top cover, amethyst	1 pc.	26781
Housing top cover, berry	1 pc.	26782
Housing top cover, petrol	1 pc.	26783

Replacement membrane filter

Description	Packaged unit	Item no.
0.2 µm, sterile (red color code)	1 pc in blister packaging	26530
0.2 µm	10 pcs in PE bag	26535
3 µm	1 pc in blister packaging	26052
3 µm	10 pcs in PE bag	26056

Power adapter

Technical specifications for power adapter, see Technical data, p. 70 .
Item no. [26606](#)

Battery pack

Technical specifications for battery, see Technical data, p. 70 . Item no. [26630](#)

Silicone adapter

Item no. [26508](#)

Adapter housing

Color: White. Item no. [26658](#)

Wall mount

Color: White. Item no. [26539](#)

Snap tape

3 sets. Item no. [26547](#)

18 Repairs

18.1 Sending for repair

NOTICE Transporting hazardous materials without approval is prohibited by law.

Clean the instrument thoroughly and decontaminate!

- When returning products, please enclose a general description of the type of malfunction and the media used. If information regarding media used is missing, the instrument cannot be repaired.
- Shipment is at the risk and the cost of the sender.

Outside USA and Canada

Complete the “Declaration on Absence of Health Hazards” and send the instrument to the manufacturer or supplier. Ask your supplier or manufacturer for the form. The form can also be downloaded from www.brand.de.

Within USA and Canada

Please clarify the requirements for the return delivery with BrandTech Scientific, Inc **before** sending the instrument in for service.

Return only cleaned and decontaminated instruments to the address provided with the Return Authorization Number. Place the Return Authorization number so that it is clearly visible on the outside of the package.

Contact addresses

Germany:

BRAND GMBH + CO KG
Otto-Schott-Straße 25
97877 Wertheim (Germany)
T +49 9342 808 0

USA and Canada:

BrandTech® Scientific, Inc.
11 Bokum Road
Essex, CT 06426-1506 (USA)
T +1-860-767 2562

F +49 9342 808 98000
info@brand.de
www.brand.de

India:

BRAND Scientific Equipment Pvt. Ltd.
303, 3rd Floor, 'C' Wing, Delphi
Hiranandani Business Park,
Powai
Mumbai-400 076 (India)
T +91 22 42957790
F +91 22 42957791
info@brand.co.in
www.brand.co.in

F +1-860-767 2563
info@brandtech.com
www.brandtech.com

China:

BRAND (Shanghai) Trading Co., Ltd.
Rm 201-202, North Tower,
No. 199 Kaibin Rd, Xuhui District, Shanghai
Shanghai 200030 (P.R. China)
T +86 21 6422 2318
F +86 21 6422 2268
info@brand.com.cn
www.brand.cn.com

19 Warranty

We shall not be liable for the consequences of improper handling, use, servicing, operating or unauthorized repairs of the device or for the consequences of normal wear and tear, especially of wearing parts such as pistons, seals, valves and the breakage of glass. The same applies for failure to follow the instructions of the operating manual. We are not liable for damage resulting from disassembly beyond that described in the operating manual or if non-original spare parts or components have been installed.

USA and Canada:

Find more warranty information on www.brandtech.com.

20 Disposal

The symbol  means that at the end of their service life, batteries/accumulators and electronic devices must be disposed of separately from household waste (unsorted municipal waste).

Electronic devices must be disposed of in accordance with Directive 2012/19/ EU of the European Parliament and of the Council from July 04, 2012 on waste from electrical and electronic equipment and in compliance with national disposal regulations.

Both batteries and accumulators (rechargeable batteries) contain materials that can be damaging to the environment and human health. Therefore, they must be properly disposed of in accordance with Regulation (EU) 2023/1542 of the European Parliament and of the Council from July 12, 2023 on batteries and waste batteries and in compliance with national disposal regulations. Only dispose of fully discharged batteries and accumulators.

⚠WARNING! Fire hazard due to heating after a short circuit

Do not short-circuit the battery to discharge it! Tape the plug with adhesive tape to prevent a short circuit in the collection container! Never disassemble the battery!

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