


CHEMISTRY DIAPHRAGM PUMPS WITH ATEX CONFORMITY

MD 4C EX VARIO

MV 10C EX VARIO

MD 4C EX VARIO + AK + EK

MV 10C EX VARIO + AK + EK

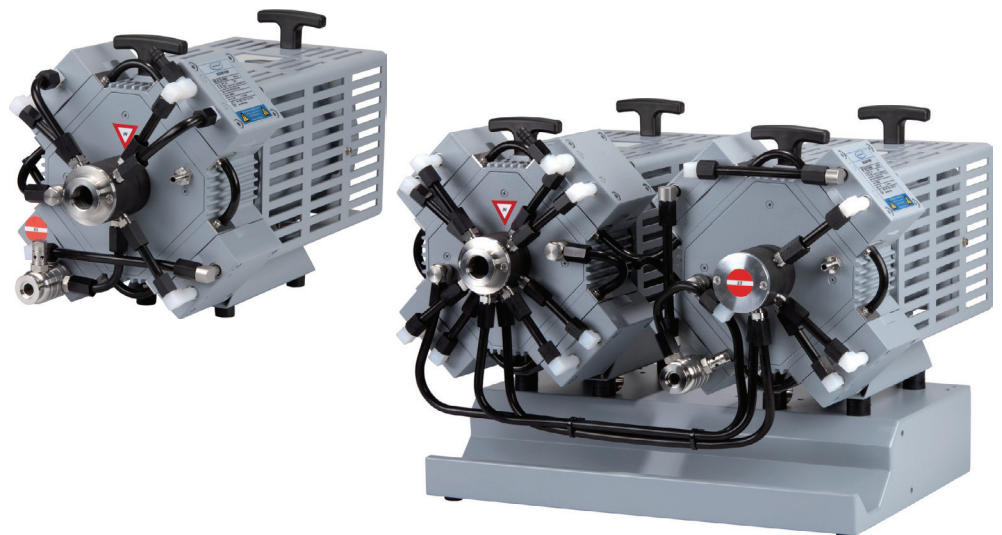

Instructions for use



Dear customer,

Your VACUUBRAND diaphragm pump shall support you at your work for a long time without any trouble and with full load output. Thanks to our large practical experience we attained much information how you could add to an efficient application and to personal safety. Please read these instructions for use prior to the initial start-up of your pump.

VACUUBRAND diaphragm pumps are the result of many years of experience in construction and practical operation of these pumps combined with the latest results in material and manufacturing technology.

Our quality maxim is the "zero fault principle":

Every delivered diaphragm pump is tested extensively including an endurance run of 14 hours. Due to this endurance run, also faults, which occur rarely, are reported and can be corrected. Every single diaphragm pump is tested on achievement of the specification after the endurance run.

Every VACUUBRAND pump leaving our factory achieves the specification. We feel obliged to this high quality standard.

We are aware that the vacuum pump should not draw a part of the real work and we hope to contribute with our products to an effective and trouble-free realisation of your work.

Yours

VACUUBRAND GMBH + CO KG

After sales service: Contact your local dealer or call +49 9342 808-5500.



Attention: The manual of the offered device with ATEX approval is available in German, English and French (DVR 3: also Dutch). The user is only permitted to operate the device if he understands one of these languages or a technically correct translation of the complete manual. The manual must be completely read and understood before operation of the device. All specified measures must be applied, or else must be replaced by equivalent measures at the user's own risk.

The device has an ATEX approval according to EC-directives and standards. Users outside of the European Community must check national regulations for hazardous areas.

Trademark index:

VACUU·LAN® (US-Reg.No 3,704,401), VACUU·BUS®, VACUU·CONTROL®, VACUU·SELECT® (US-Reg.No 5,522,260), VARIO® (US-Reg.No 3,833,788), VACUUBRAND® (US-Reg.No 3,733,388), VACUU·VIEW®, GREEN VAC® (US-Reg.No. 4,924,553), VACUU·PURE® (US-Reg No. 5,559,614) and also the shown company logos are registered trademarks of VACUUBRAND GMBH + CO KG in Germany and/or other countries.

D, A, CH, B, L

Achtung: Die Betriebsanleitung des angebotenen Geräts mit ATEX-Zulassung ist in Deutsch, Englisch und Französisch (DVR 3: auch Niederländisch) verfügbar. Der Anwender darf das Gerät nur dann in Betrieb nehmen, wenn er eine dieser Sprachen oder eine fachlich korrekte Übersetzung der vollständigen Anleitung versteht. Die Betriebsanleitung muss vor Inbetriebnahme des Geräts vollständig gelesen und verstanden werden, und alle geforderten Maßnahmen müssen eingehalten werden. Das Gerät verfügt über eine ATEX-Zulassung gemäß EU-Richtlinien und Normen. Anwender außerhalb der EU müssen die nationalen EX-Vorschriften prüfen.

GB, IRL

Attention: The manual of the offered device with ATEX approval is available in German, English and French (DVR 3: also Dutch). The user is only permitted to operate the device if he understands one of these languages or a technically correct translation of the complete manual. The manual must be completely read and understood before operation of the device and all required measures must be applied. The device has an ATEX approval according to EC-directives and standards. Users outside of the European Community must check national regulations for hazardous areas.

F, B, L

Attention : le manuel d'utilisation de l'appareil ATEX proposé est disponible en allemand, anglais et français (en néerlandais également pour le DVR 3). L'utilisateur est autorisé à utiliser l'appareil uniquement s'il comprend l'une de ces langues ou une traduction techniquement correcte du manuel complet. Le manuel doit être lu et compris entièrement avant l'utilisation de l'appareil et toutes les mesures requises doivent être appliquées. L'appareil a une certification ATEX suivant les directives CE et les standards. Les utilisateurs en dehors de la communauté européenne doivent vérifier les normes locales pour les zones dangereuses.

BG

Внимание: Ръководството за работа с предлаганото устройство, което е с ATEX одобрение е достъпно на следните езици: немски, английски и френски език (DVR 3: също холандски). Потребителят има право да работи с устройството само ако разбира много добре един от тези езици или ползва технически правилен превод на пълното ръководство. Ръководството трябва да бъде изцяло прочетено и разбрано преди да се пристъпи към работа с устройството, а всички необходими мерки трябва да бъдат спазвани. Устройството има ATEX одобрение съгласно ЕС - директивите и стандартите. Потребители извън Европейската общност трябва да се съобразят със съответните национални разпоредби за опасни зони.

CN

注意：所提供的ATEX批准设备的操作说明书有德文版、英文版和法文版（DVR 3也有荷兰语版）。客户仅允许懂得其中一种语言或者整本说明书的正确技术翻译的情况下，才可以操作该设备。该说明书在操作前必须完全阅读且能够理解，所有需要的测试必须采用。该装置根据EC指令和标准有ATEX批准。欧盟以外区域的客户必须核查各自国家针对危险区域的国家标准。

CZ



Pozor: Příručka pro nabízené zařízení s osvědčením ATEX je k dispozici v němčině, angličtině a francouzštině (DVR 3: i v holandštině). Uživatel smí dané zařízení používat pouze, pokud zná jeden z těchto jazyků nebo má k dispozici technicky správný překlad kompletního manuálu. Příručku musí kompletně přečíst a porozumět jí před zahájením provozu zařízení a musí zajistit uplatnění všech požadovaných opatření. Přístroj má schválení ATEX podle směrnic a norem EU. Uživatelé mimo území Evropského společenství, musí ověřit vnitrostátní předpisy pro prostředí s nebezpečím výbuchu.

DK



Bemærk: Betjeningsvejledningen til enheden "med ATEX-godkendelse" er til rådighed på Tysk, Engelsk og Fransk (DVR 3: også på Hollandsk)
Brugeren har kun tilladelse til at betjene enheden hvis, han/hun forstår et af de disse sprog eller har en teknisk korrekt oversættelse af hele manualen.
Betjeningsvejledningen skal læses og forstås før enheden kan anvendes og betjenes. Alle de krævede foranstaltninger skal overholdes.
Enheden har en ATEX godkendt certificering ifølge EF- direktiver og standarder.
Brugere uden for Det Europæiske Fællesskab skal kontrollere nationale regler for anvendelse.

EE



Tähelepanu: käsiraamat pakutud ATEX-i poolt heakskiidetud seadmele on saadaval saksa, inglise ja prantsuse keeles (DVR 3: puhul ka Hollandi keeles). Kasutajal on lubatud seda seadet kasutada ainult juhul, kui ta saab aru ühest nendest keeltest või on olemas tehniliselt korrektne tõlge kogu juhendist. Juhend peab olema täielikult loetud ja mõistetud enne seadmega töötamist ning kõiki vajalike meetmeid tuleb rakendada. Seadmel on ATEX-i heakskiit vastavalt EÜ-direktiividele ja standarditele. Kasutajad väljaspool Euroopa Liitu peavad kontrollima riiklikuid regulatsioone ohtlikke piirkondade kohta.

ES



Atención: Las instrucciones de funcionamiento del equipo ofrecido con homologación ATEX están disponibles en Alemán, Inglés y Francés (DVR 3: también en holandés). El usuario deberá poner en funcionamiento el equipo únicamente si comprende uno de estos idiomas o una correcta traducción técnica del manual completo en el idioma que comprenda. Las instrucciones deben ser leídas y entendidas antes de utilizar el equipo y deben cumplirse todas las medidas necesarias. El equipo posee una aprobación de conformidad ATEX con las directivas y normas de la UE. Los usuarios fuera de la UE deben verificar las normativas nacionales de protección contra explosiones.

FI



Huomio! Tuotteen käyttöohje, joka sisältää ATEX -todistuksen, on saatavilla saksaksi, englanniksi ja ranskaksi (DVR 3 myös hollanniksi). Laitetta saa käyttää vain henkilö, joka ymmärtää jotain em. kielistä tai hänellä on teknisesti tarkka käänнос alkuperäisestä käyttöohjeesta. Käyttäjän tulee lukea ja ymmärtää käyttöohje kokonaisuudessaan sekä noudattaa siinä mainittuja toimenpiteitä ennen laitteen käyttöä. Tuotteella on ATEX hyväksyntä EU-direktiivien ja standardien mukaisesti. Jos laitetta käytetään EU:n ulkopuolella, tulee noudattaa paikallisia määräyksiä.

GR

Προσοχή: Το εγχειρίδιο του προσφερόμενου οργάνου με πιστοποίηση ATEX είναι διαθέσιμο στα Γερμανικά, Αγγλικά, Γαλλικά (DVR3: επίσης Ολλανδικά). Ο χρήστης μπορεί να λειτουργεί το μηχάνημα εφόσον καταλαβαίνει μία από τις συγκεκριμένες γλώσσες ή μία τεχνικά σωστή μετάφραση του πλήρους εγχειριδίου. Το εγχειρίδιο θα πρέπει να αναγνωσθεί πλήρως και να γίνει κατανοητό πριν την λειτουργία του μηχανήματος και θα πρέπει να εφαρμόζονται όλα τα απαιτούμενα μέτρα. Το όργανο έχει πιστοποίηση ATEX σύμφωνα με τις οδηγίες και τα στάνταρ της Ευρωπαϊκής Ένωσης (Ε.Ε.). Χρήστες εκτός Ε.Ε. πρέπει να ελέγξουν τους εθνικούς τους κανονισμούς για τις τοποθεσίες επικινδυνότητας.

HR

Pažnja: Upute o rukovanju ponuđenih uređaja sa ATEX odobrenjem su dostupne na njemačkom, engleskom i francuskom jeziku (DVR 3: također i na holandskom). Korisniku je dopušteno rukovanje uređajem ako razumije jedan od ovih jezika ili ako ima prijevod tehničke dokumentacije. Uputa za rukovanje prije korištenja uređaja mora se pročitati, razumjeti te je potrebno primijeniti sve nužne mjere za ispravno rukovanje. Uređaji imaju ATEX odobrenje prema smjernicama Europske unije i važećim standardima. Korisnici izvan Europske unije moraju provjeriti nacionalnu regulativu vezanu uz rad sa potencijalno opasnim supstancama te uređajima koji se koriste u tom potencijalno opasnom području rada.

HU

Figyelem: a megajánlott ATEX minősítéssel rendelkező készülék kézikönyve németül, angolul és franciául érhető el (a DVR 3 hollandul is). Kérjük, csak akkor használja a berendezést, ha tökéletesen ért ezeken a nyelveken, vagy rendelkezésére áll egy technikailag megfelelő fordítás a teljes kézikönyvről. A kézikönyvet a készülék használatba vétele előtt kérjük végigolvasni, és betartani a benne foglaltakat. A készülék az EC-irányelveknek és előírásoknak megfelelő ATEX minősítéssel rendelkezik. Az Európai Közösségen kívüli vásárlók ellenőrizték az adott országban érvényes veszélyes helyiségekre vonatkozó előírásokat.

IT

Attenzione: il manuale della strumentazione con certificazione ATEX offerta è disponibile in Tedesco, Inglese e Francese. (DVR 3: anche in Olandese). L'utilizzatore può utilizzare l'apparecchiatura solo se capisce una di queste tre lingue o una traduzione tecnicamente corretta dell'intero manuale. Prima di utilizzare l'apparecchiatura il manuale deve essere completamente letto e tutte le indicazioni devono essere seguite. La strumentazione ha un'approvazione ATEX in conformità agli standards ed alle direttive EC. Al di fuori della Comunità Europea gli utilizzatori devono controllare la regolamentazione nazionale per le aree pericolose.

JP

注意：ATEX認証された製品に付きましては、ドイツ語、英語、フランス語（DVR 3はオランダ語も）の取扱説明書がございます。ユーザーはこれらの言語が理解できるか、技術的に正しい完全なマニュアルがある場合のみ、装置を作動することが許されます。製品使用前にマニュアルは全て読み、理解しなければなりません。また、全ての要求される対策を講じなければなりません。この製品はECの指令及び規格により承認されています。EC外のユーザーに付きましては、自国の危険場所に関する規定をご確認ください。

KR

주의 : ATEX 인증을 받은 본 장비의 매뉴얼은 독어, 영어, 불어 (DVR3 는 네덜란드어도 포함) 로 되어있습니다. 이들 언어중 하나를 독해 가능하거나 혹은 이 매뉴얼을 기술적으로 완벽하게 이해한 사용자만이 이 장비를 구동해야 합니다. 장비를 구동하기 전에 매뉴얼을 모두 읽고 이해하여야 하며, 모든 필수 요건들을 지켜야합니다. 장비는 ATEX 인증을 받았으며 이 인증은 EC의 지침과 규격에 따른 것입니다. 유럽공동체 이외 지역의 사용자들은 반드시 방폭에 대한 국가규정을 확인해야 합니다.

LT

Dėmesio: Pasiūlyto prietaiso vartotojo instrukcija yra patvirtinta ATEX ir galima vokiečių, anglų bei prancūzų kalbomis (DVR3: taip pat olandų kalba). Teisę dirbti su prietaisu turi tik vartotojas, suprantantis bent vieną iš minėtų užsienio kalbų arba turintis techniškai taisytiną vartotojo instrukciją. Prieš pradėdant dirbti su prietaisu privaloma perskaityti ir suprasti vartotojo instrukciją bei laikytis visų nurodytų reikalavimų. Prietaisas turi ATEX patvirtinimą remiantis EC-direktyvomis ir standartais. Vartotojai, nepriklausantys Europos Sąjungai, turi pasitikrinti jų šalyje galiojančius teisės aktus/reglamentus dėl pavojingų zonų.

LV

Uzmanību: Piedāvātās iekārtas ar ATEX apstiprinājumu lietošanas instrukcija ir pieejama vācu, angļu un franču valodā (DVR 3: arī holandiešu valodā). Lietotājam ir atļauts darboties ar šo iekārtu tikai tajā gadījumā, ja viņš saprot vismaz vienu no augstāk minētām valodām vai ja ir nodrošināts pilns un tehniski pareizs lietošanas instrukcijas tulkojums. Pirms iekārtas darbināšanas lietošanas instrukciju ir jāizlasa un jāizprot, kā arī jāpieņem visas nepieciešamās darbības. Iekārta ir ATEX apstiprināta, atbilstoši EK direktīvām un standartiem. Ārpus Eiropas Kopienas valstu lietotājiem ir jāpārbauda valsts nacionālos likumus par bīstamām zonām.

NL

Attentie: de handleiding voor het geoffreerde apparaat met ATEX goedkeuring is beschikbaar in het Duits, Engels en Frans (en voor de DVR 3 ook in het Nederlands). Het is de gebruiker alleen toegestaan om het apparaat te bedienen indien hij/zij een van deze talen begrijpt of indien er een technisch correcte vertaling van de complete handleiding beschikbaar is. De handleiding dient in zijn geheel gelezen en begrepen te worden en alle nodige maatregelen dienen geregeld te zijn alvorens het apparaat bedient mag worden. Het apparaat heeft een ATEX goedkeuring conform de EG-richtlijn en standaarden. Gebruikers buiten the Europese Gemeenschap moeten de nationale richtlijnen voor gevaarlijk locaties controleren.

PL

Uwaga: Instrukcja oferowanego urządzenia z dopuszczeniem ATEX jest dostępny w języku niemieckim, angielskim i francuskim (DVR 3: również holenderskim). Użytkownik może korzystać z urządzenia tylko wtedy, gdy zapoznał się z instrukcją w jednym z podanych języków, lub poprawnym technicznym tłumaczeniem pełnej instrukcji. Przed użyciem instrukcja obsługi musi być przeczytana i zrozumiana, oraz wszystkie wymagane środki bezpieczeństwa muszą być przestrzegane. Urządzenie posiada certyfikat ATEX zgodnie z dyrektywami i normami. Użytkownicy spoza Wspólnoty Europejskiej powinni sprawdzić krajowe przepisy dla stref zagrożonych wybuchem.

PT

Atenção: O manual do dispositivo oferecido com aprovação ATEX está disponível em Alemão, Inglês e Francês (DVR 3: também em holandês). O utilizador só é autorizado a operar o dispositivo se dominar um destes idiomas ou se tiver disponível um manual de operação completo, técnica e corretamente traduzido. O manual deverá ser lido e entendido na sua totalidade antes de se iniciar o funcionamento do dispositivo e todas as medidas necessárias deverão ser aplicadas.

O dispositivo tem uma aprovação ATEX de acordo com as diretivas e normas da Comunidade Europeia. Os utilizadores fora da Comunidade Europeia deverão verificar os regulamentos nacionais relativos às áreas de risco.

RO

Atentie: Manualul aparatului este disponibil in urmatoarele limbi: germana, engleza si franceza (DVR 3: de asemenea si in olandeza).

Operatorul poate utiliza acest dispozitiv doar daca intelege una din aceste limbi sau exista o traducere tehnica corecta a manualului.

Manulul trebuie citit complet si inteles inainte de punerea in functiune a aparatului si toate cerintele solicitate trebuie aplicate.

Dispozitivul are aprobare ATEX conform cu directivele si standardele EC. Utilizatorii din afara comunitatii Europene trebuie sa verifice regulile specifice tarii lor si standardele pentru utilizarea aparatului intr-un mediu periculos.

RU

Внимание: инструкция на предложенный аппарат с сертификатом ATEX доступна на немецком, английском и французском (DVR 3: также на датском). Пользователю разрешено работать с прибором только, если он понимает один из этих языков или технически корректный полный перевод инструкции.

Инструкция должна быть полностью прочитана и понята до начала работы с прибором, должны быть выполнены необходимые требования для работы прибора.

У прибора есть сертификат ATEX согласно стандартам ЕС. Пользователи вне ЕС должны проверять требования своих стран для работы в опасных областях.

SE

Observera: Manualen om den erbjudna enheten med ATEX-godkännande finns på tyska, engelska och franska (DVR 3: även nederländska). Användaren är endast tillåten att använda enheten om han förstår något av dessa språk eller en tekniskt korrekt översättning av den fullständiga manualen. Manualen måste vara helt läsas och förstås innan drift av enheten och alla nödvändiga åtgärder måste tillämpas. Enheten har en ATEX-godkännande enligt EG-direktiv och standarder. Användare utanför Europeiska gemenskapen måste kontrollera nationella regler för farliga områden.

SI

Opomba: navodila za uporabo za ponujeno napravo z odobritvijo ATEX je na voljo v nemškem, angleškem in francoskem jeziku (DVR 3: tudi nizozemščina). Uporaba naprave je dovoljena samo uporabnikom kateri so popolnoma prebrali in razumejo navodila v enem izmed ponujenih jezikov ali tehnično pravilen prevod celotnega priročnika.

Naprava ima ATEX odobritev v skladu z EC direktivami in standardi. Uporabniki zunaj Evropske skupnosti morajo preveriti nacionalne predpise za nevarna območja.

SK



Upozornenie: Návod pre ponúknutý prístroj s ATEX schválením je dostupný v nemeckom, anglickom alebo francúzskom jazyku (DVR 3: tiež v holandskom). Užívateľ môže používať prístroj iba v prípade ak rozumie jednému z týchto jazykov alebo technicky správne prekladu celého návodu na obsluhu. Je nutné si prečítať celý návod na obsluhu a pochopiť ho pred začatím práce s prístrojom a musí pritom dodržať všetky požadované opatrenia.

Prístroj má ATEX schválenie podľa požiadaviek kladených smernicami a normami Európskej únie. Užívatelia mimo Európskej únie si musia overiť platné národné predpisy kladené na priestory s nebezpečenstvom výbuchu.

TR



Dikkat : ATEX onayı ile sunulan cihazın kuulanma kitabı Almanca, İngilizce ve Fransızca dillerinde mevcuttur (DVR3: ek olarak Hollanda dilinde). Kullanıcı bu dillerden birini tam olarak anladığı yada teknik olarak bütün kullanma kitabını kendi diline çevirebildiği durumda cihazı kullanabilir. Cihazın kullanımasından önce kullanma kitabının tamamı okunmuş olmalı ve gerekli ölçümler yapılmış olmalıdır.

Cihaz EC-direktifleri ve standartlarına uygun ATEX onayına sahiptir. Avrupa Topluluğu dışında bulunan kullanıcılar, tehlikeli alanlar için kendi ülkelerinin düzenlemelerini kontrol etmelidirler.

Contents

Safety information!	12
General information	13
Intended use	13
Setting up and installing the equipment	13
Ambient conditions	17
Operating conditions	17
Safety during operation	19
Maintenance and repair	20
Notes regarding devices with ATEX conformity	21
Technical data	23
Use and operation	29
Installing in a vacuum system	29
During operation	34
Shutdown	34
Example for the build-up of a vacuum system with Connection Box 100	35
Accessories with Connection Box 100	36
Example for the build-up of a vacuum system with single components	37
Accessories for single components build-up	38
Circuit diagram for MD 4C EX VARIO	40
Circuit diagram for MV 10C EX VARIO	40
Safety analysis concerning the intrinsic safety of the vacuum sensor (pressure transducer) supply	41
Troubleshooting	42
Replacing diaphragms and valves	43
Inspecting the overpressure safety valve	52
Repair - Maintenance - Return - Calibration	55
EC Type examination certificate	57



- ➔ DANGER indicates a hazardous situation which, if not avoided, will result in death or serious injury.



- ⚠ WARNING indicates a hazardous situation which, if not avoided, could result in death or serious injury.



- CAUTION indicates a hazardous situation which, if not avoided, could result in minor or moderate injury.



NOTICE is used to address practices not related to personal injury.



Caution! Hot surface!



Isolate equipment from mains before removing the cover.



Electronic components must not be disposed of in the domestic waste at the end of their service life. Used electronic devices contain harmful substances that can cause damage to the environment or human health. End users are legally obliged to take used electric and electronic devices to a licensed collection point.

Safety information!



Attention: To operate the pump / pumping unit additional devices are required, see section “Example for the build-up of a vacuum system”.

General information

NOTICE

☞ **Read and comply with this manual before installing or operating the equipment.**

☞ Transport the pump at the provided handle.

Remove all packing material, remove the product from its packing-box, remove the protective covers from the inlet and outlet ports and keep, inspect the equipment.

If the equipment is damaged, notify the supplier and the carrier in writing within three days; state the item number of the product together with the order number and the supplier's invoice number. Retain all packing material for inspection.

Do not use the equipment if it is damaged.

If the equipment is not used immediately, replace the protective covers. Store the equipment in suitable conditions.

Intended use

! WARNING

☞ Use the equipment for the intended use only according to the ATEX conformity, e. g. for pumping gas atmosphere without aerosol in a range from 0.1 to 1000 mbar and a maximum inlet pressure (dynamic pressure) and outlet pressure of 1100 mbar (absolute).

☞ The pump and all system parts must not be used on humans or animals.

☞ Prevent any part of the human body from coming into contact with vacuum.

☞ Ensure that the individual components are only connected, combined and operated according to their design and as indicated in the instructions for use.

☞ Comply with notes on correct vacuum and electrical connections, see section "Use and operation".

! CAUTION

The pumps are marked with "X" (according to DIN EN ISO 80079-36:2016), i. e. restrictions of the operation conditions:

- **The pumps are designated for a low degree of mechanical stress and have to be installed in a way so that they can not be damaged from outside. Pumping units with glass attachment parts have to be installed protected against shocks from the outside and against glass splinters in the event of breaking (implosion). Attention: The joint clips at the glass attachment parts are not grounded. Avoid charge.**
- **The pumps are designated for an ambient and gas inlet temperature during operation of +10 to +40°C. Never exceed these ambient and gas inlet temperatures.**

NOTICE

Setting up and installing the equipment

Before system start-up of the pump or the complete system the user has to evaluate in its entirety the explosion risks and takes adequate measures in application of the principle of risk assessment.

The evaluation and the measures have to be laid down in an explosion protection document before system start-up.

The pump/system must not be operated until operation without danger in explosive atmosphere is ensured. **Provide suitable control, protection and warning systems even in the event of failure.**

The inner (pumped media) and outer part of the pump have to be evaluated separately because of their different conformity (see also section "Technical data" and "Notes regarding devices with ATEX conformity").

The pumps are **not suitable** for pumping dust and have no approval for operation below ground.

The pumps are **not suitable** for pumping self inflammable substances, for substances which are inflammable without air and explosive substances.

WARNING

- ☞ **Electrical and vacuum components might require a separate approval.**
- ☞ Include the components in the recurrent tests of tightness for pressure vessels, in the case of hits and beats an intermediate leak test is required.
- ☞ Due to the high compression ratio of the pumps, pressure at the outlet port might be generated being higher than the max. permitted pressure compatible with the mechanical stability of the system.
- ☞ Do not permit any **uncontrolled pressurizing** (e.g. make sure that the exhaust pipeline cannot become blocked). **Risk of bursting!**
- ☞ Always provide a free and pressureless exhaust pipeline.
- ☞ If there is an exhaust isolation valve, make sure that you cannot operate the equipment with the valve closed.

CAUTION

- Secure coolant hoses at the hose nozzles (e.g. with hose clip) to prevent their accidental slipping.
- Comply with **maximum permissible pressures** at inlet and outlet and pressure differences, see section "Technical data". Do not operate the pump with overpressure at the inlet.
- **Attention:** Flexible elements tend to shrink when evacuated.

NOTICE

Provide a firm level platform for the equipment and check that the system to be evacuated is mechanically stable and that all fittings are secure. Ensure a stable position of the pump without any mechanical contact except of the pump feet. Comply with all applicable **safety regulations**.

Pumping units with glass attachment parts have to be installed protected against hits from outside and protected against splinters (against implosion). Adopt suitable measures prior to the first use.

Attention: The joint clips at the glass attachment parts are not grounded. Avoid charge.

Obey all relevant requirements (regulations and guidelines) and adopt **suitable safety measures**.

WARNING

- ☞ **Pay attention to the max. permitted ambient temperature and make sure ventilation is adequate especially if pump is installed in a housing or if ambient temperature is elevated, install an external automatic ventilation system if necessary. If pumping hot process gases make sure that the maximum permitted gas inlet temperature is not exceeded.**
- ☞ If the equipment is brought from cold environment into a room for operation, allow the equipment to warm up (pay attention to water condensation on cold surfaces).
- ☞ The work clothing of the user must not lead to electrostatic discharge which may ignite the explosive atmosphere.

! WARNING

Inlet

- ☞ Connection lines at the pump inlet have to be performed **conductive grounded** and gas tight (designation of the pump connections see figures).
- ☞ The diameter of the inlet pipeline should be at the least as large as the diameter of the pump connection pipelines.
- ☞ Obey max. inlet pressure (see "Technical data").

! WARNING

Outlet

- ☞ Connection lines at the pump outlet (small flange) have to be performed **conductive grounded** and gas tight (designation of the pump connections see figures).
- ☞ The diameter of the inlet pipeline should be at the least as large as the diameter of the pump connection pipelines.
- ☞ Obey max. inlet pressure (see "Technical data").
- ☞ Lead away exhaust gas or condensate at the outlet of the pump and the overpressure safety valve under consideration of all applicable safety measures. **Avoid backflow of gases or condensate definitely.**
- ☞ Escaped and/or intentional or unintentional **released inflammable gases** or vapours which may lead to danger of explosion have to be lead away safely or lead to an appropriate place. If not possible enclose the gasses, vapours or mist safely or dispose in an other way. In case of mixtures or different gases the measures must consider the highest risk.

Check the **overpressure safety valve** at the outlet regularly.

! WARNING

- ☞ **Attention: If the exhaust pipeline is blocked the pumped gases may escape into the environment through the overpressure safety valve. If necessary take appropriate measures.**

If pumping dangerous gases install a second completely independent exhaust pipeline instead of the overpressure safety valve. If necessary assemble a hose nozzle instead of the overpressure safety valve to lead away the gases through a second exhaust pipeline. **Attention:** Never combine the exhaust pipelines, i. e. provide a sperate line for each exhaust.

Inlet and outlet of the pump

! WARNING

- ☞ To avoid damage of the pump due **ignition** of gases or gas mixtures before or behind the pump the user has to provide suitable **safety devices against incoming of flames** according to EN 12874. The user must ensure the suitability concerning gas flow, chemical resistance and safeness against clogging prior to use.

Inert gas purge

! WARNING

- ☞ The user must take suitable precautions to prevent any formation of explosive mixtures in the expansion chamber using inert gas to purge the housing. Otherwise in case of a diaphragm crack, mechanically generated sparks, hot surfaces or static electricity may ignite these mixtures.
- ☞ **Inert gas connection** lines (inlet and outlet) have to be performed conductive grounded and gas tight (designation of the pump connections see figures).
- ☞ Connect the **inert gas connections** at the crankcase not serial but parallel.
- ☞ We recommend an inert gas purge of ≥ 1 l per minute.
- ☞ Provide an external pressure limitation to a pressure of 1.1 bar absolute if **inert gas** is connected to the pump or at an air admittance valve. The inert gas has to be dry and pure. It could be recommendable to provide an appropriate control system to supervise the inert gas flow which switches off the pump in case of failure (e. g. absence of inert gas) if necessary.
- ☞ To control the faultless function of the diaphragm it could be recommendable install a **gas specific detector** at the **outlet of the inert gas purge**. A signal from the detector indicates a diaphragm crack. Switch off the pump immediately and check diaphragms.

- ☞ It is possible to leave out the inert gas purge if the surrounding of the pump is zone 2 or not an explosive atmosphere.
- ☞ Without inert gas purge the pump outer part has the specification II 3G Ex h IIB T4 Gc X.

WARNING

Gas ballast

- ☞ Use only inert gas for gas ballast.
- ☞ We recommend the use of gas ballast when pumping condensable media.
- ☞ Gas ballast connection lines (inlet and outlet) have to be performed conductive grounded and gas tight (designation of the pump connections see figures).
- ☞ Connect the **inert gas connections** at the gas ballast and the crankcase not serial but parallel.
- ☞ Provide an external pressure limitation to a pressure of 1.1 bar absolute if **inert gas** is connected to the pump or at an air admittance valve. The inert gas has to be dry and pure. It could be recommendable to provide an appropriate **control system** to supervise the inert gas flow which switches off the pump in case of failure (e. g. absence of inert gas) if necessary.

Pumping at **high inlet pressure** may lead to overpressure at the gas ballast valve.

WARNING

- ☞ Pumped gases or condensate might be pushed out in case the valve is open.
- ☞ The inert gas lines have to be chemical resistant due to possible backflow.

WARNING

Electrical connection

- ☞ **Electrical connection** of the pump must be performed only by a suitable trained and supervised personnel.
- ☞ The motor cable for pumps in 230 V version contains wires color coded as follows: green and yellow: earth; blue: neutral; brown: live.
- ☞ Provide a fuse for the mains cable (L and N) according to the current draw of the motor, see "Technical data". When using pumps with **two motors** provide a fuse for each pump motor separately. Due to a high charge current for each motor a protection is recommended, e. g., automatic fuse 10A type C.
- ☞ The **earth connection** of the device (potential equalization) has to be performed only using the green or green/yellow earth connection wire. Connect the earth connection wire of the device to the earth connection of the power supply line and the potential equalization point of the environment of the device.
- ☞ When using pumps with **two motors** absolutely connect both motors as described above. Ensure that always both motors are running.

Connection of the control cable

- ☞ **Electrical connection** of the pump must be performed only by a suitable trained and supervised personnel.
In case of pumps with two motors (MV 10C EX VARIO) connect the control cables in series.
Maximum voltage: 12 V.
Maximum current: 60 mA.
Comply with polarity!

The control cable of the pump contains four colored wires:

brown (+): 4 - 20 mA / speed control
blue (-): 4 - 20 mA / speed control
black: Status
grey: Status

Attention: Never use the antistatic connection at the housing of the device to connect the device to ground, use only the green or green and yellow connection wire of the motor. Use the antistatic connection at the housing only to perform an antistatic connection of other non-electronic devices and components, e. g. inlet and outlet vacuum hoses or something similar. Never lead away short-circuit currents using this connection. External electric devices have to be connected at the potential equalization point of the environment separately and independent of this device.

WARNING

- ☞ Equipment must be connected only to a suitable fused and protected electrical supply and a suitable earth point by suitable trained personnel. Failure to connect the motor to ground may result in deadly electrical shock.
- ☞ Check that mains voltage and current conform with the equipment (see rating plate).

The **motor** and the **motor circuit board** are protected by a thermal protection with **manual reset**.

Attention: Only manual reset is possible. Switch off the pump and disconnect from the power source. Identify and eliminate the cause of failure. Wait approx. five minutes before restarting the pump.

WARNING

- ☞ Avoid reliable high heat supply (e. g. due to hot process gases). Maximum permitted gas and ambient temperature see "Technical data".
- ☞ Make sure ventilation is adequate especially if pump is installed in a housing or if ambient temperature is elevated, install an external automatic ventilation system if necessary.
- ☞ Ensure that in case of **pumps with two motors** both motors are switched on at the same time, if not possible switch on the motor of the pump at the outlet first.

Ambient conditions

To the best of our knowledge the equipment is in compliance with the requirements of the applicable EC-directives and harmonized standards (see "Declaration of conformity") with regard to design, type and model, especially directive EN 61010-1. This directive gives in detail conditions, under which the equipment can be operated safely (see also IP degree of protection).

Adopt suitable measures in case of differences, e. g. using the equipment outdoors, installation in altitudes of more than 1000 m above mean sea level, conductive pollution or dewiness.

Pay attention to the **permissible maximum ambient and gas inlet temperatures** (see "Technical data").

Operating conditions

The pumps have a conformity **according to their specification** to be installed in areas and to pump out of areas which contain under normal operation conditions occasionally a potential explosive atmosphere as a mixture of air with inflammable gases or vapours (**category 2**). **Class of ignition and temperature see "Technical data"**.

The specification for category 2 for the outer part of the pump is only valid when purging the crankcase with inert gas. Without inert gas purging the pump has the specification II 3G Ex h IIB T4 Gc X for the outer part.

WARNING

- ☞ Adopt suitable measures to prevent the release of dangerous, explosive, corrosive or polluting fluids.
- ☞ If pumping **different substances**, it is recommended to purge the pump with air or inert gas prior to changing the pumped media in order to pump out residues and to avoid reactions of the pumped substances with each other and with the pump materials.
Take into consideration interactions and chemical reactions of the pumped media.
- ☞ Ensure that the materials of the wetted parts are compatible with the pumped substances, see section "Technical data".
- ☞ The **valves and the diaphragms** are wear parts. **Replace the diaphragms and valves at the latest at 90% of the typical lifetime or immediately at higher noise level.** The typical lifetime of a diaphragm is 15000 operation hours when pumping non-corrosive or non-condensable gases, if necessary provide an operation time counter.
- ☞ Check the **overpressure safety valve between the pump stages** (only pumps with four cylinders) regularly and replace **at the latest at 90% of the typical lifetime** (typical lifetime: 15000 operation hours).
- ☞ Take adequate precautions to protect people from the effects of dangerous substances (chemicals, thermal decomposition products of fluoroelastomers), wear appropriate safety-clothing and safety glasses.
- ☞ Obey applicable regulations when disposing of chemicals. Take into consideration that chemicals may be polluted.
- ☞ Adopt suitable measures to prevent contamination of the pumped substances or the environment.

Electronic equipment is never 100% fail-safe. This may lead to an indefinite status of the equipment. Provide protective measures against malfunction and failure.

WARNING

- ☞ Operating the pump, stand still of the pump or operating the air admittance valve must not lead to a critical dangerous situation under any circumstances.

Ensure that in case of failure the pump and the vacuum system always will turn into a safe status.

WARNING

- ☞ In case of leaks in the manifold or diaphragm cracks pumped substances might be released into the environment or into the pump housing. To reduce the risk of leaks, ask for a diaphragm pump with additional safety diaphragm.
- ☞ Obey especially notes on operation and use and maintenance.
- ☞ Failure of the pump (e. g. by power failure) or connected components, parts of the supply (e. g. coolant) or change of parameters (e. g. increase of pressure in the coolant system) must not lead to a critical dangerous situation under any circumstances.

Safety during operation



- ➔ In case install an appropriate collecting and disposal system and take protective action for pump and environment.
- ➔ Potentially explosive mixtures at the outlet of the pump have to be drained appropriately, sucked off or diluted with inert gas to non-explosive mixtures.
- ☞ Pay attention to the symbol "hot surfaces" on the equipment. Adopt suitable measures to prevent any danger arising from the formation of hot surfaces or electric sparks. Provide a suitable protection against contact if necessary.
- ☞ Check liquid level in both catchpots regularly and drain condensate in time.
- Comply with applicable regulations when disposing of chemicals. Take into consideration that chemicals may be polluted. Take adequate precautions to protect people from the effects of dangerous substances (chemicals, thermal decomposition products of fluoroelastomers), wear appropriate safety-clothing and safety glasses.



Use only genuine spare parts and accessories.

- ☞ Otherwise safety and performance of the equipment as well as the electromagnetic compatibility of the equipment might be reduced.
- ☞ **If not using genuine spare parts the ATEX conformity becomes invalid.**

Do not start the pump if the pressure difference between inlet and outlet port exceeds 1.1 bar at maximum.

Prevent the backpressure of gases and the backflow of condensates.
Never suck liquids or dust into the pump.

Ensure that in case of failure the pump and the vacuum system always will turn into a safe status.

Provide appropriate protective measures (i.e. precautions which allow for the requirements of the respective application) even for the case of failure and **malfunction**.

Comply especially with notes on operation and use and maintenance.

Due to the residual **leak rate of the equipment**, there might be an exchange of gas, albeit extremely slight, between the environment and the vacuum system. Adopt suitable measures to prevent contamination of the pumped substances or the environment.

Maintenance and repair

Wear parts have to be replaced regularly. In normal use, the lifetime of the diaphragms and valves is typically 15,000 operating hours. Bearings have a typical durability of 40000 h.

Use only genuine spare parts and accessories.

☞ Otherwise safety and performance of the equipment as well as the electromagnetic compatibility of the equipment might be reduced.

☞ **If not using genuine spare parts the ATEX conformity becomes invalid.**

Ensure that maintenance is done only by suitable trained and supervised technicians.

➔ **Isolate equipment from mains and wait two minutes** before starting maintenance to allow the capacitors to discharge.

- Before starting maintenance vent the pump, isolate the pump and other components from the vacuum system. Allow sufficient cooling of the pump. Drain condensate, if applicable.

☞ **Ensure that the pump cannot be operated accidentally. Never operate the pump if covers or other parts of the pump are disassembled. Never operate a defective or damaged pump.**

☞ **Attention:** The pump might be contaminated with process chemicals which have been pumped during operation. Ensure that the pump is decontaminated before maintenance and take adequate precautions to protect people from the effects of dangerous substances if contamination has occurred.

☞ Take adequate precautions to protect people from the effects of dangerous substances (chemicals, thermal decomposition products of fluoroelastomers), wear appropriate safety-clothing and safety glasses.

☞ To avoid danger due to electrostatic ignition wipe the device or parts of the device only with a humid cloth.

In order to comply with law (occupational, health and safety regulations, safety at work law and regulations for environmental protection) vacuum pumps, components and measuring instruments returned to the manufacturer can be repaired only when certain procedures (see section "**Repair - Maintenance - Return - Calibration**") are followed.



Notes regarding devices with ATEX conformity

This information sheet does not replace the instructions for use!

In directive 89/391/EEC the European Union has defined requirements for safety at work, which later have been detailed for explosive atmospheres in the directive 1999/92/EC.

The requirements to equipment for operation in explosive atmospheres have been defined in the directive 94/9/EC, since April 20th 2016 in directive 2014/34/EU (also known as ATEX 100a or ATEX 95).

Concerning technology these directives only state general requirements. Technical details and test procedures are "recommended" in European standards (e.g. for equipment with ATEX conformity: EN 1127, E50014ff, EN 60079, DIN EN ISO 80079-36:2016).

Since July 1st 2003 equipment intended for use in potentially explosive atmospheres shall only be sold if the equipment is in accordance with directive 94/9/EG (ATEX), since April 20th 2016 in accordance with directive 2014/34/EU. This applies only to new equipment. The repair of equipment sold prior to July 1st, 2003 is not affected. EX approvals according to directive 94/9/EC remain valid under 2014/34/EU.

Requirements to users and manufacturers of equipment according to ATEX:

The **user of a system according to ATEX** has to analyse the overall explosion risks according to directive 1999/92/EC and has to take adequate measures by means of a risk assessment. The evaluation and validation of measures has to be documented in an explosion protection document, specifying among other things the areas classified as potentially explosive (see below). The inner part of the pump (pumped gases) and the environment of the pump (outer explosion protection) have to be evaluated separately. Often the requirements will be different.

The responsibility for the classification of areas with potentially explosive atmospheres into zones and for the specification of requirements to the equipment (category of equipment, class of ignition, temperature class etc.) is accountable strictly to the operator of the system. The equipment used has to comply with these requirements. VACUUBRAND does not advise customers concerning requirements to the equipment under the specific aspects of the customer's applications.

Equipment with ATEX-conformity is to be used by appropriately instructed personnel only.

ATEX marking, intended use:

The marking of the equipment corresponding with the standards (see above) gives the areas in which the equipment may be used according to its intended use.

**Technical data, safety notes and intended use:
See instructions for use!**

The equipment must not be used if the instructions for use are not read, fully understood and obeyed.

Clean, check and maintain equipment regularly.

Do never operate a defective equipment.

Electrical connection:

VACUUBRAND diaphragm pumps with ATEX-conformity (device cat. 2) are equipped with a 230 V / 50 Hz single-phase motor and are protected by a thermal cutout with manual reset combined with an overcurrent protection.

See instructions for use concerning notes on electrical connection.

Where applicable the motors can be connected directly to an appropriate power supply by suitably trained and authorised personnel (electrician). It is the responsibility of the user to meet the requirements of explosion protection in the power supply and connection area.

For operation of VACUUBRAND ATEX VARIO pumps a motor speed control signal (4 – 20 mA) is required.

Vacuum and outlet connections:

**See instructions for use: Make sure that all mechanical connections
are always gas tight and electrically conducting.**

The exhaust pipeline must be designed in a way that it cannot become blocked.

If the exhaust pipeline becomes blocked, the overpressure safety device at the pump releases potentially dangerous gases into the environment. Take appropriate safety measures (see instructions for use).

Evaluate the risk of explosion for the release of pumped gases at the outlet of the pump.

Use and connection of inert gas supply:

For external ATEX equipment category 2G (e.g. for use in Ex-zone 1) the crank case of the pump has to be flushed with inert gas (e.g., nitrogen). A flow of approx. 1 l/min (at atmospheric pressure up to max. 0.1 bar over pressure) is sufficient. If necessary the flow should be monitored using a flow meter. Without inert gas purge the pump has external equipment category 3G. This means the pump has to be used in this case only in a surrounding with Ex - zone 2 (or no Ex-zone).

It is recommended to install a gas detector (specific to the pumped gas) at the outlet of the inert gas tubing. The maximum permitted overpressure at the inert gas outlet is 0.1 bar (check the flow rate through the gas detector!). A positive signal from the detector indicates a diaphragm failure. In this event switch off the pump immediately and check the pump.

**In the case of pumping condensable vapours, use gas ballast. Connect inert gas with
maximum overpressure of 0.1 bar to the gas ballast inlet (see instructions for use).**

**Inlet of air into the gas ballast must never lead to formation of an explosive mixture inside the
pump
or at the outlet of the pump!**

Technical data

Type		MD 4C EX VARIO MD 4C EX VARIO + AK + EK	MV 10C EX VARIO MV 10C EX VARIO + AK + EK
ATEX approval Inner part (pumped gases) Outer part Surrounding of the pump with inert gas purge Surrounding of the pump without inert gas purge Motor		II 2G Ex h IIC T3 Gb X II 2G Ex h IIB T4 Gb X II 3G Ex h IIB T4 Gc X II 2G Ex db IIB T4 Gb	
Max. pumping speed* according to ISO 21360	m ³ /h	3.7	8.1
Ultimate vacuum* (absolute) without gas ballast	mbar	3	2
Ultimate vacuum* (absolute) with gas ballast	mbar	10	10
Max. permitted pressure at the inlet and the outlet (absolute)	bar	1.1	
Max. permitted pressure difference between inlet and outlet	bar	1.1	
Max. permitted pressure at inert gas connection (absolute)	bar	1.1	
Max. permitted ambient temperature during operation	°C	+10 to +40	
Max. permitted ambient temperature during storage	°C	-10 to +60	
Permitted gas inlet temperature	°C	+10 to +40	
Max. surface temperature of the pump (outer part)	°C	110	
Max. permitted atmospheric moisture during operation (no condensation)	%	30 to 85	
Degree of protection of the pump IEC 529 Pump Pump + AK + EK		IP 54** IP 54**	
Nominal current draw	A	3.2	2 x 3.2
Maximum charge current / duration	A / ms	150 / 2.5	2 x 150 / 2.5
Max. permitted range of voltage supply / frequency		230V ±10%	
Motor speed min./max.	min ⁻¹	30/1500	
Motor power	kW	0.25	2 x 0.25
Max. voltage control cable	V	12	
Max. current control cable	mA	60	
Impedance control cable (per motor) max. / typical	Ω	200 / 170	
Max. integral leak rate	mbar·l/s	0.1	

* Technical data according to EN 61010-1 and EN 1012-2 recommendation. The pump achieves its ultimate pumping speed and ultimate vacuum only at operating temperature (after approx. 15 min.).

** Only when positioning the pump horizontal.

We reserve the right for technical modifications without prior notice!

Documents are only to be used and distributed completely and unchanged. It is strictly the users' responsibility to check carefully the validity of this document with respect to his product.

Type		MD 4C EX VARIO MD 4C EX VARIO + AK + EK	MV 10C EX VARIO MV 10C EX VARIO + AK + EK
A-weighted emission sound pressure level*** (uncertainty K_{pA} : 3 dB(A))	dB(A)	42	42
Inlet Pump Pump + AK + EK		KF 25 KF 25	
Outlet Pump Pump + AK + EK		KF 16 hose nozzle 10 mm	
Gas ballast connection		hose nozzle 10 mm	
Inert gas purge connection		hose nozzle 8 mm	
Recommended flow of inert gas	l/min	≥ 1	
Coolant connection at exhaust waste vapor condenser		hose nozzle 6-8 mm	
Overall dimensions L x B x H, approx. Pump Pump + AK + EK	mm mm	469 x 265 x 305 600 x 365 x 420	560 x 457 x 410 651 x 555 x 452
Mass, approx. Pump Pump + AK + EK	kg kg	28 36	61 63

*** Measurement according to EN ISO 2151:2004 and EN ISO 3744:1995 at 230V/50Hz and ultimate vacuum with exhaust tube at outlet.

We reserve the right for technical modifications without prior notice!

Wetted parts

Components	Wetted parts
Inlet	stainless steel*
Outlet Pump Pump + AK + EK	stainless steel** PBT
Hoses	PTFE, antistatic
Fittings	ETFE / stainless steel*
Housing cover insert	PTFE, carbon reinforced
Head cover	ETFE, carbon reinforced
O-ring in head cover	FPM
Diaphragm clamping disc	ETFE, carbon reinforced
Valve	FFKM or PTFE
Diaphragm	PTFE
Overpressure safety relief device	stainless steel***, PTFE, carbon reinforced, FFKM
Separator cover plate	PTFE, carbon reinforced
Catchpot, collecting flask, vapor condenser	borsilicate glass

* stainless steel 1.4404

** stainless steel 1.4541

*** stainless steel 1.4404 and 1.4310

Pump parts


Position	Component	Position	Component
1	inlet (vacuum connection)	10	exhaust waste vapor condenser
2	outlet (gas!)	11	coolant inlet (hose nozzle 6-8 mm)
3	rating plate	12	coolant outlet (hose nozzle 6-8 mm)
4	handle	13	catchpot
5	gas ballast connection (only inert gas)	14	collecting flask
6	inlet inert gas purge	15	mains cable
7	outlet inert gas purge	16	control cable
8	overpressure valve		
9	antistatic connection		

We reserve the right for technical modifications without prior notice!


MD 4C EX VARIO



MV 10C EX VARIO

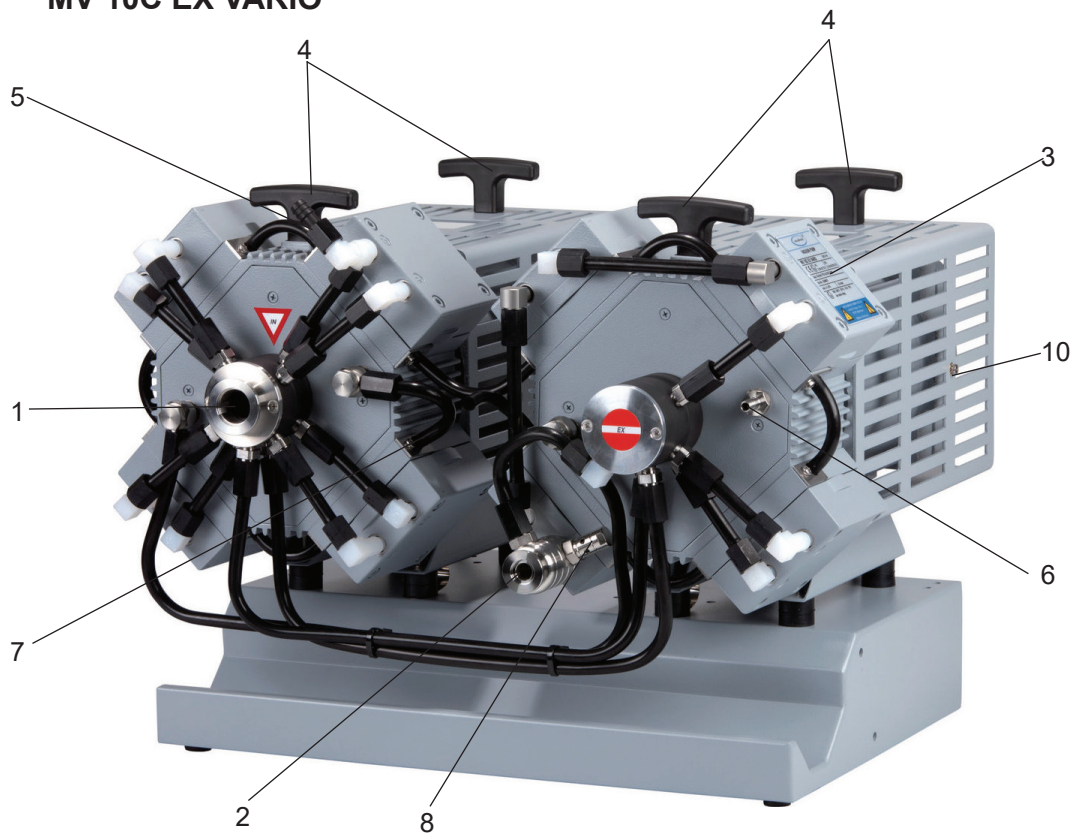


MD 4C EX VARIO + AK + EK

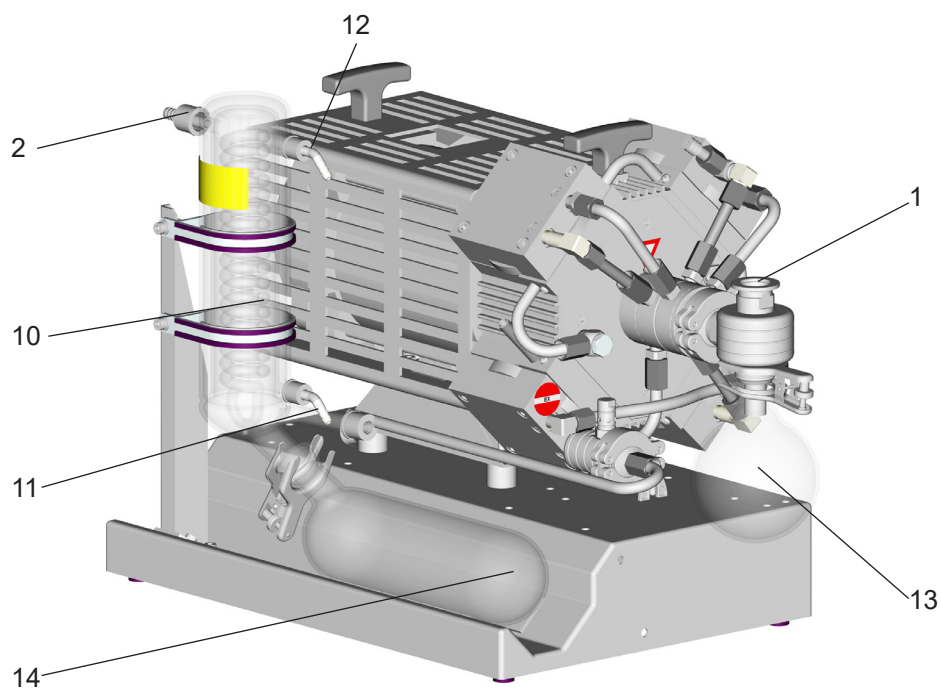


Attention: Further designation of the pump components see figure MD 4C EX VARIO

MV 10C EX VARIO + AK + EK

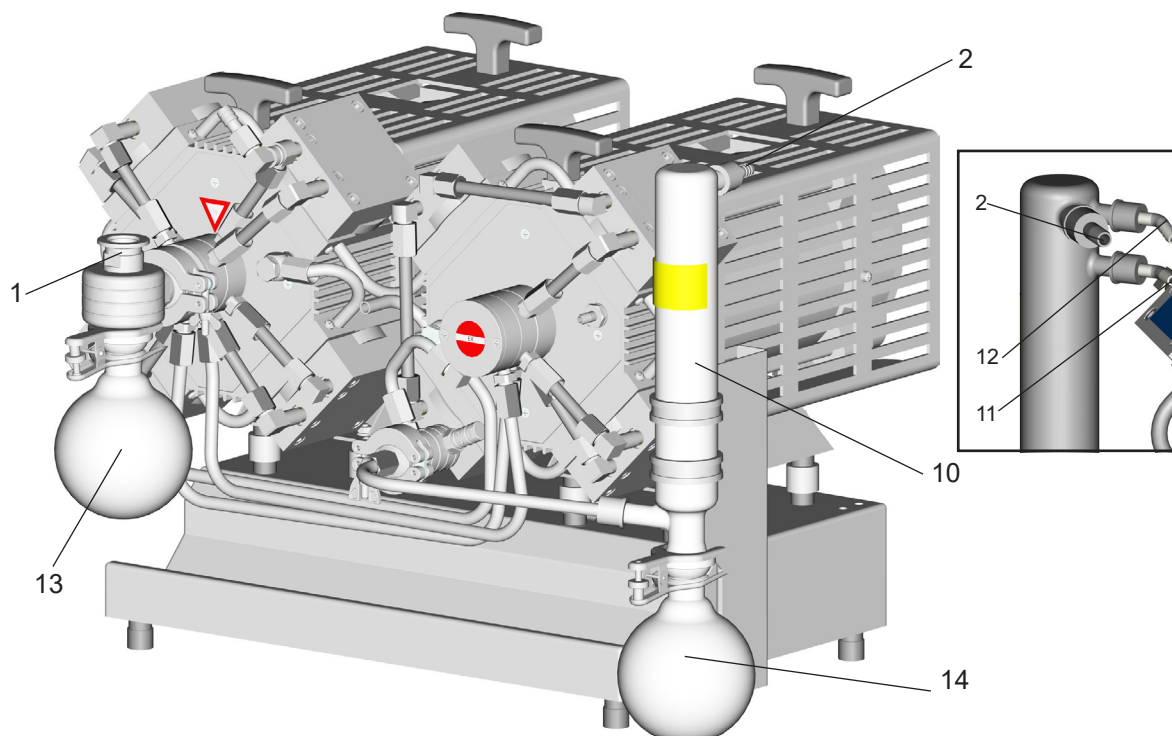

Achtung: Weitere Bezeichnungen der Pumpenteile siehe Abb. MV 10C EX VARIO

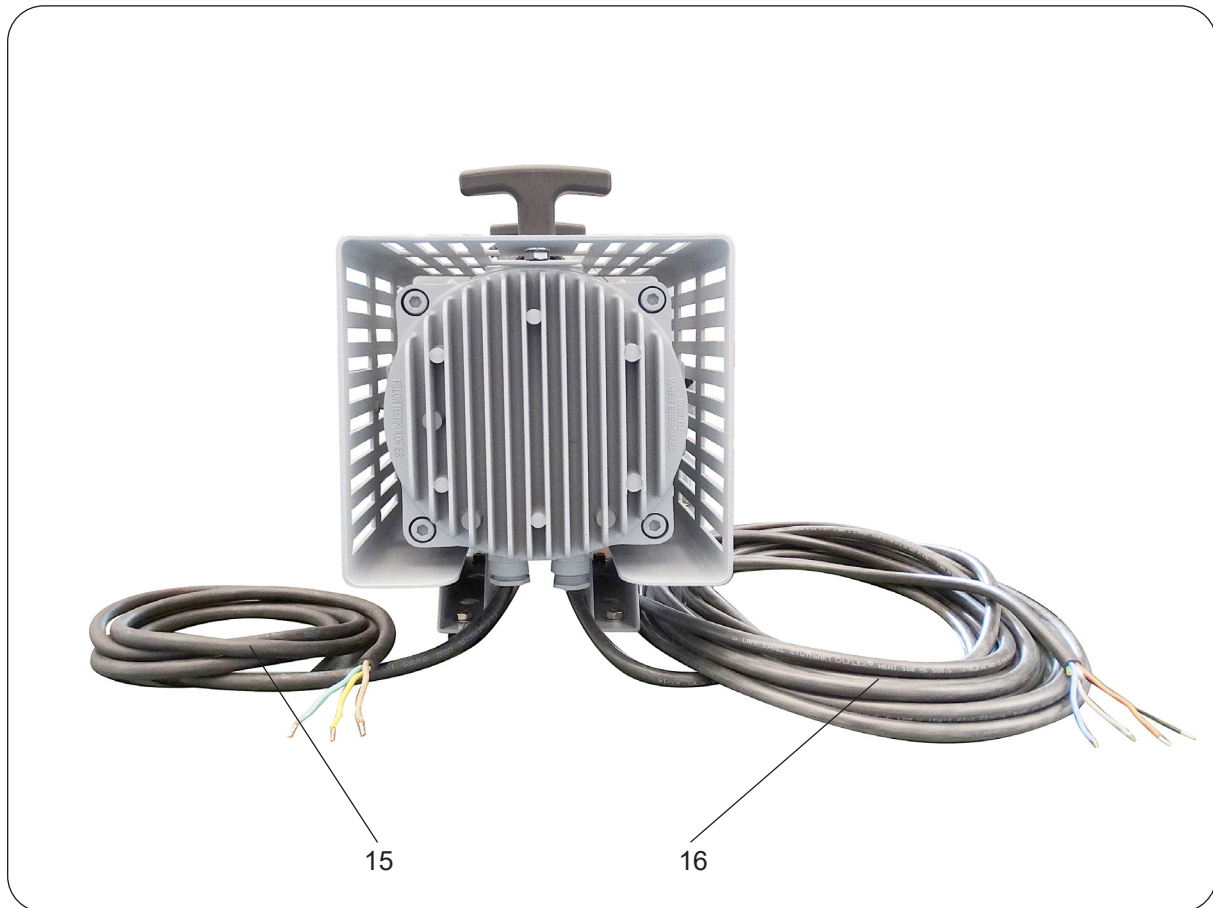

Mains cable

green and yellow: earth
blue: neutral (N)
brown: live (L)

Control cable

brown (+): 4 - 20 mA / speed control
blue (-): 4 - 20 mA / speed control
black: status
grey: status

Use and operation

Installing in a vacuum system

CAUTION

- All connection lines at the pump (e. g. inlet, outlet, gas ballast connection, inert gas connection) have to be performed **conductive grounded**.
- Connect pump and all connected components to ground.
- Avoid contact of the pump with **rusty metal parts**, e. g. by positioning the pump with sufficient distance to surrounding parts.
- The pump is designated for a low degree of mechanical stress. Protect the pump from damage and beats. In case of damage switch off the pump immediately, check leak rate if necessary.
- Connection lines at the pump inlet have to be gas tight. Particles and dust must not be aspirated, the user has to provide appropriate filters if necessary. The user must ensure their suitability concerning gas flow, chemical resistance and safety against clogging prior to use.
- Connect an exhaust line gas tight at the pump outlet if necessary. Always dispose of exhaust gases appropriately (e.g. into a fume hood). If there is risk of release of dangerous or polluting fluids, install an appropriate system to catch and dispose of those fluids.
- Reduce the transmission of vibration and prevent mechanical load due to rigid pipelines. Insert elastic hoses or flexible elements as couplings between the pump and rigid pipes. **Attention:** Flexible elements tend to shrink when evacuated.
- Use of a suitable valve to isolate the pump from the vacuum system is recommended to allow the pump to warm up before pumping condensable vapors or to clean the pump before it is switched off.
- The gas outlet must never be blocked. The exhaust line has always to be free (pressureless) to ensure an unimpeded discharge of gas.
- Especially if the gas ballast valve is open, a power failure may cause accidental ventilation of the pump. In case this constitutes a potential source of danger, take appropriate safety measures.
- Make sure ventilation is adequate especially if the pump is installed in a housing or if the ambient temperature is elevated. Provide external ventilation if necessary. Keep a distance of minimum 20 cm between fan and ambient parts.

NOTICE

Avoid throttling losses by using connecting pipes with large diameter and by keeping them as short as possible.

In case of perturbing exhaust noise connect an exhaust hose.

Install outlet pipelines always falling to avoid backflow of condensate towards the pump.

When assembling, ensure **vacuum-tightness**. After assembly, check the whole system for leaks.

Secure hose connections at the pump appropriately against accidental detaching.

NOTICE

Attach the pipelines of the coolant circuit to the respective hose nozzles (hose nozzles 6-8 mm, see image) at the vapor condenser. Check hose connections prior to starting operation of the cooling system.

Secure coolant hoses at the hose nozzles (e.g. with hose clip) to prevent their accidental slipping.

⚠ CAUTION

- **Attention:** Install hoses of the cooling system in a way to avoid flow / dropping of condensed water onto the pumping unit (especially cables and electronic parts).

⚠ WARNING

Connection of the mains cable

- ☞ Electrical connection of the pump must be performed only by a suitable trained and supervised personnel.
When using pumps with two motors (MV 10C EX VARIO) connect mains cable parallel.
The motor cable for pumps in 230 V version contains wires color coded as follows:
green and yellow: earth
blue: neutral (N)
brown: live (L)
- ☞ Provide a fuse for the mains cable (L and N) according to the current draw of the motor, see "Technical data". When using pumps with **two motors** provide a fuse for each pump motor separately. Due to a high charge current for each motor a protection is recommended, e. g., automatic fuse 10A type C.
- ☞ The **earth connection of the device** (potential equalization) has to be performed only using the green or green/yellow earth connection wire. Connect the earth connection wire of the device to the earth connection of the power supply line and the potential equalization point of the environment of the device.
Attention: Never use the antistatic connection at the housing of the device to connect the device to ground, use only the green or green and yellow connection wire of the motor. Use the antistatic connection at the housing only to perform an antistatic connection of other non-electronic devices and components, e. g. inlet and outlet vacuum hoses or something similar. Never lead away short-circuit currents using this connection. External electric devices have to be connected at the potential equalization point of the environment separately and independent of this device.
- ☞ In case of **power failure** unintentional venting of the system is possible. Adopt suitable measures if this might lead to a dangerous situation.

⚠ WARNING

Connection of the control cable

- ☞ **Electrical connection** of the pump must be performed only by a suitable trained and supervised personnel.
In case of pumps with two motors (MV 10C EX VARIO) connect the control cables in series, load per motor $\leq 200 \Omega$.
Maximum voltage: 12 V.
Maximum current: 60 mA.
Comply with polarity!

The **control cable** of the pump contains four colored wires:

brown (+): 4 - 20 mA / speed control
blue (-): 4 - 20 mA / speed control
0 - 4 mA: Motor stops
4 - 20 mA: Linear increase of the motor speed from 30 - 1500 rpm
20 mA: Maximum motor speed of 1500 rpm

black: Status of control electronics
grey: Status of control electronics
contact closed: ok
contact open: fault
Max. 250 V, 2 A

⚠ CAUTION

The **motor** and the **motor circuit board** are protected by a thermal protection with **manual reset**.

Attention: Only manual reset is possible. Switch off the pump and disconnect from the power source. Identify and eliminate the cause of failure. Wait approx. five minutes before restarting the pump.

WARNING

Inlet

- ☞ If there is a danger of **deposits** in the pump chamber (check inlet and outlet of the pump) control the pump chamber regularly and clean if necessary.

WARNING

Connection of the inert gas purge

- ☞ If the surrounding of the pump is not an explosive atmosphere it is possible to leave out the inert gas purge.

WARNING

Connection of the inert gas gas ballast

- ☞ When using air rather than inert gas, risk of significant damage to equipment and/or facilities, risk of personal injury or even loss of life exists due to the formation of hazardous and/or explosive mixtures if air and pumped media react inside or at the outlet of the pump.

WARNING

Outlet of the pump

- ☞ **Always connect an exhaust lines at the pump outlet (small flange or hose nozzle 10 mm).** Perform the outlet line **conductive grounded** and gas tight (designation of the pump connections see figures). Always connect the exhaust to a **suitable treatment plant** (e. g. hood).

WARNING

Overpressure safety valve at the outlet of the pump

The **overpressure safety valve** at the outlet prevents reliable high pressures in the system. Check overpressure safety valve regularly and replace if necessary.

DANGER

Attention: Deposits and condensate may lead to increased temperature and to excess of the max. permitted temperatures! Increased temperatures possibly may lead to ignition of flammable mixtures inside the pump.
Use inert gas gas ballast if necessary and check pump regularly for deposits.
Install a gas washing bottle, condenser, filter, separator, etc. at the inlet of the pump if necessary.

Pumps with two motors

WARNING

- ☞ Ensure that in case of pumps with two motors both motors are switched on at the same time, if not possible switch on the motor of the pump at the outlet first.
- ☞ Connect the **inert gas connections** at the crankcase not serial but parallel.



Attention: Glass parts (flasks and condensers) of pumping units (pumps with AK (IK) and EK) have no protection against splinters (against implosion) or against leakage in case of mechanical damage or hits from outside.

The user must ensure an appropriate protection against splinters / implosion or leakage prior to the first use.

Pumps with glass parts




- ☞ To avoid electrostatic charging, friction against glass parts in potentially explosive atmospheres is not permitted.

Adhesive labels on the pump




- ☞ To avoid electrostatic charging, the removal of rating plates and adhesive labels in potentially explosive atmospheres is not permitted.

MD 4C EX VARIO + AK + EK

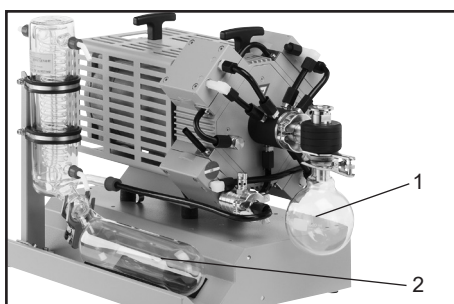


- ➔ Assemble **catchpot at the inlet (1)** using joint clips.
- ☞ The **collecting flask at the outlet (2)** is already assembled.

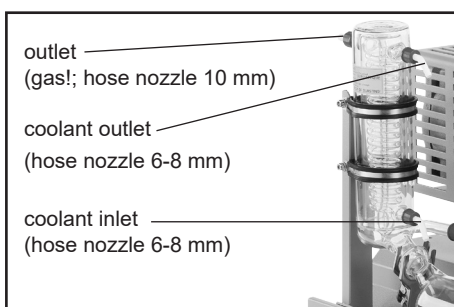


- ➔ Assemble **hose nozzles** for cooling water inlet and outlet at the vapor condenser.

MV 10C EX VARIO + AK + EK

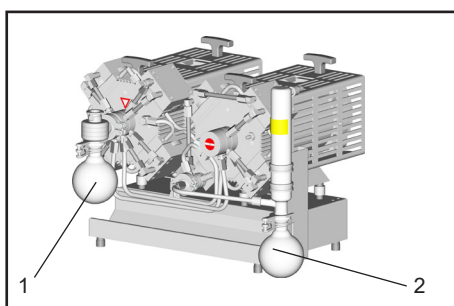

- ➔ Assemble **catchpot at the inlet (1) and the outlet (2)** using joint clips.

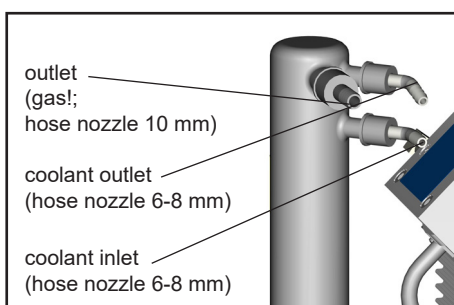

- ➔ Assemble **hose nozzles** for cooling water inlet and outlet at the vapor condenser.

Separation of condensate:

Attention: Due to electrostatics reasons the vapor condenser has no protective layer to avoid separation of condensed water! Risk of bursting!

The exhaust waste vapor condenser enables an efficient condensation of the pumped vapors at the outlet.

WARNING

- ☞ Next to 100% solvent recovery.
- ☞ Install hoses for cooling water at the inlet and outlet (hose nozzle 6-8 mm), secure hose connections against gliding down.

CAUTION

- Maximum permissible coolant pressure at the vapor condenser: 6 bar (absolute). Backflow always pressureless.
- Assemble an optional coolant valve only in the inlet line of the exhaust waste vapor condenser.
- Comply with the maximum permissible coolant pressures of additional components in the coolant circuit (e.g. coolant valve).
- Avoid overpressure in the coolant circuit (e.g. caused by blocked or squeezed coolant hoses).

Permissible range of coolant temperature at the vapor condenser:
-15°C to +20°C

Check hose connections prior to starting operation of the cooling system.
Check coolant hoses regularly during operation.

Do not allow the catchpots to get overfilled. Maximum liquid level approx. 80% to avoid problems when removing the catchpots.
Check liquid level in both catchpots regularly and drain catchpots in time.

Removing the catchpots:

Catchpot / collecting flask at outlet:

Remove joint clip, remove catchpot and drain condensate.

Catchpot at inlet:

Admit air or inert gas (via pump inlet) to atmospheric pressure. Remove joint clip, remove catchpot and drain condensate. Reassemble drained catchpots.

Important: Comply with regulations when disposing of solvents/condensates. Re-use if possible, purify if contaminated.

WARNING

For **condensable vapors** (water vapor, solvents, ...):

- ☞ Do not pump vapor before pump has reached its operating temperature and without inert gas ballast.
- ☞ Use inert gas at the gas ballast to avoid the formation of explosive mixtures.
- ☞ With gas ballast ultimate vacuum will be reduced
- ☞ Do not pump vapor before pump has reached its operating temperature

In case of low boiling solvents when the formation of condensate is unlikely, the use of gas ballast might be unnecessary.

WARNING

- ☞ Operating the pump without gas ballast increases the solvent recovery rates at the vapor condenser.

! WARNING

☞ The pumped gases at the outlet of the pump or at the exhaust waste vapor condenser have to be drained under consideration of all applicable safety regulations.

The user has to check the possibility of an explosive atmosphere at the outlet of the pump or the exhaust waste vapor condenser. Therefore for example calculate the solvent partial pressure at the present cooling water temperature at the exhaust waste vapor condenser and compare with the explosion limits of the pumped solvent. If there is an explosive mixture, drain the gas at the outlet in antistatic lines according to the valid explosion protection guidelines.

During operation

! CAUTION

- **Maximum ambient temperature:** 40 °C
- If the pump is installed in altitudes of more than 1000 m above mean sea level check compatibility with applicable safety requirements, especially IEC 60034 (motor might overheat due to insufficient cooling).
- **If the pump has been exposed to increased ambient temperature (> 40°C) check the pump for damage and perform a leak test if necessary (integral leak rate < 0.1 mbar*l/s).**

The pump achieves its pumping speed, ultimate total vacuum and vapor pumping rate only at operating temperature (after approx. 15 minutes).

! WARNING

☞ Prevent internal condensation, transfer of liquids or dust. The diaphragm and valves will be damaged, if liquids are pumped in significant amounts. Check the pump regularly at the outside for **pollution or deposits**, clean the pump if necessary to avoid increase of the operation temperature of the pump.

Shutdown

NOTICE

Short-term:

Has the pump been exposed to condensate?

Allow the pump to continue to run at atmospheric pressure for a few minutes.

Has the pump been exposed to media which may damage the pump materials or form **deposits**?

Check and clean pump heads if necessary.

Long-term:

Take measures as described in section short-term shutdown.

Separate pump from the apparatus.

Close inlet and outlet port (e. g. with transport caps).


Close the gas ballast valve.

Store the pump in dry condition


Example for the build-up of a vacuum system with Connection Box 100

Attention: Electrical connection of the pump must be performed only by electrically qualified person with additional qualification in explosion protection taking into account all relevant standards and regulations.

Build-up with MD 4C EX VARIO

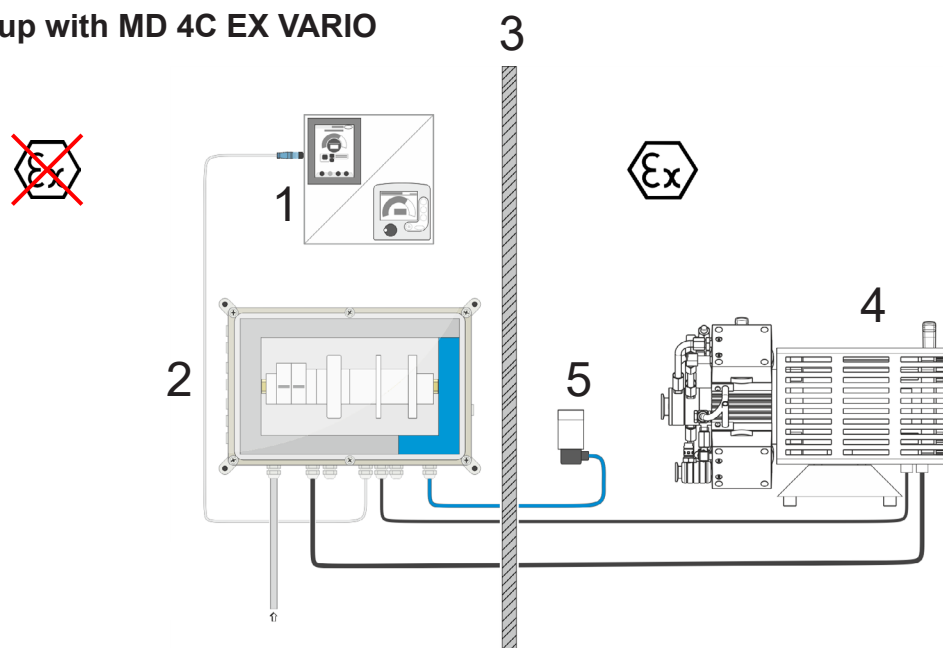

Build-up with MV 10C EX VARIO

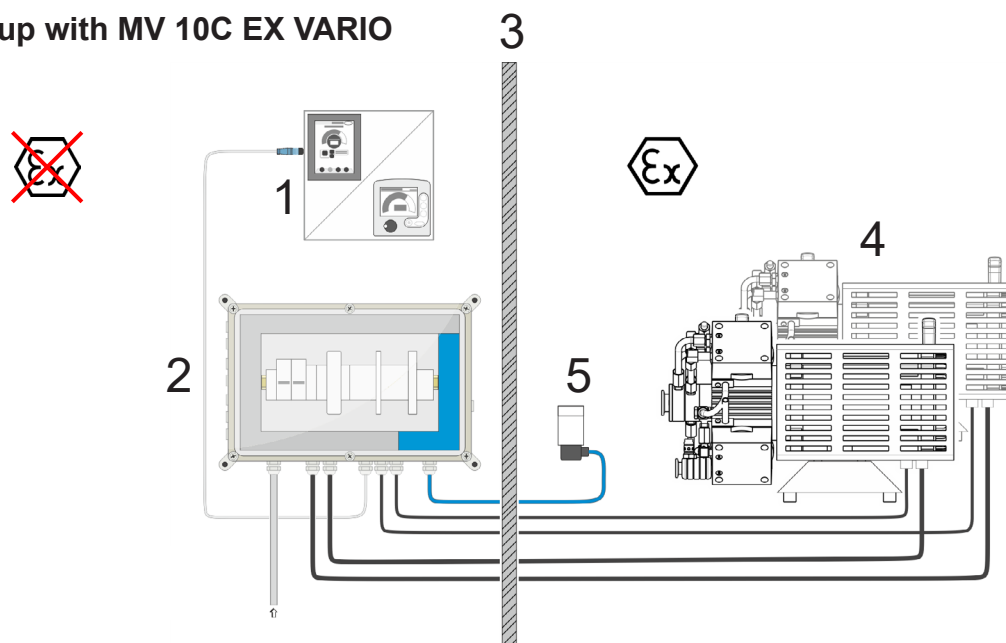

1	VACUUBRAND vacuum controller VACUU·SELECT / CVC 3000
2	Connection Box 100
3	Separating wall (spatial separation of explosive and non-explosive atmosphere)
4	VACUUBRAND VARIO diaphragm pump with ATEX approval
5	Vacuum sensor with ATEX approval

Accessories with Connection Box 100

Build-up of a system with Connection Box 100

Connection Box 100	20635435
Vacuum controller VACUU·SELECT operating panel..... without sensor, with wall plug power supply	20700050
alternatively: Vacuum controller CVC 3000	20636559
without sensor, without wall plug power supply	
Wall plug power supply VACUU·SELECT vacuum controller, VACUU·VIEW / extended, CVC / DCP 3000	20612089
Supply voltage 100-230 V~ / 50-60 Hz, Including interchangeable plugs: CEE, CH, UK, US, AUS	
ATEX vacuum sensor DMP 331i	20635423
BDSSENSORS, with KF DN 16, measuring range: 1000-1 mbar accuracy 1 mbar, with ATEX approval: II 1G Ex ia IIC T4 Ga, output signal 4 - 20 mA, for operation with EX i - power / input isolating amplifier 24V alternatively:	
ATEX vacuum sensor Cerabar PMP21	20635424
Endress+Hauser, with KF DN 16, measuring range: 1000-1 mbar, accuracy: 3 mbar, with ATEX approval: II 2G Ex ia IIC T4 Gb, output signal 4 - 20 mA, for operation with EX i - power / input isolating amplifier 24V	
Extension cable VACUU·BUS, 2 m	20612552

NOTICE

Several components are installed in the Connection Box 100 that are needed to evaluate the sensor signal. In this measurement chain, the measurement uncertainty of the vacuum value displayed on the vacuum controller increases to ± 3 mbar (for sensor 20635423 at 0.1 % FSO) and ± 4 mbar (for sensor 20635424 at 0.3 % FSO).


Further information concerning accessories on www.vacuubrand.com and on request.

PTFE tubing antistatic, KF DN 16/500 mm	20686030
PTFE tubing antistatic, KF DN 16/1000 mm	20686031
PTFE tubing antistatic, KF DN 25/500 mm	20686032
PTFE tubing antistatic, KF DN 25/1000 mm	20686033
Centring ring KF DN 20/25C..... with FEP coated seal ring (chemically resistant)	20635722

Example for the build-up of a vacuum system with single components

Attention: Electrical connection of the pump must be performed only by electrically qualified person with additional qualification in explosion protection taking into account all relevant standards and regulations.

For installation the single components have to be arranged by the customer in a suitable switch cabinet and wired correctly. For installation the customer himself is responsible.

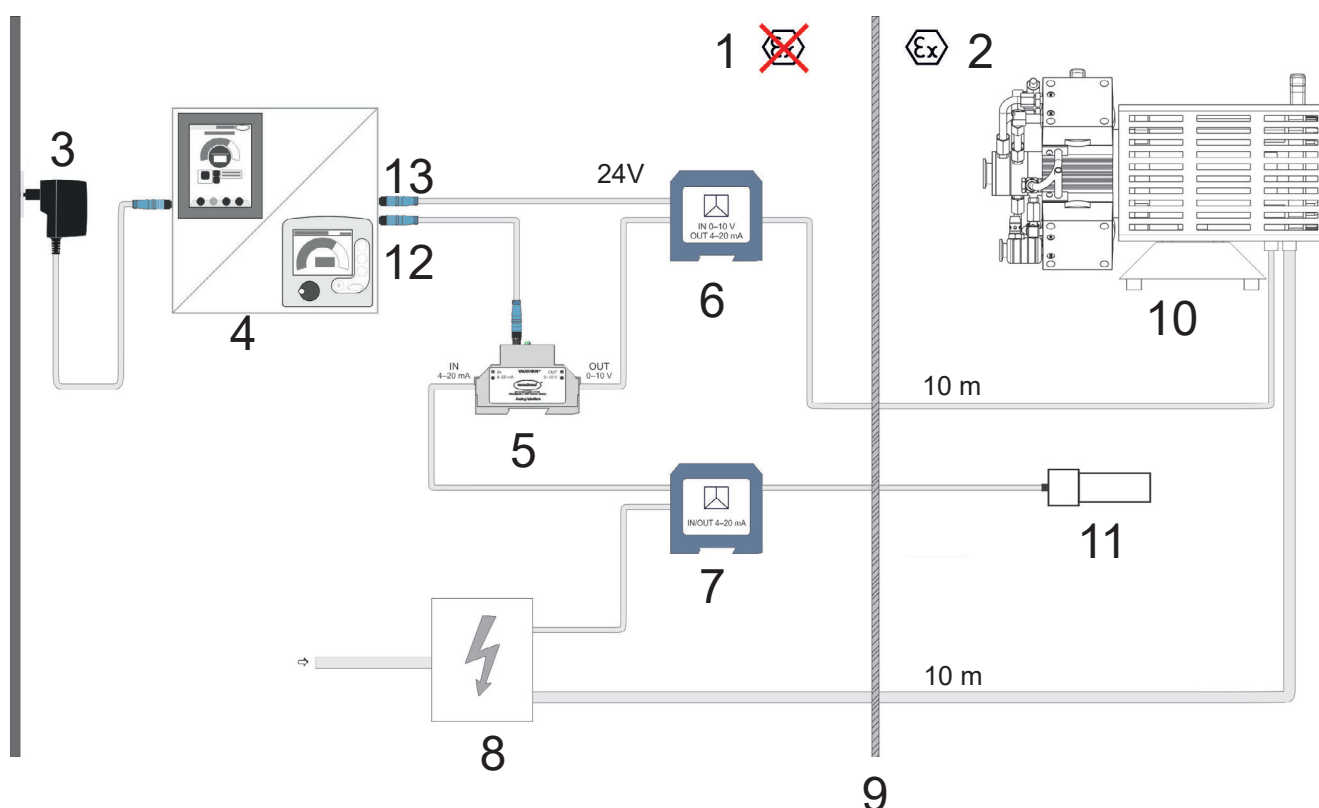

1	Zone without explosive atmosphere
2	Zone with explosive atmosphere
3	Power supply VACUUBRAND vacuum controller
4	VACUUBRAND vacuum controller VACUU·SELECT / CVC 3000
5	VACUUBRAND Analog I/O Module (4-20 mA / 0-10 V)
6	Isolation amplifier (signal converter)
7	Power/input isolating amplifier
8	Power supply pump
9	Separating wall (spatial separation of explosive and non-explosive atmosphere)
10	VACUUBRAND VARIO diaphragm pump with ATEX approval
11	Vacuum sensor with ATEX approval
12	Extension cable VACUU·BUS
13	VACUU·BUS cable with male plug and with litz wires

Accessories for single components build-up

Build-up of a system with single components

For installation the single components have to be arranged by the customer in a suitable switch cabinet and wired correctly.

Vakuum Controller VACUU·SELECT Bedieneinheit..... **20700050**
ohne Sensor, mit Steckernetzteil

alternatively:

Vacuum controller CVC 3000 **20636559**
without sensor, without wall plug power supply

Wall plug power supply VACUU·SELECT vacuum controller,
VACUU·VIEW / extended, CVC / DCP 3000 **20612089**
Supply voltage 100-230 V~ / 50-60 Hz,
Including interchangeable plugs: CEE, CH, UK, US, AUS

Analog I/O interface module VACUU·BUS **20635425**
converts an input signal 4 - 20 mA into a VACUU·BUS signal,
converts a VACUU·BUS signal into an output signal 0 - 10 V,
with 2 m VACUU·BUS cable

Isolating amplifier, for the connection of the ATEX VARIO pump/s **20635426**
for electrical isolation of analog signals, with screw connection, input signal: 0 - 10 V,
output signal: 4 - 20 mA, supply voltage 24V DC

Ex-i power/input isolating amplifier, for the connection of an ATEX vacuum sensor **20635427**
sends fed or active 0/4-20 mA signals from the Ex area to a load (active or passive)
to the safe area, electrical 3-way isolation, SIL 2, wide range power supply

ATEX vacuum sensor DMP 331i **20635423**
BDSSENSORS, with KF DN 16, measuring range: 1000-1 mbar
accuracy 1 mbar, with ATEX approval: II 1G Ex ia IIC T4 Ga, output signal 4 - 20 mA,
for operation with EX i - power / input isolating amplifier 24V
alternatively:

ATEX vacuum sensor Cerabar PMP21 **20635424**
Endress+Hauser, with KF DN 16, measuring range: 1000-1 mbar, accuracy: 3 mbar,
with ATEX approval: II 2G Ex ia IIC T4 Gb, output signal 4 - 20 mA,
for operation with EX i - power / input isolating amplifier 24V

Extension cable VACUU·BUS, 2 m **20612552**

VACUU·BUS cable 2 m, with male plug and with litz wires **20612462**


NOTICE

Several components are installed in this structure that are needed to evaluate the sensor signal. In this measurement chain, the measurement uncertainty of the vacuum value displayed on the vacuum controller increases to ± 3 mbar (for sensor 20635423 at 0.1 % FSO) and ± 4 mbar (for sensor 20635424 at 0.3 % FSO).

Further information concerning accessories on www.vacuubrand.com and on request.



PTFE tubing antistatic, KF DN 16/500 mm	20686030
PTFE tubing antistatic, KF DN 16/1000 mm	20686031
PTFE tubing antistatic, KF DN 25/500 mm	20686032
PTFE tubing antistatic, KF DN 25/1000 mm	20686033
Centring ring KF DN 20/25C.....	20635722
with FEP coated seal ring (chemically resistant)	

Circuit diagram for MD 4C EX VARIO

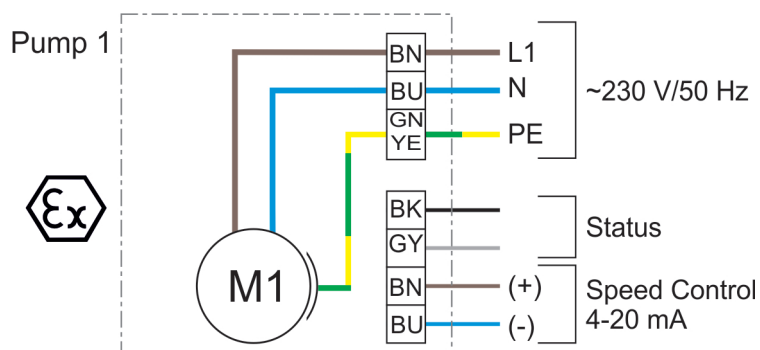

BU	blue
BN	brown
GN YE	green-yellow
GY	grey
BK	black

Circuit diagram for MV 10C EX VARIO

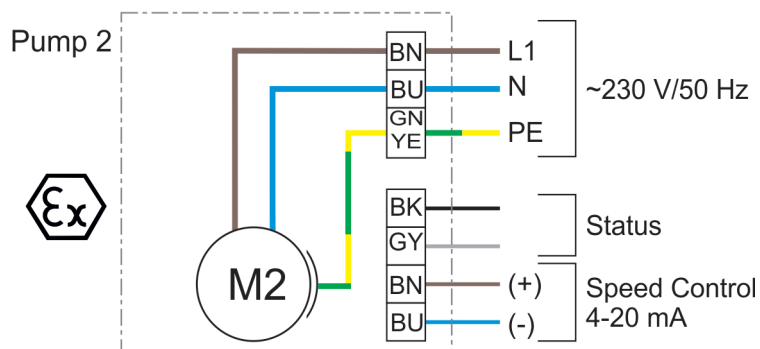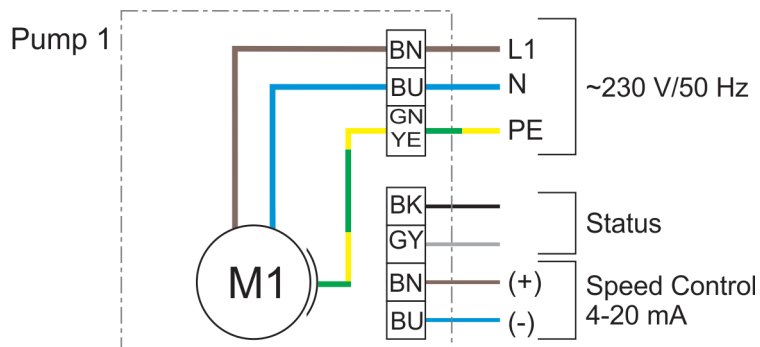

BU	blue
BN	brown
GN YE	green-yellow
GY	grey
BK	black

Safety analysis concerning the intrinsic safety of the vacuum sensor (pressure transducer) supply

Vacuum sensor (pressure transducer)

Vacuum sensor DMP 331i

Manufacturer: BD-Sensors, order no.: 111-1001-E-1-100-200-2-111
VACUUBRAND order no.: 20635423

Vacuum sensor Cerabar PMP21

Manufacturer: Endress+Hauser, order no.: PMP21-BA1U2HBWBJ
VACUUBRAND order no.: 20635424

Intrinsically safe power and current supply

Ex-i power/input isolating amplifier 4-20 mA

Manufacturer: Phoenix Contact, order no.: MACX MCR-EX-SL_RPSSI-I-UP-2865793
VACUUBRAND order no.: 20635427

The vacuum sensors are, according to their instructions for use, designed for operation with an intrinsically safe current supply, whereby the measuring signal output is carried out via current 4 - 20 mA. The power supply is carried out via the Ex-i power/input isolating amplifier.

Safety-related technical data according to ATEX (manufacturer's data)

Specification	Vacuum sensor DMP 331i	Vacuum sensor Cerabar PMP21	Ex-i power/input isolating amplifier 4-20 mA MACX MCR-EX-SL_RPSSI-I-UP-2865793	Rating
Voltage	U _{max} = 28 V	U < 30 V	Max. 25.2 V	fulfilled
Current	I _{max} = 93 mA	I < 100 mA	Max. 93 mA	fulfilled
Power output	P _{max} = 660 mW	P < 800 mW	Max. 587 mW	fulfilled
Inductance	L = 0	L = 0	Max. 2 mH	fulfilled
Capacity	C _{max} = 27 nF	C = 11.6 nF	Max. 107 nF	fulfilled

For safe operation the maximum permitted values for voltage, current and power output of the vacuum sensors must not be exceeded. Additionally the maximum permitted values for inductance and capacity at the power/input isolating amplifier must be observed. These conditions are met.

Troubleshooting

Never operate a defective pump!

Read and obey the instructions for use!

Fault	Possible cause	Remedy
☐ Pump does not start or stops immediately.	➔ Mains not plugged in, electrical supply failure? ➔ Pressure in outlet pipeline too high? ➔ Motor / circuit board overloaded? ➔ Control signal missing or control signal ≤ 4 mA?	✓ Check power supply and fuse in the building installation. ✓ Remove blockade in line, open valve. ✓ Allow the device to cool down, identify cause of failure and eliminate. Then switch on the pump. In case of recurrence contact customer service. ✓ Connect control signal.
☐ Pump does not achieve its ultimate vacuum or usual pumping speed.	➔ Centring ring at small flange connection not correctly positioned or leak in the pipeline or vacuum system? ➔ Long, narrow line? ➔ Pump has been exposed to condensate? ➔ Deposits have been formed inside the pump? ➔ Diaphragms or valves damaged? ➔ Outgassing substances or vapour generated in the process?	✓ Check pump directly - connect vacuum gauge directly at pump inlet - then check connection, pipeline and vacuum system if necessary. ✓ Use lines with larger diameter, length as short as possible. ✓ Allow pump to run for some minutes with atmospheric pressure at the inlet. ✓ Clean and inspect the pump heads. ✓ Replace diaphragms and/or valves. ✓ Check process parameters.
☐ Pump too noisy. Attention: Switch off pump immediately and inspect!	➔ Loud exhaust noise? ➔ Diaphragm crack or diaphragm clamping disc loose? ➔ Motor bearing defective? ➔ Deposits have been formed inside the pump? ➔ Diaphragms or valves defective? ➔ Other than above mentioned causes?	✓ Connect hose or silencer to pump outlet. ✓ Perform maintenance. ✓ Contact local distributor. ✓ Clean and inspect pump heads. ✓ Replace diaphragms and/or valves. ✓ Contact local distributor.
☐ Pump seized.		✓ Contact local distributor.

Replacing diaphragms and valves

NOTICE

Maintenance of the pump must be performed only by a suitable trained and supervised personnel.

All bearings are encapsulated and are filled with long-life lubricant and run under normal operating conditions maintenance free. Replacement of the bearing must be performed at the latest at 90% of the nominal lifetime of the bearing (nominal lifetime of the bearing: 40000 operation hours, if necessary provide an operation time counter). If the noise level is enhanced switch off the pump immediately and inspect. Replacement of the bearing must only performed at the factory. Never operate a defective pump.

NOTICE

The valves and the diaphragms are wear parts. Replace the diaphragms at the latest at 90% of the typical lifetime or immediately at higher noise level. The typical lifetime of a diaphragm is 15000 operation hours when pumping non-corrosive or non-condensable gases, if necessary provide an operation time counter. Pumping condensable media, operation at high temperature, great gas flow or deposits reduce the lifetime of a diaphragm significantly. Therefore check the diaphragm regularly for faultless condition.

☞ Prevent internal condensation, transfer of liquids or dust. The diaphragm and valves will be damaged, if liquids are pumped in significant amount.

NOTICE

To control the faultless function of the diaphragm install a gas specific detector at the outlet of the inert gas flush. A signal from the detector indicates a diaphragm crack. Switch off the pump immediately and check diaphragms.

If the rated ultimate vacuum is no longer achieved, the pump interior, the hose connections, the diaphragms and the valves must be cleaned and the diaphragms and valves must be checked for cracks or other damage. Replace defective parts immediately.

NOTICE

Inspect the fittings of the pump regularly, the maximum permitted leak rate is 0.1 mbar l / s.

When replacing diaphragms and valves check the overpressure safety valve at the outlet of the pump.

Remove possible deposits in the crank chamber of the pump when replacing diaphragms and valves.

If the pump is exposed to corrosive gases or gases and vapors which may form deposits (e. g. crystallization) or deposits due to pumped particles, maintenance should be carried out frequently (according to the experience of the user).

☞ Regular maintenance will improve the lifetime of the pump and also protect both man and environment.



! WARNING

! CAUTION

Before starting maintenance vent the system, isolate the pump and other components from the vacuum system and the electrical supply. Drain condensate if applicable, avoid the release of pollutants. Allow sufficient cooling of the pump. Before starting maintenance, wait two minutes after isolating the equipment from mains to allow the capacitors to discharge.

Never operate the pump if covers or other parts of the pump are disassembled. Ensure that the pump cannot be operated accidentally. Never operate a defective or damaged pump.

Ensure that the maintenance technician is familiar with the safety procedures which relate to the products processed by the pumping system.

Attention: The pump might be contaminated with the process chemicals that have been pumped during operation. Ensure that the pump is decontaminated before maintenance and take adequate precautions to protect people from the effects of dangerous substances if contamination has occurred.

☞ Wear appropriate safety-clothing when you come in contact with contaminated components.

NOTICE

☞ **Please read section "Replacing diaphragms and valves" completely before starting maintenance.**

Partially the pictures show pumps in other versions. This doesn't influence replacing diaphragms and valves of the pump.

NOTICE

If not using genuine spare parts the ATEX conformity becomes invalid.

Attention: Open the pump only outside of explosive atmospheres and only by suitably trained personnel.

Always replace both diaphragms of a pump head!

Always disassemble and assemble only one pump head before opening the next pump head. The single parts of a pump head are measured in a way so to avoid that the diaphragm clamping disc hits the head cover. Never mix parts of different pump heads or add or leave out washers.

We recommend to replace all diaphragms and valves of a pump at the same time.

Spare parts

Set of seals (diaphragms, valves, O-rings)

MD 4C EX VARIO / MD 4C EX VARIO + AK + EK **2 x 20696837**

MV 10C EX VARIO / MV 10C EX VARIO + AK + EK **4 x 20696837**

Diaphragm key SW 66 (special tool) **20636554**

Sealing bond (PTFE) **20637514**

Valve (one piece)..... **20637225**


(internal overpressure valve in MD 4C EX VARIO , MD 4C EX VARIO + AK + EK,
MV 10C EX VARIO, MV 10C EX VARIO + AK + EK)

O-Ring 28 x 2.5 **20635628**

at the spherical ground joint of the catchpot at the inlet

Flat seal (overpressure safety valve at the outlet)..... **20637081**


Spring (overpressure safety valve at the outlet)..... **20637065**

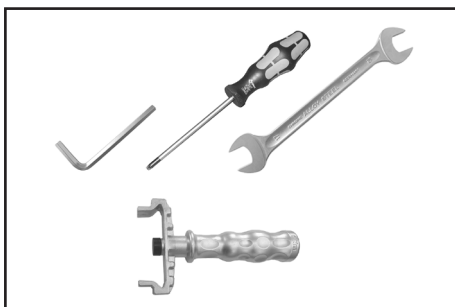


Tools required (metric):

- Phillips screw driver size 2
- Open-ended wrench size 10/15/17
- Open-ended wrench size 19, 4.5 mm thick
- Torque wrench size 19
- Hex key size 5
- Diaphragm key SW 66 (cat. no.: 20636554)
- Open-ended wrench size 19 with torque indicator


Cleaning and inspecting the pump heads

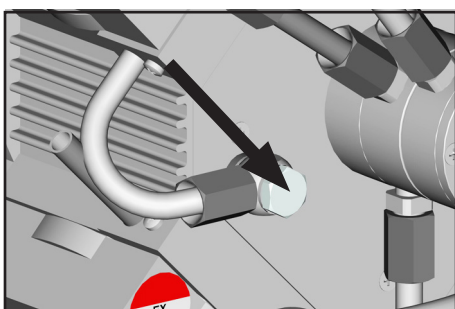


- ➔ Use open-ended wrench to unscrew the inert gas fitting at the housing cover.

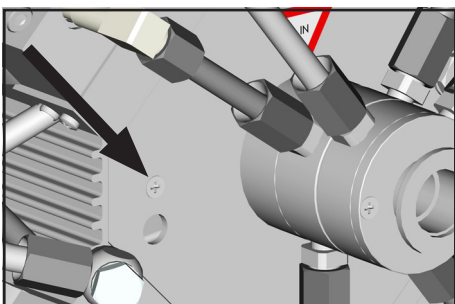


- ➔ Unscrew the housing plate from the pump.

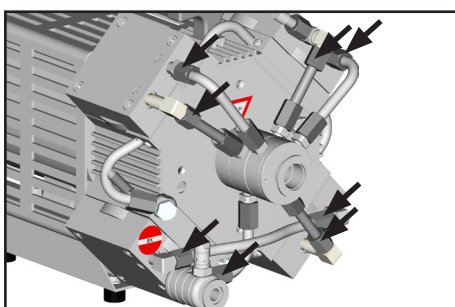

- ➔ Unscrew the union nut at the elbow fitting of the pump head and remove hose.
- ⚠ Turn the fittings as little as possible and do not remove the fittings from the pump head. Through reassembly a leak may result.

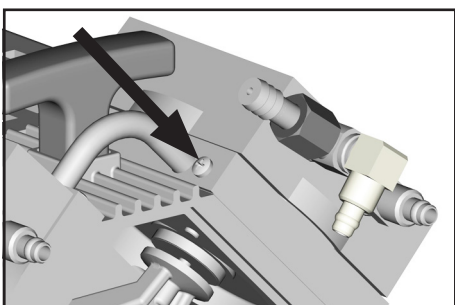


- ➔ Position the pump on the motor. Unscrew the locking plate at the inert gas connection and remove hose.

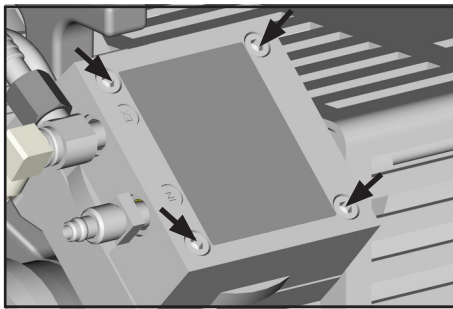


- ➔ Use hex key to remove four socket head screws from pump head and remove upper housing (housing cover with housing cover insert and head cover).
- ⚠ Never remove parts by using a spiky or sharp-edged tool (e. g. screw driver), we recommend to use a rubber mallet or compressed air (to be blown carefully into port).
- ➔ Carefully remove head cover from housing cover insert and check valves. Note position of valves and remove.
- ⚠ Replace valves if necessary.
- ⚠ Use petroleum ether or industrial solvent to remove deposits. Do not inhale.

Replacing the diaphragm

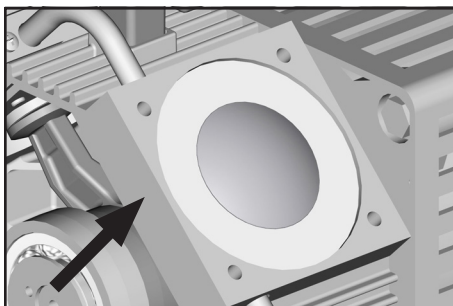


- ➔ Remove intermediate flange by hand.

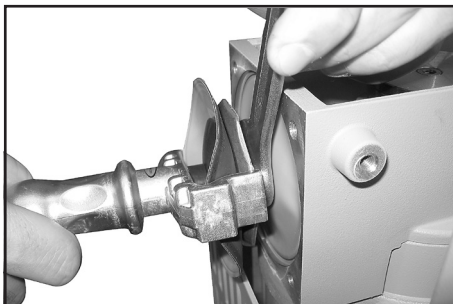

- ⚠ Check diaphragm for damage and replace if necessary.
- ➔ Lift diaphragm carefully sidewise.
- ⚠ Never use a spiky or sharp-edged tool to lift the diaphragm.
- ➔ Use the diaphragm key to grip the diaphragm support disc below the diaphragm.
- ➔ Unscrew diaphragm support disc with diaphragm and diaphragm clamping disc.


- ➔ Never increase the amount of washers.
Risk of temperature increase!
Risk of explosion!

⚠ DANGER

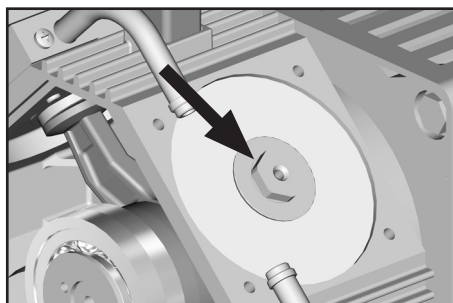


- ➔ Remove clamping disc of the safety diaphragm with an open-ended wrench and remove together with safety diaphragm and support disc.

View of the disassembled pump head parts (fig. MD 4C EX VARIO)

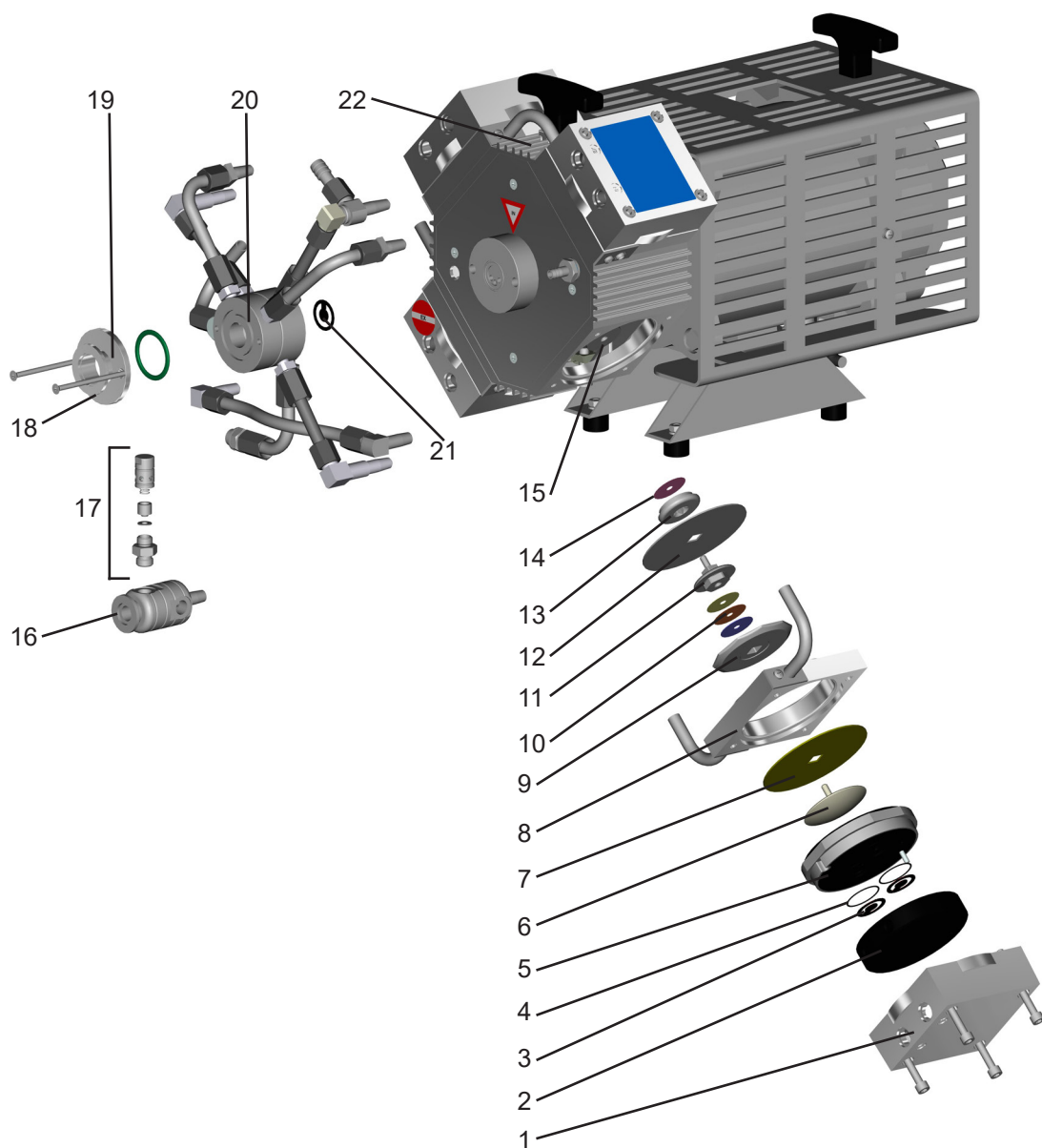


Position	Component	Position	Component
1	housing cover	12	safety diaphragm
2	housing cover insert	13	diaphragm support disc
3	valve	14	washer
4	O-ring	15	rod
5	head cover	16	outlet
6	diaphragm clamping disc	17	overpressure safety valve
7	work diaphragm	18	inlet
8	intermediate flange	19	distributor cover
9	diaphragm support disc	20	distributor
10	washer	21	internal overpressure safety valve
11	safety diaphragm clamping disc	22	housing

View of the disassembled pump head parts (fig. MV 10C EX VARIO)

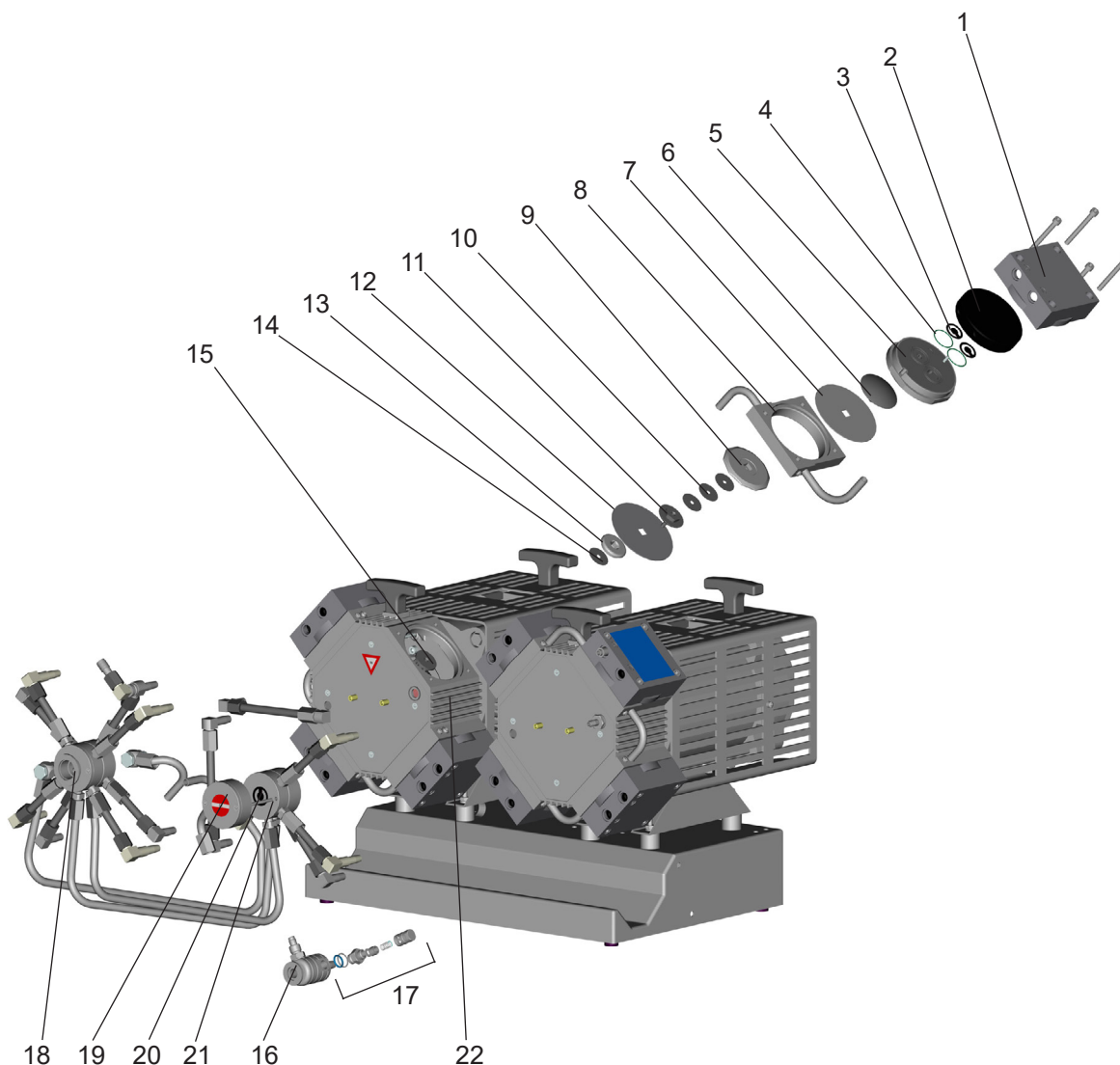


Position	Component	Position	Component
1	housing cover	12	safety diaphragm
2	housing cover insert	13	diaphragm support disc
3	valve	14	washer
4	O-ring	15	rod
5	head cover	16	outlet
6	diaphragm clamping disc	17	overpressure safety valve
7	work diaphragm	18	inlet
8	intermediate flange	19	distributor cover
9	diaphragm support disc	20	distributor
10	washer	21	internal overpressure safety valve
11	safety diaphragm clamping disc	22	housing

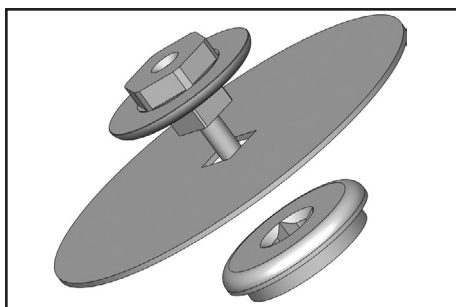

Before reassembling ensure that all parts are clean, dry and lint free, especially the valve seat.

- ➔ Position new safety diaphragm between safety diaphragm clamping disc and safety diaphragm support disc.
- ☞ Attention: Position diaphragm with pale side towards diaphragm clamping disc.
- ☞ Make sure that the square head screw of the diaphragm clamping disc is correctly seated in the guide hole of the diaphragm support disc.

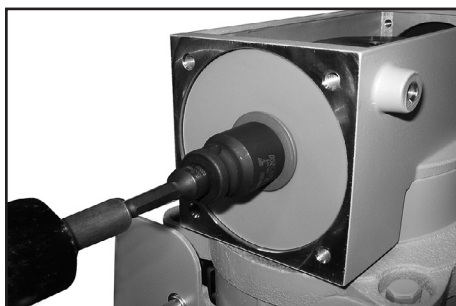


Use torque wrench w/f 19 to screw safety diaphragm support disc, safety diaphragm and safety diaphragm clamping disc to rod.

- ☞ **Attention: Absolutely assemble all washers correctly if available (between safety diaphragm support disc and rod). Never assemble more washers!**


! DANGER

- ☞ Make sure that the square head screw of the diaphragm clamping disc is correctly seated in the guide hole of the diaphragm support disc.

Optimum torque for the safety diaphragm clamping disc: **6 Nm.**

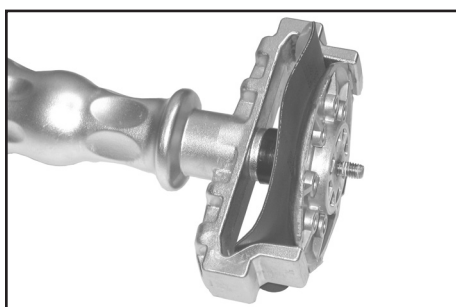


- ➔ Position new work diaphragm between diaphragm clamping disc with square head screw and diaphragm support disc.
- ☞ Attention: Position diaphragm with pale side towards diaphragm clamping disc (to pump chamber).
- ☞ Make sure that the square head screw of the diaphragm clamping disc is correctly seated in the guide hole of the diaphragm support disc.
- ➔ Lift diaphragm at the side and position carefully together with diaphragm clamping disc and diaphragm support disc in the diaphragm key.
- ☞ Avoid damage of the diaphragm: Do not bend diaphragm too much.

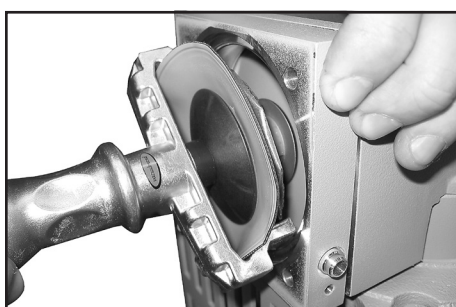

- ➔ Position intermediate flange to housing.
- ➔ Screw diaphragm clamping disc, diaphragm, diaphragm support disc and washers to connecting rod.
- ➔ Optimum torque for the diaphragm support disc: 6 Nm, it is recommended to use a torque key. Attach hex key to diaphragm key (hexagonal bolt 6 mm wide).
- Attention:** Never use the diaphragm key with any additional tools like tongs or hex keys without torque limitation.


Assembling the pump head

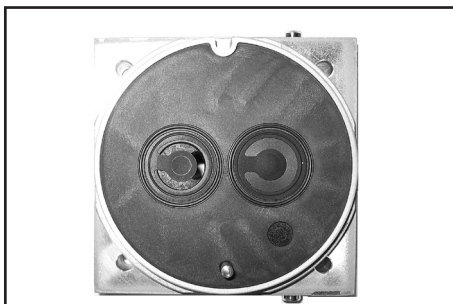


- ➔ Position pump in a way so that the disassembled pump head is above, support if necessary.
- ➔ By turning eccentric bushing (front of connecting rod), bring connecting rod into a position in which diaphragm is in contact with housing and centred with respect to bore.

Reassemble in reverse order.

- ➔ Install head cover with O-ring, valves and housing cover with housing cover insert.

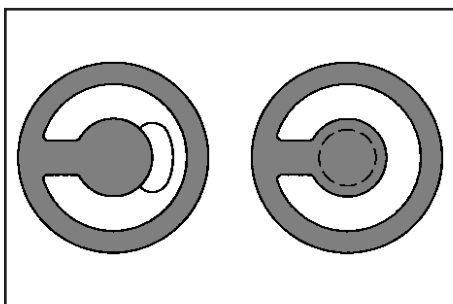


- ☞ Make sure that the valves are correctly seated: Valves at the outlet with round centred opening under valve, valves at the inlet with kidney-shaped opening beside valve.
- ☞ Ensure that the diaphragm is positioned centrally so that it will become clamped uniformly between housing and head cover.

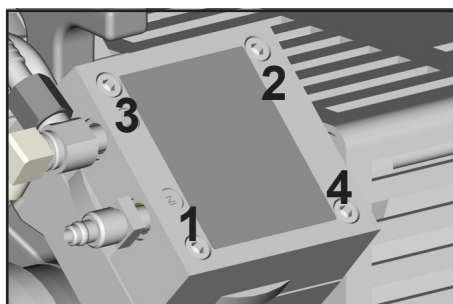

- ➔ Screw in four socket head screws fixing housing cover crosswise (e. g. in the sequence 1,2,3,4) first slightly, then tighten.
- ☞ Do not tighten until head cover is in contact with housing, torque 12 Nm.


Attention: Ensure that the diaphragm clamping disc never hits the head cover! In case of loud knocking switch off the pump immediately.

Inspect pump or replace diaphragms and valves at the other pump heads analogue.

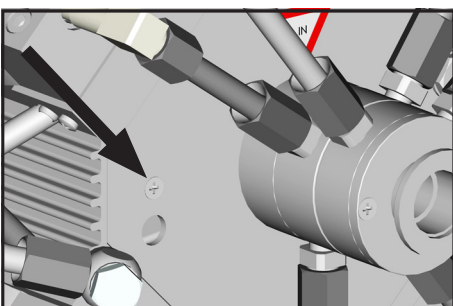


- ➔ Screw housing plate to pump.

Individual performance check of a pump head

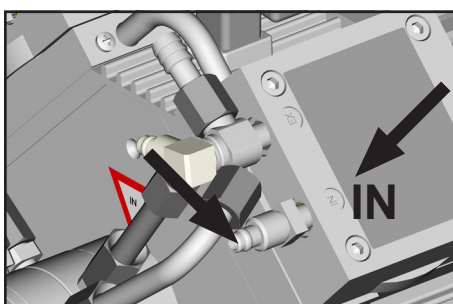


- ➔ Measure the pressure at the inlet port of the individual head: Use a suitable vacuum gauge (e. g. DVR 2pro, cat. no.: 20682906), make sure that it is correctly calibrated, and measure the pressure at the inlet port. A vacuum of less than 120 mbar should be indicated.
- ☞ If the reading is higher, recheck the pump chamber and make sure that the valves and the diaphragms are correctly seated (diaphragms concentric with bore).


Assembling fittings

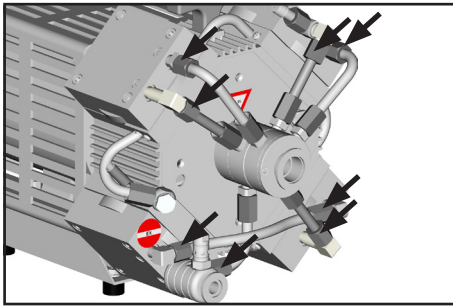

- ➔ Position pump on pump feet.
- ➔ Connect hose to elbow fitting.
- ➔ Tighten union nuts first by hand and then tighten one full turn using open ended wrench.

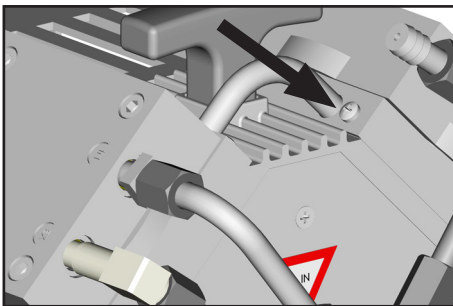

- ➔ Position the pump on the motor.
Attach hose to the inert gas connection and screw the locking plate.

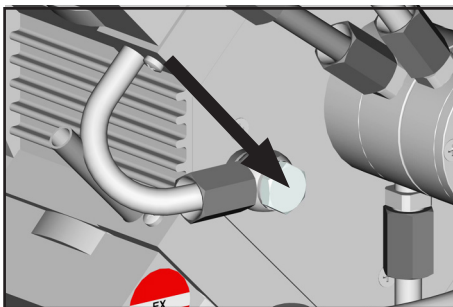

- ➔ Use open-ended wrench to screw the inert gas fitting at the housing cover.

DANGER

Attention: Perform a leak test always after opening the pump by using an appropriate leak detector (e. g. helium leak detector)!

Ensure that the diaphragm clamping disc never hits the head cover while pump is running! Noise check!

If the pump does not achieve the ultimate pressure:

- ☞ In case the diaphragms and valves have been replaced, a run-in period of several hours is required before the pump achieves its ultimate vacuum.


If all pump heads achieve a vacuum below 120 mbar but pump does not achieve the ultimate total pressure:

Check hose connectors between pump heads and manifolds for leaks. If necessary recheck pump chamber.


Inspecting the overpressure safety valve

Inspecting the overpressure safety valve at the outlet:


- ➔ Unscrew the overpressure safety valve at the outlet of the pump.

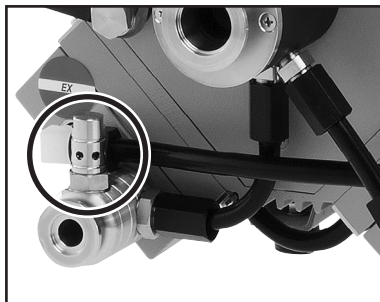

MD 4C EX VARIO

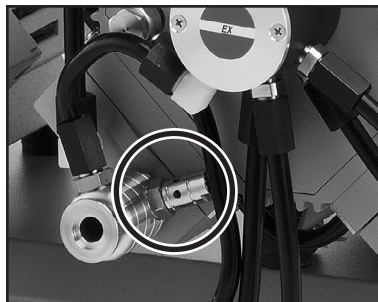


MV 10C EX VARIO

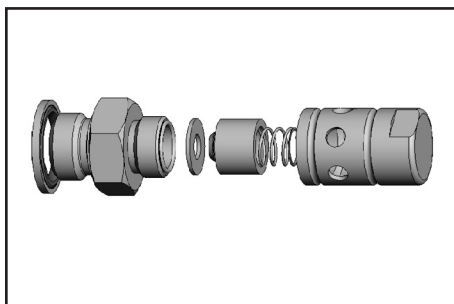

- ➔ Disassemble the overpressure safety valve and inspect the parts for faultless condition, replace damaged parts.
- ➔ Reassemble the valve correctly.
- ➔ Screw the overpressure safety valve to the outlet of the pump.

Inspection and replacement of the internal overpressure safety valve at the manifold: (only MD 4C EX VARIO / MD 4C EX VARIO + AK + EK)

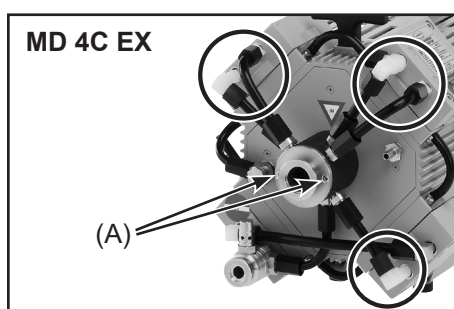

- ➔ Use an open-ended wrench (width 17 mm) to unscrew at the pump head the union nuts of the tubings which are connected to the manifold (5x).


Elbow fitting (3x):

- ➔ Turn the fitting with an open-ended wrench (width 14 mm) to detach the tubing from the pump head (1/4 of a turn at maximum).
- ⚠ Do not remove the elbow fitting from the pump head.

Straight fitting (2x):

- ➔ Pull the tubing off the fitting.

- ➔ Unscrew the two countersunk screws (A) at the cover plate and remove cover plate together with distributor and tubing.

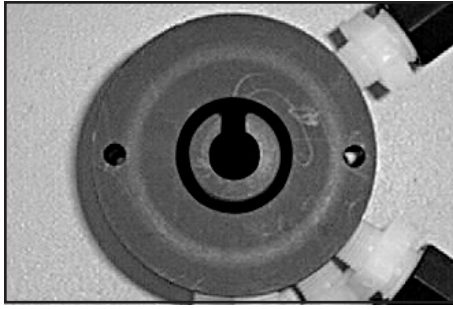

- ➔ Note position of valve and remove.
- ☞ Check valve for damage and replace if necessary. Make sure that the valve is correctly seated.
- ➔ Reassemble distributor and cover plate.

Elbow fitting (3x):


- ➔ Slip the tubing onto the elbow fitting by turning the fitting with an open-ended wrench (width 15 mm).

Straight fitting (2x):


- ➔ Slip the tubing onto the fitting.

- ➔ Tighten union nuts first by hand and then tighten **one** full turn using the open ended wrench (width 17 mm).

Inspection and replacement of the internal overpressure safety valve at the manifold at the outlet: (only MV 10C EX VARIO / MV 10C EX VARIO + AK + EK)

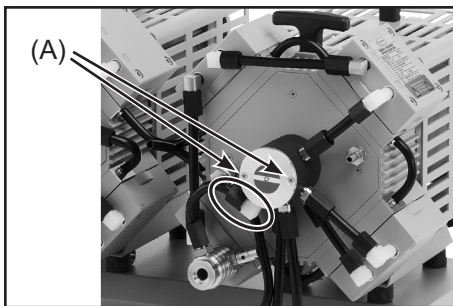

- ➔ Use an open-ended wrench (width 17 mm) to unscrew at the pump head the union nut of the tubing which is connected to the outlet.
- ➔ Turn the fitting with an open-ended wrench (width 14 mm) to detach the tubing from the fitting (1/4 of a turn at maximum).
- ☞ Do not remove the elbow fitting from the distributor cover.
- ➔ Unscrew the two countersunk screws (A) at the cover plate and remove cover plate together with distributor cover.

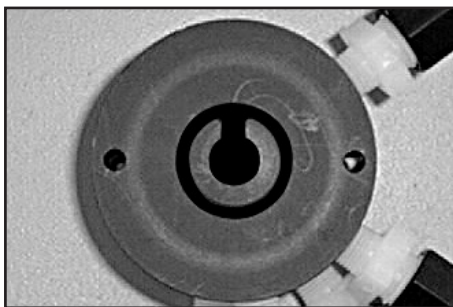

- ➔ Note position of valve and remove.
- ☞ Check valve for damage and replace if necessary. Make sure that the valve is correctly seated.
- ➔ Reassemble distributor cover and cover plate.
- ➔ Slip the tubing onto the elbow fitting by turning the fitting with an open-ended wrench (width 15 mm).
- ➔ Tighten union nut first by hand and then tighten **one** full turn using the open ended wrench (width 17 mm).

Notes on assembling fittings


The threads of the fittings at the pump heads might be sealed with PTFE bond (width 10 mm or 5 mm, thickness 0.1 mm).

If the fittings are turned or have been removed, use new PTFE bond when reassembling.

If not using new fittings clean the thread of the fittings from PTFE bond.


Wind the thread of the fitting two times against the direction of turning in (anticlockwise) with PTFE bond. Tighten the PTFE bond firmly. Ensure that the PTFE bond ends flush with the fittings. Ensure that the PTFE bond not covers the openings in the elbow fittings.

Screw in elbow fittings (clockwise) until only half of a thread course is visible. Do not overtighten. Avoid canting of the thread.

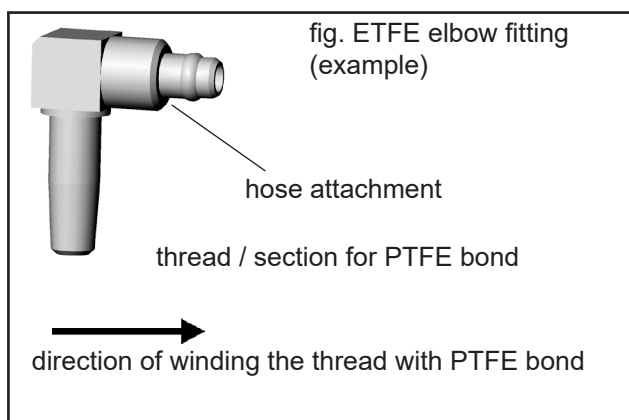

Seal only the lower half of the thread of the stainless steel fittings with PTFE bond (width 5 mm), see figure. Sealing the complete thread with PTFE bond leads to interruption of the conductive connection!

Never seal the hose attachment of metal fittings with PTFE bond.

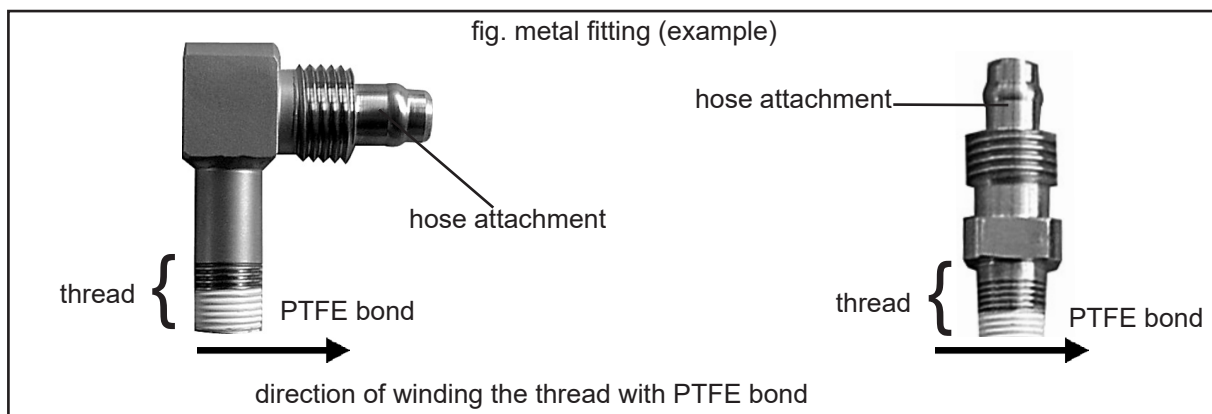

Repair - Maintenance - Return - Calibration

IMPORTANT

Every employer (user) is held responsible for the health and safety of his employees. This also applies to service personnel performing repair, maintenance, return or calibration.

The **health and safety clearance form** informs the contractor about any possible contamination of the device and forms the basis for the risk assessment.

In case of devices which have been in contact with biological substances of risk level 2 contact the VACUUBRAND service absolutely before dispatching the device. These devices have to be completely disassembled and decontaminated by the user prior to shipment. Do not return devices which have been in contact biological substances of risk level 3 or 4. These devices cannot be checked, maintained or repaired. Also decontaminated devices must not returned to VACUUBRAND due to a residual risk.

The same conditions apply to on-site work.

No repair, maintenance, return or calibration is possible unless the correctly completed health and safety clearance form is returned. Devices sent are rejected if applicable. Send a completed copy of the **health and safety clearance form** to us in advance. The declaration must arrive before the equipment. Enclose a second completed copy with the product.

Remove all components from the device that are no original VACUUBRAND components. VACUUBRAND will not be responsible for lost or damaged components that are no original components.

Drain the device completely of fluids and residues. Decontaminate the device.

Close all openings airtight especially if using substances hazardous to health.

To expedite repair and to reduce costs, please enclose a detailed description of the problem and the product's operating conditions with every product returned.

If you do not wish a repair on the basis of our **quotation**, the device may be returned to you disassembled and at your expense.

In many cases, the components must be cleaned in the factory prior to repair.

For cleaning we use an environmentally friendly water based process. Unfortunately the combined attack of elevated temperature, cleaning agent, ultrasonic treatment and mechanical stress (from pressurised water) may result in damage to the paint. Please mark in the health and safety clearance form if you wish a repaint at your expense just in case such a damage should occur. We will also replace parts for cosmetic reasons at your request and at your expense.

Before returning the device

Pack the device properly, if necessary, please order original packaging materials at your costs.

Mark the package completely

Enclose the completed health and safety clearance form.

Notify the carrier of any possible contamination if required.

Scrapping and waste disposal

Dispose of the equipment and any components removed from it safely in accordance with all local and national safety and environmental requirements. Particular care must be taken with components and waste oil which have been contaminated with dangerous substances from your processes. Do not incinerate fluoroelastomer seals and O-rings. You may authorize us to dispose of the equipment **at your expense**. Otherwise we return the device at your expense.

Health and safety clearance form

1. Device (Model):
2. Serial no.:
3. Reason for return / malfunction:
4. Has the device been used in a copper process step (e.g., semiconductor production): ☐ yes ☐ no
5. Substances (gases, liquids, solids, biological material, e. g. bacteria, viruses) in contact with the device / which have been pumped:
.....
.....
.....
6. Risk level of the used biological material: ☐ none ☐ 1 ☐ 2* ☐ 3** ☐ 4**
 * Contact the VACUUBRAND service absolutely before dispatching the device.
 ** Devices which have been in contact with biological substances of risk level 3 or 4 cannot be checked, maintained or repaired. Also decontaminated devices must not returned to VACUUBRAND due to a residual risk.
7. Radioactive contamination: ☐ yes ☐ no
8. Prior to return to the factory the device has been decontaminated: ☐ yes ☐ no
 Description of the decontamination method and the test / verification procedure:

9. All parts of the device are free of hazardous, harmful substances: ☐ yes ☐ no
10. Protective measures required for service staff:
.....
11. If the paint is damaged, we wish a repaint or a replacement of parts for reason of appearance (repaint and replacement at customer's expense): ☐ yes ☐ no

12. Legally binding declaration

We assure for the returned device that all substances, which have been in contact with the device are listed in section 5 and that the information is complete and that we have not withheld any information. We declare that all measures - where applicable - have been taken listed in section "Repair - Maintenance - Return - Calibration". By our signature below, we acknowledge that we accept liability for any damage caused by providing incomplete or incorrect information and that we shall indemnify VACUUBRAND from any claims as regards damages from third parties. We are aware that as expressed in § 823 BGB (Public Law Code of Germany) we are directly liable for injuries or damages suffered by third parties, particularly VACUUBRAND employees occupied with handling/repairing the product. Shipping of the device must take place according to regulations.

Name: Signature:

Job title: Company's seal:

Date:

Release for repair grant by VACUUBRAND (date / signature):

Protective measures: ☐ Protective gloves, safety goggles ☐ Hood ☐ External cleaning

VACUUBRAND GMBH + CO KG
Alfred-Zippe-Straße 4
97877 Wertheim - Germany

T +49 9342 808-5660 F +49 9342 808-5666
E-Mail: service@vacuubrand.com
www.vacuubrand.com



EC Type examination certificate



APRAZ

a.s.b.l.

Your inspecting authority



- (1) **EU TYPE EXAMINATION CERTIFICATE**
- (2) **Equipment or Protective system intended for use
in potentially explosive atmospheres
Directive 2014/34/EU**
- (3) EU type examination certificate number: **APRAZ 20ATEX0209 X**
- (4) Equipment or protective system: Three-phase and single-phase asynchronous motors
Single-phase: BA(V or X) 3xy C or E or CP
Three-phase: BA(V or X) 3xy TR
- (5) Applicant - Manufacturer: Elnor Motors NV Belgium
- (6) Address: De Costerstraat 45
3150 Haacht (Wespelaar)
Belgium
- (7) This equipment or protective system and any acceptable variation thereto is specified in the schedule to this certificate and the documents therein referred to.
- (8) APRAZ, notified body n° 0029 in accordance with article 17 of the directive 2014/34/EU of 26 February 2014, certifies that this equipment or protective system has been found in compliance with the essential health and safety requirements relating to the design and construction of equipment and protective systems intended for use in potentially explosive atmospheres, given in annexe II of the directive.

The examination and test results are recorded in the confidential file n° 17566/006.
- (9) The conformity of the essential health and safety requirements has been assured by compliance with:
EN 60079-0: 2018
EN 60079-1: 2014
- (10) If the sign "X" is placed after the certificate number, it indicates that the equipment or protective system is subject to special conditions for safe use specified in the schedule to this certificate.
- (11) This EU TYPE EXAMINATION CERTIFICATE relates only to the design and construction of this specified equipment or protective system in accordance with the directive 2014/34/EU. Further requirements of the directive apply to the manufacturing and supply of this equipment or protective system. These are not covered in this certificate.
- (12) The marking of the equipment or protective system shall include the following:
⊕ II 2 G Ex db IIB T4...T3 Gb

Brussels, 07 December 2020

APRAZ
Chaussée de Vilvorde 156
1120 Bruxelles
tel 02/264 03 60 fax 02/268 89 58
mail : info@apraz.com

Ir. B. Nève
General Manager

This certificate can only be reproduced in its entirety, schedule included, without any modification

Rev. 0

1/3

R.A. 001 rev2 - 08-2020



APRAGAZ
a.s.b.l.
Your inspecting authority



(13)

SCHEDULE

(14)

EU TYPE EXAMINATION CERTIFICATE N° APRAGAZ 20ATEX0209 X

(15) Description of the equipment or protective system

Three-phase and single-phase asynchronous motors:

Single-phase: BA(V or X) 3xy C or E or CP

Three-phase: BA(V or X) 3xy TR

BA: a naturally cooled explosion proof motor

BAV: a fan cooled explosion proof motor

BAX: a motor cooled by the application

The motor consists of a flameproof enclosure which is made of an aluminium or a steel sheet metal carcass and has aluminium or cast iron end shields.

The fastenings are made by hex cap screw of property class 4.6, nuts of property class 5.0 and socket cap screws of property class 8.8.

The motor may be equipped with two ends shafts and can be electrically directly fed by a single cable or through a connection in the terminal box.

Temperature classes

The motors on S1 duty service cooled by the process have the temperature class T3. The motors have a temperature class T4 if they are equipped with a thermal protection suitable to the temperature class.

The motors with natural cooling, the motors energized by converters and the motors on S2 to S9 duty service are provided with a direct thermal protection or a PTC resistance suitable to the temperature class.

Electrical characteristics

Power range: from 0,25 kW to 2,5 kW

Voltage supply: up to 690 Volt – For VAC motors: 50/60 Hz

Insulation class: F

Routine tests

The manufacturer shall make the routine verifications and tests necessary to ensure that the electrical apparatus produced complies with the specifications submitted to the testing station together with the prototype or sample.

Every part of the motor legible for routine testing whether it's made from aluminium, iron cast or steel that is tested at 4*Pref is exempted from overpressure pressure testing, otherwise the routine tests at 11,01 bar need to be executed.

The control of the weld of the steel carcass is always to be tested at 11,01 bar.

A motor of every fabrication number is additionally tested on the test bench. This to verify the maximum surface temperatures on different parts on the motor for the specific type of motor.

Cables resisting up to 120°C should be selected for a T3 and cables resisting up to 100°C should be selected for a T4 motor.



This certificate can only be reproduced in his entirely, schedule included, without any modification

Rev. 0

2/3

R.A. 001 rev2 – 08-2020



APRAGAZ

a.s.b.l.
Your inspecting authority



- (13) **SCHEDULE**
- (14) **EU TYPE EXAMINATION CERTIFICATE N° APRAGAZ 20ATEX0209 X**

- (15) Marking:

Elnor Motors NV Belgium
De Costerstraat 45
3150 Haacht (Wespelaar)
Belgium

Year CE + Identification n° of NOBO production control phase
Type: BA(X/V) 3xy
Serial number: 650xxxxx
Manufacturer date, fabrication number and batch number : JJWW/NFNF xxx



II 2 G Ex db IIB T4...T3 Gb
APRAGAZ 20ATEX0209 X

Warning:

Do not open under voltage.

Eventual prescriptions

Choice of cable needs to be compatible with a temperature of X °C.
Ambient temperature range: -40°C up to +70°C.
An adequate warning is to be printed when the motor is delivered without cable glands, plugs or cable.

Note:

This X °C is in function of the temperature class of the motor. X=120°C for a T3 motor and X=100°C for a T4 motor.
Depending on the foreseen duty class of the motor the temperature range can be narrowed.

- (16) The examination and test results are recorded in confidential report n° U194718.

- (17) Special conditions for safe use:

An appropriate connection of the free end of the cable shall be foreseen in case the motor is provided with a permanently connected cable.

If operated with a frequency drive, or when a temperature class better than a T3 has to be respected a thermal protection device has to be installed directly in the motor. The following table details the maximum set points of the direct thermal protection device per temperature class:

Temperature class	Set point windings in °C	Set point end shields/bearings in °C
T3	150	135
T4	135	120

- (18) Essential health and safety requirements:

The essential health and safety requirements are covered by the conformity to European standards listed at (9).

This certificate can only be reproduced in his entirely, schedule included, without any modification

Rev. 0

R.A. 001 rev2 – 08-2020

3/3

EG-Konformitätserklärung EC Declaration of Conformity Déclaration CE de conformité



Hersteller / Manufacturer / Fabricant:

VACUUBRAND GMBH + CO KG · Alfred-Zippe-Str. 4 · 97877 Wertheim · Germany

Hiermit erklärt der Hersteller, dass das Gerät konform ist mit den Bestimmungen der Richtlinien:

Hereby the manufacturer declares that the device is in conformity with the directives:

Par la présente, le fabricant déclare, que le dispositif est conforme aux directives:

- 2014/34/EU
- 2011/65/EU, 2015/863
- 2006/42/EG (gilt nur für Motor / applies only for motor / ne s'applique qu'au moteur)

Membranvakuumpumpe / Diaphragm vacuum pump / Pompe à membrane:

Typ / Type / Type: MD 4C EX VARIO / MD 4C EX VARIO + AK + EK / MV 10C EX VARIO / MV 10C EX VARIO + AK + EK

Artikelnummer / Order number / Numéro d'article: 20696936 / 20696937 / 20696946 / 20696947

Seriennummer / Serial number / Numéro de série: Siehe Typenschild / See rating plate / Voir plaque signalétique

Angewandte harmonisierte Normen / Harmonized standards applied / Normes harmonisées utilisées:

EN 1127-1:2019, EN ISO 80079-36:2016 (ISO 80079-36:2016), EN IEC 63000:2018 (IEC 63000:2016)

Motor / motor / moteur: EN 60204-1:2018 (IEC 60204-1:2016), EN IEC 60079-0:2018 (IEC 60079-0:2017),

EN 60079-1:2014 (IEC 60079-1:2014)

Bevollmächtigter für die Zusammenstellung der technischen Unterlagen / Person authorised to compile the technical file /
Personne autorisée à constituer le dossier technique:

Dr. Constantin Schöler · VACUUBRAND GMBH + CO KG · Germany

Die technische Dokumentation gemäß der Richtlinie 2014/34/EU (ATEX) wird im Auftrag von VACUUBRAND GMBH + CO KG bei der entsprechend beauftragten Hinterlegungsstelle hinterlegt:

The technical documentation in accordance with directive 2014/34/EU (ATEX) is deposited on behalf of VACUUBRAND GMBH + CO KG with the relevant authorized depository:

La documentation technique conformément à la directive 2014/34/EU (ATEX) est déposée au nom de VACUUBRAND GMBH + CO KG auprès de l'institution de dépôt mandatée à cet effet:

IBExU Institut für Sicherheitstechnik GmbH · Fuchsmühlenweg 7 · 09599 Freiberg · Germany

Ort, Datum / place, date / lieu, date: Wertheim, 21.07.2025

(Dr. Constantin Schöler)

Geschäftsführer / Managing Director / Gérant

ppa.

(Jens Kaibel)

*Technischer Leiter / Technical Director /
Directeur technique*

VACUUBRAND GMBH + CO KG

Alfred-Zippe-Str. 4
97877 Wertheim

Tel.: +49 9342 808-0

Fax: +49 9342 808-5555

E-Mail: info@vacuubrand.com

Web: www.vacuubrand.com

VACUUBRAND®

Einbauerklärung für Maschinen
Declaration of Incorporation of the Machinery
Déclaration d'incorporation des machines

Hersteller / Manufacturer / Fabricant:

VACUUBRAND GMBH + CO KG · Alfred-Zippe-Str. 4 · 97877 Wertheim · Germany

Hiermit erklärt der Hersteller, dass die unvollständige Maschine konform ist mit den Bestimmungen der Richtlinie:

Hereby the manufacturer declares that the incomplete machinery is in conformity with the directive:

Par la présente, le fabricant déclare, que la quasi-machine est conforme à directive:

■ 2006/42/EG

Membranvakuumpumpe / Diaphragm vacuum pump / Pompe à membrane:

Typ / Type / Type: MD 4C EX VARIO / MD 4C EX VARIO + AK + EK / MV 10C EX VARIO / MV 10C EX VARIO + AK + EK

Artikelnummer / Order number / Numéro d'article: 20696936 / 20696937 / 20696946 / 20696947

Seriennummer / Serial number / Numéro de série: Siehe Typenschild / See rating plate / Voir plaque signalétique

Angewandte harmonisierte Normen / Harmonized standards applied / Normes harmonisées utilisées:

EN ISO 12100:2010 (ISO 12100:2010), EN 1012-2:1996 + A1:2009

Weitere angewandte Normen / Other standards applied / Autres normes utilisées:

EN 61010-1:2010 + A1:2019 + A1:2019/AC:2019 (IEC 61010-1:2010 + COR:2011 + A1:2016, modifiziert / modified / modifié + A1:2016/COR1:2019)

Die technische Dokumentation nach Anhang VII B wurde erstellt. Der Hersteller verpflichtet sich, die technische Dokumentation zur unvollständigen Maschine den zuständigen Stellen in Papierform auf Verlangen zu übermitteln.

The technical documentation in accordance with annex VII B has been compiled. The manufacturer undertakes to submit the technical documentation relating to the incomplete machine to the relevant national authorities as paper mold on request.

La documentation technique selon l'annexe VII B a été établie. Le fabricant s'engage à remettre la documentation technique concernant la quasi-machine aux services compétents sous forme papier à leur demande.

Bevollmächtigter für die Zusammenstellung der technischen Unterlagen / Person authorised to compile the technical file / Personne autorisée à constituer le dossier technique:

Dr. Constantin Schöler · VACUUBRAND GMBH + CO KG · Germany

Die Inbetriebnahme dieser unvollständigen Maschine ist so lange untersagt, bis festgestellt wurde, dass die Maschine, in die sie eingebaut werden soll, den Bestimmungen der EG-Richtlinie Maschinen, den harmonisierten Normen, europäischen Normen oder den entsprechenden nationalen Normen entspricht.

This incomplete machine must not be put into service until the machinery into which it is to be incorporated has been declared in conformity with the provisions of the EC Machinery Directive, the harmonized standards, European standards, or the relevant national standards.

La mise en service de cette quasi-machine est interdite tant qu'il n'a pas été constaté que la machine dans laquelle elle doit être incorporée est conforme aux dispositions de la directive CE Machines, aux normes harmonisées, aux normes européennes ou aux normes nationales correspondantes.

Ort, Datum / place, date / lieu, date: Wertheim, 21.07.2025



(Dr. Constantin Schöler)

Geschäftsführer / Managing Director / Gérant

ppa.



(Jens Kaibel)

Technischer Leiter / Technical Director /
Directeur technique

VACUUBRAND GMBH + CO KG

Alfred-Zippe-Str. 4
97877 Wertheim

Tel.: +49 9342 808-0
Fax: +49 9342 808-5555
E-Mail: info@vacuubrand.com
Web: www.vacuubrand.com

VACUUBRAND®

DECLARATION OF CONFORMITY – China RoHS 2

VACUUBRAND GMBH + CO KG has made reasonable efforts to ensure that hazardous materials and substances may not be used in its products.

In order to determine the concentration of hazardous substances in all homogeneous materials of the subassemblies, a “Product Conformity Assessment” (PCA) procedure was performed. As defined in GB/T 26572 the “Maximum Concentration Value” limits (MCV) apply to these restricted substances:

- Lead (Pb): 0.1%
- Mercury (Hg): 0.1%
- Cadmium (Cd): 0.01%
- Hexavalent chromium (Cr(+VI)): 0.1%
- Polybrominated biphenyls (PBB): 0.1%
- Polybrominated diphenyl ether (PBDE): 0.1%

Environmentally Friendly Use Period (EFUP)

EFUP defines the period in years during which the hazardous substances contained in electrical and electronic products will not leak or mutate under normal operating conditions. During normal use by the user such electrical and electronic products will not result in serious environmental pollution, cause serious bodily injury or damage to the user's assets. The Environmentally Friendly Use Period for VACUUBRAND products is 40 years.



MATERIAL CONTENT DECLARATION FOR VACUUBRAND PRODUCTS						
有毒有害物质或元素 Hazardous substances						
部件名称 Part name	铅 Pb	汞 Hg	镉 Cd	六价铬 Cr(+VI)	多溴联苯 PBB	多溴二苯醚 PBDE
包装 Packaging	O	O	O	O	O	O
塑料外壳 / 组件 Plastic housing / parts	O	O	O	O	O	O
真空油 Vacuum oil	O	O	O	O	O	O
电池 Battery	O	O	O	O	O	O
玻璃 Glass	X	O	O	O	O	O
电子电气组件 Electrical and electronic parts	X	O	O	O	O	O
控制器 / 测量设备 Controller / measuring device	X	O	O	O	O	O
金属外壳 / 组件 Metal housing / parts	X	O	O	O	O	O
电机 Motor	X	O	O	O	O	O
配件 Accessories	X	O	O	O	O	O
此表格是按照SJ/T 11364-2014中规定所制定的。 This table is created according to SJ/T 11364-2014.						

VACUUBRAND®

- O: 表示该有毒有害物质在该部件所有均质材料中的含量均在GB/T 26572规定的限量要求以下。
O: Indicates that the above mentioned hazardous substance contained in all homogeneous materials of the part is below the required limit as defined in GB/T 26572.
- X: 表示该有毒有害物质至少在该部件某一均质材料中的含量超出GB/T 26572规定的限量要求。
X: Indicates that the above mentioned hazardous substance contained in at least one of the homogeneous materials of this part is above the required limit as defined in GB/T 26572.

电池、玻璃器皿和配件可能不属于所附设备所包含的内容，它们可能有各自单独的EFUP标记和/或可能正在维护其部件EFUP标记的更新。

Batteries, glassware and accessories might not be content of the enclosed device and may have its own EFUP-marking and/or might be maintaining parts with changing EFUP-marking.

除上表所示信息外，还需声明的是，这些部件并非是有意图用铅（Pb）、汞（Hg）、铬（Cd）、六价铬（Cr(+VI)）、多溴联苯（PBB）或多溴二苯醚（PBDE）来制造的。

Apart from the disclosures in the above table, the subassemblies are not intentionally manufactured or formulated with lead (Pb), mercury (Hg), cadmium (Cd), hexavalent chromium (Cr+VI), polybrominated biphenyls (PBB), and polybrominated diphenyl ethers (PBDE).

Products manufactured by VACUUBRAND may enter into further devices (e.g., rotary evaporator) or can be used together with other appliances (e.g., usage as booster pumps).

With these products and appliances in particular, please note the EFUP labeled on these products.


VACUUBRAND will not take responsibility for the EFUP of those products and appliances.

Place, date: Wertheim, 06 September 2022



(Dr. Constantin Schöler)
Managing Director

ppa.

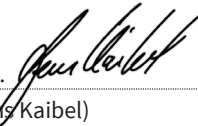

(Jens Kaibel)

Technical Director

VACUUBRAND GMBH + CO KG

Alfred-Zippe-Str. 4
97877 Wertheim
Germany

Tel.: +49 9342 808-0

Fax: +49 9342 808-5555


E-Mail: info@vacuubrand.com

Web: www.vacuubrand.com

Declaration of Conformity – China RoHS 2

V5_September 2022

Copyright 2022

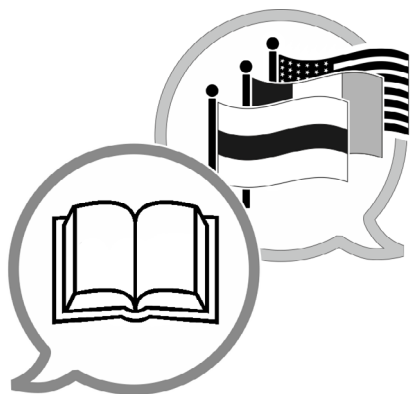

www.vacuubrand.com/manuals

Manufacturer:

VACUUBRAND GMBH + CO KG
Alfred-Zippe-Str. 4
97877 Wertheim
GERMANY

Phone:

Head office +49 9342 808-0
Sales +49 9342 808-5550
Service +49 9342 808-5660

Fax: +49 9342 808-5555

Email: info@vacuubrand.com

Web: www.vacuubrand.com