ELEKTROMAGNETISCHE SAUGLEITUNGSVENTILE SOLENOID OPERATED IN-LINE VALVES ELECTROVANNES D'ISOLEMENT

VV 6 - VV 6C - VV 15C - VV 25

VV-B 6 - VV-B 6C - VV-B 15C

Instructions for use

Inhalt / Contents / Contenu

Unbedingt beachten!	11
⊕ Hinweise zur Gerätekennzeichnung (ATEX).....	12
Technische Daten	13
Medienberührte Werkstoffe	13
Ersatzteile	14
Safety information!	15
⊕ Important information: Equipment marking (ATEX)	16
Technical Data	17
Wetted parts	17
Spare parts	18
Notes importantes!	19
⊕ Notes importantes concernant le marquage des appareils (ATEX).....	20
Données techniques	21
Matériaux exposés au gaz	21
Pièces de rechange	22

Trademark Index:

VACUU-LAN® (US-Reg.No 3,704,401), VACUU-BUS®, VACUU-CONTROL®, VACUU® (US-Reg. No 5,522,262), VACUU-SELECT® (US-Reg.No 5,522,260), VARIO® (US-Reg.No 3,833,788), VACUUBRAND® (US-Reg.No 3,733,388), VACUU-VIEW®, GREEN VAC® (US-Reg.No. 4,924,553), VACUU-PURE® (US-Reg No. 5,559,614) und die gezeigten Firmenlogos sind registrierte Trademarks der Firma VACUUBRAND GMBH + CO KG in Deutschland und/oder anderen Ländern.

Trademark index:

VACUU-LAN® (US-Reg.No 3,704,401), VACUU-BUS®, VACUU-CONTROL®, VACUU-SELECT® (US-Reg.No 5,522,260), VARIO® (US-Reg.No 3,833,788), VACUUBRAND® (US-Reg.No 3,733,388), VACUU-VIEW®, GREEN VAC® (US-Reg.No. 4,924,553), VACUU-PURE® (US-Reg No. 5,559,614) and also the shown company logos are registered trademarks of VACUUBRAND GMBH + CO KG in Germany and/or other countries.

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VACUU-LAN® (US-Reg.No 3,704,401), VACUU-BUS®, VACUU-CONTROL®, VACUU® (US-Reg. No 5,522,262), VACUU-SELECT® (US-Reg.No 5,522,260), VARIO® (US-Reg.No 3,833,788), VACUUBRAND® (US-Reg.No 3,733,388), VACUU-VIEW®, GREEN VAC® (US-Reg.No. 4,924,553), VACUU-PURE® (US-Reg No. 5,559,614) et les logos de l'entreprise sont des marques déposées ou des marques de VACUUBRAND GMBH + CO KG en Allemagne et/ou dans d'autres pays.

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➔ DANGER indicates a hazardous situation which, if not avoided, will result in death or serious injury.

➔ WARNING indicates a hazardous situation which, if not avoided, could result in death or serious injury.

• CAUTION indicates a hazardous situation which, if not avoided, could result in minor or moderate injury.

NOTICE is used to address practices not related to personal injury.

Electronic components must not be disposed of in the domestic waste at the end of their service life. Used electronic devices contain harmful substances that can cause damage to the environment or human health. End users are legally obliged to take used electric and electronic devices to a licensed collection point.

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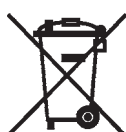
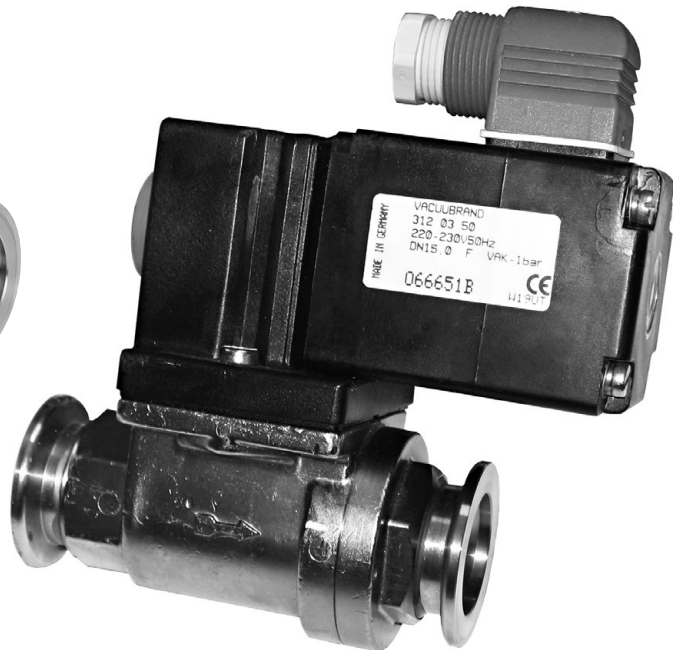
EN

Attention: This manual is not available in all languages of the EU. The user must not operate the device if he does not understand this manual. In this case a technically correct translation of the complete manual has to be available. The manual must be completely read and understood before operation of the device and all required measures must be applied.  "Safety instructions for vacuum equipment"

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Safety information!

NOTICE

The valves VV-B 6 / VV-B 6C and VV-B 15C are to be connected only to the VACUUBRAND BUS system VACUU•BUS!

Remove all packing material, remove the product from its packing-box, remove the protective covers, inspect the equipment.

If the equipment is damaged, notify the supplier and the carrier in writing within three days; state the item number of the product together with the order number and the supplier's invoice number. Retain all packing material for inspection.

Comply with **national safety regulations and safety requirements** concerning the use of vacuum and electrical equipment. Comply with all **applicable and relevant safety requirements** (regulations and guidelines), **implement the required actions and adopt suitable safety measures**.

! DANGER

➤ The in-line valve (230 V~) must be connected only to a **suitable electrical supply** and a suitable earth point. Failure to connect to ground may result in deadly electrical shock.

➤ Adopt suitable measures to prevent the release of **dangerous, toxic, explosive, corrosive, noxious or polluting fluids, vapours and gases**.

➤ This device is not approved for operation in potentially explosive atmospheres.

➤ Valves without the "Ex" mark on the rating plate of the control unit are not approved for operation with dangerous or explosive gases or with potentially explosive or inflammable substances. Do not operate the valve with dangerous or explosive gases or with potentially explosive or inflammable substances.

➤ VACUU•BUS valves bearing the "Ex" mark on the rating plate of the control unit are approved for operation with potentially explosive atmospheres according to their ATEX classification imprinted on their rating plate, but they are **not approved for operation in potentially explosive atmospheres** (see section „Ex Important information: Equipment marking (ATEX)“, pg. 16).

➤ **Note:** The equipment may be contaminated with chemicals, which have been processed during operation. Ensure that the equipment is completely decontaminated before maintenance commences. Take adequate precautions to protect people from the effects of dangerous substances if contamination has occurred. Use appropriate protective clothing, safety goggles and protective gloves.

! WARNING

📖 **Read and comply with this manual.**

📖 **Do not use the equipment if it is damaged.**

📖 **Do not disassemble any valve if under pressure or vacuum!**

📖 **Use only OEM spare parts and accessories.**

! CAUTION

- Check that mains voltage and current conform with the equipment (see rating plate).
- **If idle, the in-line valves are closed. If energised, the in-line valves are open.**
- Comply with max. permissible temperatures and pressures (see "Technical data").
- Check that the system to be evacuated and all hose connections are mechanically stable and that all fittings are secure.
- Check that the connections between in-line valve and tubings are leak-tight. Check the valve diaphragm regularly for leak-tightness and correct seating, especially prior to initial operation.
- Ensure that the materials of the wetted parts are compatible with the pumped substances, see section "Technical data".

NOTICE

Check that the gas is flowing in the correct direction (shown by the arrow on the valve housing).

Ensure that maintenance is done only by a suitably trained and supervised technician. In order to comply with law (occupational, health and safety regulations, safety at work law and regulations for environmental protection) components returned to the manufacturer can be repaired only when certain procedures are followed.

 **Important information: Equipment marking (ATEX)**

Only valid for products with ATEX marking. If the ATEX marking is shown on the rating plate of the respective product, VACUUBRAND GMBH + CO KG assures, that the device complies with the provisions of the directive 2014/34/EU. The applied harmonized standards are indicated in the EC Declaration of Conformity of the Machinery (see instructions for use).

**VACUUBRAND equipment bearing the ATEX mark
(see rating plate)**

The classification according to ATEX is only valid for the inner part (wetted part, pumped gas or vapor) of the equipment. The equipment is not suitable for use in external, potentially explosive atmospheres (environment).

The overall category of the equipment depends on the connected components. If the connected components do not comply with the classification of the VACUUBRAND equipment, the specified category of the VACUUBRAND equipment is no longer valid.

Vacuum pumps and vacuum gauges in category 3 are intended for connection to equipment in which during normal operation explosive atmospheres caused by gases, vapors or mists normally don't occur; or, if they do occur, are likely to do so only infrequently and for a short period only.

Equipment in this category ensures the requisite level of protection during normal operation.

The use of gas ballast or the operation of venting valves is only permitted if thereby explosive atmospheres normally don't occur in the interior of the equipment or, if they do occur, are likely to do so only infrequently and for a short period.

The pumps are marked with "X" (according to DIN EN ISO 80079-36:2016), i.e., restrictions of the operation conditions:

- The equipment is designated for a low degree of mechanical stress and has to be installed in a way so that it cannot be damaged from outside.
Pumping units have to be installed so that they are protected against shocks from the outside and against glass splinters in the event of breakage (implosion).
- The equipment is designated for an ambient and gas inlet temperature during operation of +10 to +40°C. Never exceed these ambient and gas inlet temperatures. If pumping / measuring gases which are not potentially explosive, extended gas inlet temperatures are permissible. See instructions for use, section "Gas inlet temperatures" or "Technical data".

After any intervention at the equipment (e.g., repair / maintenance) the ultimate vacuum of the pump has to be checked. Only if the pump achieves its specified ultimate vacuum is the pump's leak rate low enough to ensure that no explosive atmospheres will occur in the interior of the equipment.

After any intervention at the vacuum sensor, the leak rate of the equipment has to be checked.

Attention: This manual is not available in all languages of the EU. The user must not operate the device if he does not understand this manual. In this case a technically correct translation of the complete manual has to be available. The manual must be completely read and understood before operation of the device. All specified measures must be applied, or else must be replaced by equivalent measures at the user's own risk.

Technical Data

Type	VV 6 VV-B 6	VV 6C VV-B 6C	VV 15C VV-B 15C	VV 25
Max. permissible pressure range	0 - 1.5 bar	0 - 1.5 bar	0 - 1 bar	0 - 1 bar
Leak rate	1*10 ⁻⁵ mbar*l/s	1*10 ⁻² mbar*l/s	1*10 ⁻² mbar*l/s	1*10 ⁻⁴ mbar*l/s
Max. permissible gas temperature*	80 °C	100 °C	100 °C	80 °C
Max. permissible range of ambient temperature	0 - 50 °C	0 - 40 °C	0 - 40 °C	0 - 50 °C
Electrical supply	24 V= or** 230 V~ 50/60 Hz	24 V=	24 V=	230 V~ 50/60 Hz
Power draw (operation) 24 V= 230 V~ 50/60 Hz	8 W 8 W / 15 VA	6 W -	14 W -	- 16 W / 32 VA
Degree of protection IEC 60529	IP 65			
Max. switching frequency per minute	120	50	50	120
Connection	small flange KF DN 16 or hose nozzle DN 6/10 mm		small flange KF DN 16 or DN 25	small flange KF DN 25
Nominal width	6 mm		15 mm	25 mm
Weight approx.	0.53 kg	0.35 kg	1.2 kg	1.4 kg

* if processing potentially explosive atmospheres (only VV-B 6 / VV-B 6C / VV-B 15C) : +10°C to +40°C

** 24V= : Cat.-No. 20674090 / 20674290

230 V~ 50/60 Hz: Cat.-No. 20674094

Wetted parts

Component	VV 6 VV-B 6	VV 6C VV-B 6C	VV 15C VV-B 15C	VV 25
Housing	PP	PVDF	PVDF	brass, nickel plated
Diaphragm	FKM	PTFE	PTFE	FKM
Valve seal	FKM	fluoroelastomer	PTFE	FKM
Hose nozzle	PPS	PPS	-	-
Small flange	stainless steel			

We reserve the right for technical modification without prior notice!

Spare parts

VV 6 / VV-B 6	Quantity	Cat.-No.
Seal ring 1.5 mm, R 1/4"	2	20639729
Small flange KF DN 16	2	20642462
Hose nozzle G 1/4 6/10 mm with O-ring 11x3	2	20642470
Solenoid-operated valve 230 V~ 50/60 Hz	1	23120361
Solenoid-operated valve 120 V~ 60 Hz	1	23120360

VV 6C / VV-B 6C	Quantity	Cat.-No.
Seal ring 1.5 mm, R 1/4"	2	20639729
Small flange KF DN 16	2	20642462
Hose nozzle G 1/4 6/10 mm with O-ring 11x3	2	20642470
Solenoid-operated valve 24 V= complete	1	20638424
O-ring 25 x 2	1	23120808
Valve seal (PTFE, white)	1	on demand
Valve seal (fluoroelastomer, black)	1	on demand

VV 15C / VV-B 15C	Quantity	Cat.-No.
Seal ring 1.5 mm, G 1/2"	2	20638882
Small flange KF DN 16	2	20643809
Small flange KF DN 25	2	20638817
Solenoid-operated valve 24 V=	1	20638881

VV 25	Quantity	Cat.-No.
Small flange KF DN 25	2	20640660

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Hereby the manufacturer declares that the device is in conformity with the directives:

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Alfred-Zippe-Str. 4 · 97877 Wertheim

T +49 9342 808-0 · F +49 9342 808-5555

info@vacuubrand.com - www.vacuubrand.com

VACUUBRAND®

DECLARATION OF CONFORMITY – China RoHS 2

VACUUBRAND GMBH + CO KG has made reasonable efforts to ensure that hazardous materials and substances may not be used in its products.

In order to determine the concentration of hazardous substances in all homogeneous materials of the subassemblies, a “Product Conformity Assessment” (PCA) procedure was performed. As defined in GB/T 26572 the “Maximum Concentration Value” limits (MCV) apply to these restricted substances:

- Lead (Pb): 0.1%
- Mercury (Hg): 0.1%
- Cadmium (Cd): 0.01%
- Hexavalent chromium (Cr(+VI)): 0.1%
- Polybrominated biphenyls (PBB): 0.1%
- Polybrominated diphenyl ether (PBDE): 0.1%

Environmentally Friendly Use Period (EFUP)

EFUP defines the period in years during which the hazardous substances contained in electrical and electronic products will not leak or mutate under normal operating conditions. During normal use by the user such electrical and electronic products will not result in serious environmental pollution, cause serious bodily injury or damage to the user’s assets. The Environmentally Friendly Use Period for VACUUBRAND products is 40 years.

MATERIAL CONTENT DECLARATION FOR VACUUBRAND PRODUCTS						
有毒有害物质或元素 Hazardous substances						
部件名称 Part name	铅 Pb	汞 Hg	镉 Cd	六价铬 Cr(+VI)	多溴联苯 PBB	多溴二苯醚 PBDE
包装 Packaging	0	0	0	0	0	0
塑料外壳 / 组件 Plastic housing / parts	0	0	0	0	0	0
真空油 Vacuum oil	0	0	0	0	0	0
电池 Battery	0	0	0	0	0	0
玻璃 Glass	X	0	0	0	0	0
电子电气组件 Electrical and electronic parts	X	0	0	0	0	0
控制器 / 测量设备 Controller / measuring device	X	0	0	0	0	0
金属外壳 / 组件 Metal housing / parts	X	0	0	0	0	0
电机 Motor	X	0	0	0	0	0
配件 Accessories	X	0	0	0	0	0
此表格是按照SJ/T 11364-2014中规定所制定的。 This table is created according to SJ/T 11364-2014.						

Declaration of Conformity – China RoHS 2**V5_September 2022**

Copyright 2022

VACUUBRAND®

- O: 表示该有毒有害物质在该部件所有均质材料中的含量均在GB/T 26572规定的限量要求以下。
O: Indicates that the above mentioned hazardous substance contained in all homogeneous materials of the part is below the required limit as defined in GB/T 26572.
- X: 表示该有毒有害物质至少在该部件某一均质材料中的含量超出GB/T 26572规定的限量要求。
X: Indicates that the above mentioned hazardous substance contained in at least one of the homogeneous materials of this part is above the required limit as defined in GB/T 26572.

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Batteries, glassware and accessories might not be content of the enclosed device and may have its own EFUP-marking and/or might be maintaining parts with changing EFUP-marking.

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Apart from the disclosures in the above table, the subassemblies are not intentionally manufactured or formulated with lead (Pb), mercury (Hg), cadmium (Cd), hexavalent chromium (Cr+VI), polybrominated biphenyls (PBB), and polybrominated diphenyl ethers (PBDE).

Products manufactured by VACUUBRAND may enter into further devices (e.g., rotary evaporator) or can be used together with other appliances (e.g., usage as booster pumps).

With these products and appliances in particular, please note the EFUP labeled on these products.

VACUUBRAND will not take responsibility for the EFUP of those products and appliances.

Place, date: Wertheim, 06 September 2022

(Dr. Constantin Schöler)
Managing Director

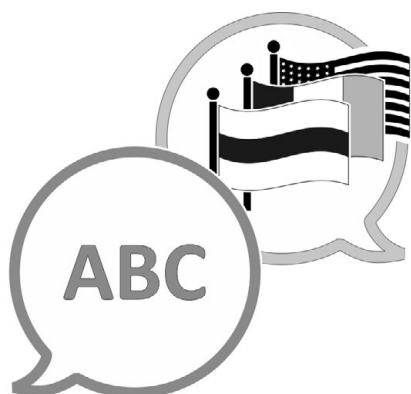
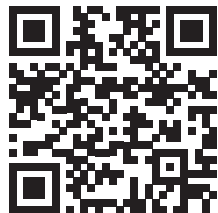
ppa. (Jens Kaibel)
Technical Director

VACUUBRAND GMBH + CO KG

Alfred-Zippe-Str. 4
97877 Wertheim
Germany

Tel.: +49 9342 808-0
Fax: +49 9342 808-5555
E-Mail: info@vacuubrand.com
Web: www.vacuubrand.com

[VACUUBRAND > Support > Manuals](#)

Manufacturer:

VACUUBRAND GMBH + CO KG
Alfred-Zippe-Str. 4
97877 Wertheim
GERMANY

Phone:

Head office +49 9342 808-0
Sales +49 9342 808-5550
Service +49 9342 808-5660

Fax: +49 9342 808-5555

Email: info@vacuubrand.com

Web: www.vacuubrand.com